

# Minitom



**Small, automatic precision cut-off machine for sectioning all materialographic and ceramic specimens**



*The transparent protection screen allows visual examination during the cutting process without the risk of splashing*

- **Simple operation**
- **Powerful motor**
- **High precision**
- **Functional design**
- **Easy to clean**

Minitom is a small, low-speed, precision cut-off machine for sectioning all types of materials. Maximum specimen size: 30 mm dia. The cutting speed is continuously variable and the motor is designed to ensure that the selected speed remains constant at any load. The required specimen size can be set with an accuracy of + 0.01 mm, by means of a micrometer screw, and the cutting pressure can be selected independent of the weight of the specimen. The motor is equipped with a thermal overload protection although prior to actually switching out an overload-indicator provides a visual warning that either the cutting procedure should be stopped or the cutting pressure should be reduced. Minitom can be adjusted to stop cutting at any stage of the operation, e.g. when the material has just been sliced through. The cutting process is liquid-cooled

with the cooling liquid draining back continuously into the coolant tank.

## Operation

Minitom is very easy to operate, with fixation of the specimen and the setting of specimen size, cutting speed, and cutting pressure taking only a few minutes. The workpiece is fastened in a holder mounted on a movable arm. This arm acts as a steelyard with two weights, one of which balances the specimen holder with specimen, the other being used to set the exact cutting pressure. During the cutting operation the arm is guided down over the cut-off wheel. When the cutting operation is finished the cut-off wheel stops and the specimen falls into a basket in the coolant tank thus protecting the specimen from drying and making retrieval easy. The coolant tank is very easy to remove and both the retrieval of the finished specimens and the cleaning of the coolant tank are therefore simple and straightforward. Minitom's practical and attractive design is the product of a very extensive programme of design development.

*Minitom is small, compact and extremely user-friendly*




## Specimen holders

Minitom can be supplied with a variety of specimen holders to suit a wide range of specimen types.

## Cut-off wheels

Minitom uses CBN and diamond cut-off wheels with a diameter of 75 mm (3") to 127 mm (5") and a center hole diameter of 12.7 mm (1/2"). The cut-off wheels are available with different grain size and concentration, thus ensuring that the correct cut-off wheel is available for any type of material.

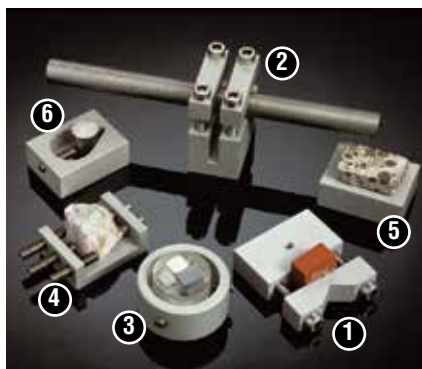

Minitom offers holders for all specimen shapes

## Technical specifications

	Voltage: 100-240 V / 50-60 Hz Output: 38 W
<b>Motor shaft</b>	Speed: 100-420 rpm Diameter: 12.7 mm (1/2") Wheel diameter: 75-127 mm (3"-5")
<b>Specimen holder arm</b>	Balances specimen holder with workpieces up to 200 g in weight Weight 0-350 g, which is used to set the cutting pressure Axial movement: 0-25 mm
<b>Coolant tank</b>	Volume: 250 ml
<b>Dimensions</b>	Width: 280 mm Depth: 400 mm Height: 400 mm Weight: 8 kg

Struers' equipment is in conformity with the provisions of the applicable International Directives and their appurtenant Standards. (Please contact your local supplier for details)

## Specifications

### Minitom

Low-speed precision cut-off machine. Complete with universal specimen holder (04436901) and flange set 65 mm dia. as well as 42 mm dia. Cut-off wheels are ordered separately. 100-240 V / 50-60 Hz

	Cat. no.
<b>1 Specimen holders</b> For general use. Vice type with max. opening 25 mm / 1".	04436901
<b>2</b> For long specimens. Double parallel vice with max. opening 25 mm / 1".	04436902
<b>3</b> For round specimens. Max. opening 30 mm dia.	04436903
<b>4</b> For irregularly-shaped specimens. With 5 screws. Max. opening 30 mm.	04436904
<b>5</b> For adhesive mounting of specimens. 30 x 40 mm.	04436905
<b>6</b> For round specimens. Teardrop type with max. opening 20 mm dia.	04436906

Struers' products are subject to constant product development. Therefore, we reserve the right to introduce changes in our products without notice.

### Struers ApS

Pederstrupvej 84  
DK-2750 Ballerup, Denmark  
Phone +45 44 600 800  
Fax +45 44 600 801  
struers@struers.dk  
www.struers.com

### AUSTRALIAN & NEW ZEALAND

**Struers Australia**  
27 Mayneview Street  
Milton QLD 4064  
Australia  
Phone +61 7 3512 9600  
Fax +61 7 3369 8200  
info.au@struers.dk

### BELGIUM (Wallonia)

**Struers S.A.S.**  
370, rue du Marché Rollay  
F- 94507 Champigny  
sur Marne Cedex  
Téléphone +33 1 5509 1430  
Télécopie +33 1 5509 1449  
struers@struers.fr

### BELGIUM (Flanders)

**Struers GmbH Nederland**  
Zomerdijk 34 A  
3143 CT Maassluis  
Telefoon +31 (10) 599 7209  
Fax +31 (10) 5997201  
netherlands@struers.de

### CANADA

**Struers Ltd.**  
7275 West Credit Avenue  
Mississauga, Ontario L5N 5M9  
Phone +1 905-814-8855  
Fax +1 905-814-1440  
info@struers.com

### CHINA

**Struers Ltd.**  
No. 1696 Zhang Heng Road  
Zhang Jiang Hi-Tech Park  
Shanghai 201203, P.R. China  
Phone +86 (21) 6035 3900  
Fax +86 (21) 6035 3999  
struers@struers.cn

### CZECH REPUBLIC & SLOVAKIA

**Struers GmbH Organizační složka**  
v dekoltechnický park  
Plešská 1920,  
CZ-252 63 Roztoky u Prahy  
Phone +420 233 312 625  
Fax +420 233 312 640  
czechrepublic@struers.de  
slovakia@struers.de

### GERMANY

**Struers GmbH**  
Carl-Friedrich-Benz-Straße 5  
D- 47877 Willich  
Phone +49 (0) 2154 486-0  
Fax +49 (0) 2154 486-222  
verkauf@struers.de

### FRANCE

**Struers S.A.S.**  
370, rue du Marché Rollay  
F-94507 Champigny  
sur Marne Cedex  
Téléphone +33 1 5509 1430  
Télécopie +33 1 5509 1449  
struers@struers.fr

### HUNGARY

**Struers GmbH**  
Magyarország Fióktelep  
Tatai út 53  
2821 Gyermely  
Phone +36 (34) 880546  
Fax +36 (34) 880547  
hungary@struers.de

### IRELAND

**Struers Ltd.**  
Unit 11 Evolution@ AMP  
Whittle Way, Catcliffe  
Rotherham S60 5BL  
Tel. +44 0845 604 6664  
Fax +44 0845 604 6651  
info@struers.co.uk

### ITALY

**Struers Italia**  
Via Monte Grappa 80/4  
20020 Arese (MI)  
Tel. +39-02/38236281  
Fax +39-02/38236274  
struers.it@struers.it

### JAPAN

**Marumoto Struers K.K.**  
Takanawa Muse Bldg. 1F  
3-14-13 Higashi-Gotanda,  
Shinagawa  
Tokyo  
141-0022 Japan  
Phone +81 3 5488 6207  
Fax +81 3 5488 6237  
struers@struers.co.jp

### NETHERLANDS

**Struers GmbH Nederland**  
Zomerdijk 34 A  
3143 CT Maassluis  
Telefoon +31 (10) 599 7209  
Fax +31 (10) 5997201  
netherlands@struers.de

### NORWAY

**Struers ApS, Norge**  
Sjøskogveien 44C  
1407 Vinterbro  
Telefon +47 970 94 285  
info@struers.no

### AUSTRIA

**Struers GmbH**  
Zweigniederlassung Österreich  
Betriebsgebiet Puch Nord 8  
5412 Puch  
Telefon +43 6245 70567  
Fax +43 6245 70567-78  
austria@struers.de

### POLAND

**Struers Sp. z o.o.**  
Oddział w Polsce  
ul. Jasnogórska 44  
31-358 Kraków  
Phone +48 12 661 20 60  
Fax +48 12 626 01 46  
poland@struers.de

### ROMANIA

**Struers GmbH, Sucursala Bucuresti**  
Str. Preciziei nr. 6R  
062203 sector 6, Bucuresti  
Phone +40 (31) 101 9548  
Fax +40 (31) 101 9549  
romania@struers.de

### SWITZERLAND

**Struers GmbH**  
Zweigniederlassung Schweiz  
Weissenbrunnstrasse 41  
CH-8903 Birmensdorf  
Telefon +41 44 777 63 07  
Fax +41 44 777 63 09  
switzerland@struers.de

### SINGAPORE

**Struers Singapore**  
627A Aljunied Road,  
#07-08 BizTech Centre  
Singapore 389842  
Phone +65 6299 2268  
Fax +65 6299 2661  
struers.sg@struers.dk

### SPAIN

**Struers España**  
Camino Cerro de los Gamos 1  
Building 1 - Pozuelo de Alarcón  
CP 28224 Madrid  
Teléfono +34 917 901 204  
Fax +34 917 901 112  
struers.es@struers.es

### FINLAND

**Struers ApS, Suomi**  
Hietalahdenranta 13  
00180 Helsinki  
Puhelin +358 (0)207 919 430  
Faksi +358 (0)207 919 431  
finland@struers.fi

### SWEDEN

**Struers Sverige**  
Box 20038  
161 02 Bromma  
Telefon +46 (0)8 447 53 90  
Telefax +46 (0)8 447 53 99  
info@struers.se

### UNITED KINGDOM

**Struers Ltd.**  
Unit 11 Evolution @ AMP  
Whittle Way, Catcliffe  
Rotherham S60 5BL  
Tel. +44 0845 604 6664  
Fax +44 0845 604 6651  
info@struers.co.uk

### USA

**Struers Inc.**  
24766 Detroit Road  
Westlake, OH 44145-1598  
Phone +1 440 871 0071  
Fax +1 440 871 8188  
info@struers.com