DuroCit-3

#5020065

01.07.2016

Instructions

Anleitung

Mode d'emploi

English

Description

DuroCit-3 is a yellow three component acrylic mounting system with filler, consisting of one powder and two liquids. DuroCit-3 is used when good edge retention and good workability properties are required.

DuroCit-3 cures at room temperature.

Material Safety Data

Please read the following carefully:

DuroCit Powder

WARNING

Contains: Dicyclohexyl phthalate. Dibenzoyl peroxide.

H317: May cause an allergic skin reaction.

P261: Avoid breathing dust/fume/gas/mist/vapours/spray.

P280: Wear protective clothing and gloves.

P302/352: IF ON SKIN: Wash with plenty of soap and water.

P333/313: If skin irritation or rash occurs: Get medical advice/attention.

P363: Wash contaminated clothing before reuse.

P501: Dispose of contents/container in accordance with local regulations.

DuroCit-3 Liquid I

Danger

Contains: Methyl 2-methylprop-2-enoate. Styrene.

H226: Flammable liquid and vapour.

H315: Causes skin irritation.

H317: May cause an allergic skin reaction.

H319: Causes serious eye irritation.

H361d: Suspected of damaging the unborn child.

H372H: Causes damage to organs (hearing organs) through prolonged or repeated exposure.

P210: Keep away from heat/sparks/open flames/hot surfaces. – No smoking.

P280: Wear protective clothing, gloves, eye and face protection.

P260: Do not breathe gas, fume, vapours or spray.

P302/352: IF ON SKIN: Wash with plenty of soap and water.

P305/351/338 IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

P337/313: If eye irritation persists: Get medical advice/attention.

DuroCit-3 Liquid II

Danger

Contains: Methyl 2-methylprop-2-enoate. Styrene.

H225: Highly flammable liquid and vapour.

H315: Causes skin irritation.

H317: May cause an allergic skin reaction.

H335: May cause respiratory irritation.

H361d: Suspected of damaging the unborn child.

H373H: May cause damage to organs (hearing organs) through prolonged or repeated exposure.

P210: Keep away from heat/sparks/open flames/hot surfaces. – No smoking.

P280: Wear protective clothing, gloves, eye and face protection.

P260: Do not breathe gas, fume, vapours or spray..

P302/352: IF ON SKIN: Wash with plenty of soap and water.

P501: Dispose of contents/container in accordance with local regulations.

Mounting of Specimens

Measuring of Components

The ratio for DuroCit Powder:DuroCit-3 Liquid I: DuroCit-3 Liquid II is **14:8:4** by weight and **15:10:5** by volume.

Measuring of Components – premixed liquid

Ratio for Durocit Powder:Premixed liquid is 14:12 by weight and 15:15 by volume.

Weighing is recommended to obtain the best mounts and for reproducibility.

Important!

Do NOT pour excess liquid back into the bottle as this will significantly reduce the shelf-life.

Mixing of Components

- Measure out the amount of DuroCit-3 Liquid I into a mixing cup (max. 32 g / 30 ml).
- Measure out the amount of DuroCit Powder into a mixing cup (max. 56 g / 45 ml).
- Pour the Powder into Liquid I and stir thoroughly for at least 1-2 min., until the mixture is homogeneous and without lumps.
- Carefully add the correct amount of Liquid II into the Liquid I and Powder mixture.
- Stir thoroughly for 1 min. Mix carefully to avoid trapping air in the mixture.

Mixing of Components – premixed liquid

Alternatively liquid I and liquid II can be premixed in the correct ratio in advance. Storing it for 5°C the mixture will keep for 4 weeks.

- Measure out the amount of premixed liquid into a mixing cup (max. 48 g / 45 ml).
- Measure out the amount of DuroCit Powder into a mixing cup (max. 56 g / 45 ml).
- Pour the Powder into the premixed liquid and stir thoroughly for at least 1-2 min., until the mixture is homogeneous and without lumps.

For best results, use the stirrer to press the mixture against the wall of the mixing cup to break any lumps.

Selecting the Mounting Cup

All Struers mounting cups can be used. FlexiForm is recommended for easy unmounting of the mount. If possible, choose a mounting cup where the distance between the specimen and cup wall is approx. 5 mm/0.2".

Pre-treatment of specimens

To promote the best possible adhesion of DuroCit-3 to the specimen ensure that the specimen is thoroughly cleaned and dry prior to mounting. Each surface must be free from grease, moisture and other contaminants.

Mounting

- Place the cleaned specimen in the centre of the mounting cup.
- Mix the required amount of DuroCit-3.
- Pour the mixture carefully over the specimen; run the mixture slowly along the stirrer or the mixing cup handle.
- Ensure that no air remains trapped within cracks or cavities in the specimen by moving the specimen or by tapping the mounting cup on the table.

Curing

DuroCit-3 is a room temperature curing system. For a 30 mm dia. mount, the curing time is about 30 minutes. We recommend keeping the mount in the mounting cup until the temperature is below 40°C/104°F.

After Curing

After removal from the mounting cup, the mount is immediately ready for preparation.

Technical Data

	Components	Ratio by weight	Ratio by volume
DuroCit-3	DuroCit Powder	14	15
	DuroCit-3 Liquid I	8	10
	DuroCit-3 Liquid II	4	5
	Premixed liquid	12	15

NB For the best mount use weighing method.

Density

DuroCit Powder: 1.26 g/ml
 DuroCit-3 Liquid I: 1.05 g/ml
 DuroCit-3 Liquid II: 1,10 g/ml
 Premixed Liquid 1.07 g/ml

Soluble in DuroCit-3 Liquid I and II are soluble in Acetone

Filler The filling material in DuroCit is a form of aluminium silicate.

Curing

Can be prepared (ground and polished) after:	30 min.
Peak temperature *:	138°C / 280 °F
Pot life:	4 min.

*30 mm dia. mounting cup without specimen at 21°C / 70°F.

Properties after Curing

Hardness: 85 Shore D
 Density: 1,6 g/ml
 Chemical resistance: DuroCit-3 is resistant to the most common etching materials.

Tips

- It can be an advantage to prolong the first polishing step a little to get the optimal planeness between specimen and DuroCit-3 mounting material.
- DuroCit-3 can be cured in a pressure chamber to minimize possible air trapped in the mount.

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