# RotoPol-21

# RotoPol-22

# RotoPol-25

Instruction Manual

Manual No.: 15157001

Date of Release 15.03.2002

<b>Table of Contents</b>	<b>Page</b>
User's Guide .....	1
Reference Guide.....	21
Quick Reference Guide.....	35

---

Always state *Serial No* and *Voltage/frequency* if you have technical questions or when ordering spare parts. You will find the Serial No. and Voltage on the type plate of the machine itself. We may also need the *Date* and *Article No* of the manual. This information is found on the front cover.

The following restrictions should be observed, as violation of the restrictions may cause cancellation of Struers legal obligations:

**Instruction Manuals:** Struers Instruction Manual may only be used in connection with Struers equipment covered by the Instruction Manual.

**Service Manuals:** Struers Service Manual may only be used by a trained technician authorised by Struers. The Service Manual may only be used in connection with Struers equipment covered by the Service Manual.

Struers assumes no responsibility for errors in the manual text/illustrations. The information in this manual is subject to change without notice. The manual may mention accessories or parts not included in the present version of the equipment.

The contents of this manual are the property of Struers. Reproduction of any part of this manual without the written permission of Struers is not allowed.

All rights reserved. © Struers 2002.

**Struers A/S**  
Valhøjs Allé 176  
DK-2610 Rødovre/Copenhagen  
Denmark  
Telephone +45 36 70 35 00  
Fax +45 38 27 27 01

---

## **RotoPol -21/-22/-25**

### **Safety Precaution Sheet**

#### **To be read carefully before use**

1. The operator(s) should be fully instructed in the use of the machine, any connected equipment and accessories and the applied consumables according to the relevant Instruction Manuals.
2. The machine must be placed on a safe and stable table with an adequate working height. All functions on the machine and any connected equipment must be in working order.
3. Operators should ensure that the actual voltage corresponds to the voltage on the back of the machine. The machine must be earthed. Follow the local regulations. Always turn the power off and remove the plug or the cable before opening the machine.
4. Connect only to cold water. Make sure that the water connections are leak-proof and that the water outlet is working.
5. The connection to the mains water supply should be turned off if the machine is left unattended.
6. Establish efficient exhaust from the working area if consumables containing alcohol, for example, are being used.
7. Make sure that the grinding or polishing discs are correctly mounted, the driving pins engaged with the holes in the turntable.
8. Keep clear of the rotational disc and the specimen mover if these are mounted during operation. Make sure that any screws for clamping of specimens in specimen holders protrude as little as possible. While grinding manually, be careful not to touch the grinding paper. Do not drop the specimen while grinding.
9. If you observe malfunctions or hear unusual noises - stop the machine and call technical service.

---

The equipment is designed for use with consumables supplied by Struers. If subjected to misuse, improper installation, alteration, neglect, accident or improper repair, Struers will accept no responsibility for damage(s) to the user or the equipment.

Dismantling of any part of the equipment, during service or repair, should always be performed by a qualified technician (electromechanical, electronic, mechanical, pneumatic, etc.)

---

# User's Guide

Table of Contents	Page
<b>1. Getting Started</b>	
Checking Contents of Packing .....	3
RotoPol-21 .....	3
RotoPol-22 .....	3
RotoPol-25 .....	3
Recirculation Cooling Unit Optional (TRECI) .....	4
Recirculation Cooling Unit Connector, Optional (ROPRE)...	4
Getting Acquainted with RotoPol .....	5
Unpacking RotoPol .....	5
Placing RotoPol .....	5
Supplying Power .....	6
Direction of Turntable .....	6
Supplying Water .....	6
Connection to Water Mains .....	6
Connection to Water Outlet .....	7
Software Settings.....	7
Setting the Language .....	8
Language Setting for RotoPol-21 and -25 .....	8
Language Setting for RotoPol-22.....	9

## **2. Basic Operations**

Using the Controls.....	10
Front Panel of RotoPol-21.....	10
Front Panel of RotoPol-22.....	12
General Use of Keys.....	16
Reading the Display.....	16
Routine Operations.....	16
Mounting a Disc.....	16
Starting the Disc.....	16
Stopping the Disc.....	16
User Select on RotoPol-21 and -22.....	17
Selecting Water and Speed.....	17
Setting to User Select.....	17
Setting to GRINDING.....	17
Setting to POLISHING.....	17
User Select on RotoPol-25.....	18
Selecting Water.....	18
Setting to User Select.....	18
Setting to GRINDING.....	18
Setting to POLISHING.....	18
Settings with a RotoCom Memory Unit.....	19

## **1. Getting Started**

### **Checking Contents of Packing**

#### *RotoPol-21*

In the packing case you should find the following parts:

- 1 RotoPol-21
- 2 Splash rings
- 1 Inlet hose 3/4" x 2 m
- 1 Filter gasket
- 1 Gasket  $\varnothing 11/\varnothing 24$  x 1.5
- 1 Reduction ring with gasket
- 1 Outlet hose  $\varnothing 32$  x 3 m
- 1 Hose clamp
- 1 Template
- 1 Set of Instruction Manuals

#### *RotoPol-22*

- 1 RotoPol-22
- 2 Splash rings
- 1 Inlet hose 3/4" x 2 m
- 1 Filter gasket
- 1 Gasket  $\varnothing 11/\varnothing 24$  x 1.5
- 1 Reduction ring with gasket
- 1 Outlet hose  $\varnothing 32$  x 3 m
- 1 Hose clamp
- 1 Template
- 1 Set of Instruction Manuals

#### *RotoPol-25*

- 1 RotoPol-25
- 1 Power cable
- 2 Splash rings
- 1 Inlet hose 3/4" x 2 m
- 1 Filter gasket
- 1 Gasket  $\varnothing 11/\varnothing 24$  x 1.5
- 1 Reduction ring with gasket
- 1 Outlet hose  $\varnothing 32$  x 3 m
- 1 Hose clamp
- 1 Template
- 1 Set of Instruction Manuals

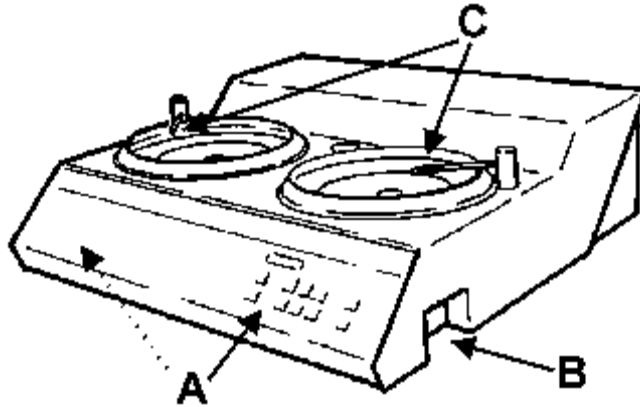
*RotoPol-21/-22/-25*  
*Instruction Manual*

<i>Recirculation Cooling Unit</i>	1	TRECI complete
<i>Optional (TRECI)</i>	1	Outlet angle
	1	Funnel
	2	Disposable plastic containers (TREPO)
<i>Recirculation Cooling Unit</i>	1	Connection box
<i>Connector, Optional (ROPRE)</i>	1	Elbow 90°
	1	Reduction piece, large
	1	Reduction piece, small
	1	Hose clamp
	1	Additive for Cooling Fluid, 1 l (ADDUN)
	1	Diagram


**Getting Acquainted with RotoPol**

Take a moment to familiarise yourself with the location and names of all the RotoPol components:

- A Front panel control(s)
- B Main switch
- C Water taps with sprinkler tube and turntables

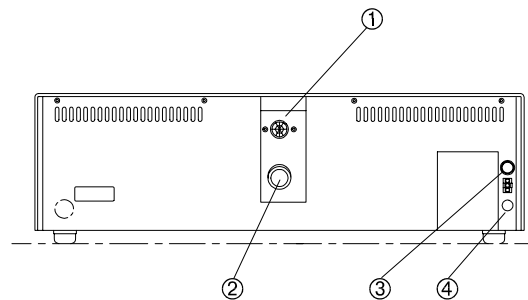*Back of RotoPol*

- ① Water inlet tube
- ② Water outlet tube
- ③ Socket for specimen mover
- ④ Socket for recirculation cooling unit

**Unpacking RotoPol**

All RotoPol models are detached from the bottom of the packing case by removing the four screws from below.

**Placing RotoPol**

RotoPol should be placed on a steady table with an adequate working height. The machine must be close to the power supply, water mains and water outlet facilities. If water recirculation is used, there must be room under the table for the Recirculation Cooling Unit (TRECI).

## Supplying Power

**IMPORTANT**  
Check that the mains voltage corresponds to the voltage stated on the type plate on the side of the machine.

### *RotoPol-21 and -22*

RotoPol-21 and -22 are factory mounted with electric cable. Mount a plug on the cable or connect to mains:

Black and brown: phase.  
Yellow/green: earth.

### *RotoPol-25*

Plug the power cable into the rear panel of RotoPol-25 and connect the other end to a grounded electrical outlet.

## Direction of Turntable

### *RotoPol-21 and -22*

Check that the turntable rotates in the direction indicated by the label (anti-clockwise). If the direction of rotation is incorrect, switch two of the phases.

## Supplying Water

Water for wet grinding and disc cooling may be supplied from the water mains or a Recirculation Cooling Unit.

### *Connection to Water Mains*

- Mount the straight end of the pressure hose onto the water inlet tube (upper) in the hose box on the back of RotoPol-21/22/25 (see drawing on the previous page):
  - Insert the filter gasket in the coupling nut with the flat side against the pressure hose.
  - Tighten the coupling nut completely.
- Mount the other end of the pressure hose on the water mains tap for cold water:
  - Mount the reduction ring with gasket on the water mains tap, if necessary.
  - Introduce the gasket and tighten the coupling nut completely.

**IMPORTANT**  
Only connect to cold water.

*Connection to Water Outlet*

- Locate the end of the drain hose where a short length of the steel reinforcement has been stripped away. Mount the stripped end of the hose onto the water outlet tube. Use hose clamp for fastening.
- Lead the other end of the drain hose to the water outlet. Arrange the tube so that it slopes throughout its length. Shorten the hose, if necessary. The drain hose should be attached to the water outlet system if possible.

**REMEMBER**  
Avoid sharp bends in the drain hose and make sure that the hose slopes throughout its length.

**Software Settings**

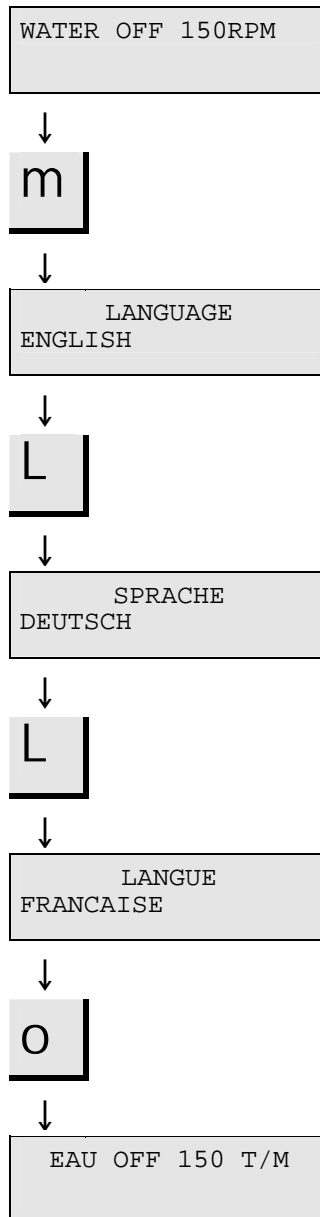
- Switch on the power at the main switch. The following display will appear briefly:

- RotoPol-2X -  
VERSION X.Y

- The operation menu will then appear:

WATER OFF 150RPM

**Setting the Language**  
*Language Setting for  
RotoPol-21 and -25*

Example screen (operation menu).

The MENU **m** key activates the setup menu.

Press SELECT **L** to choose English, German, French or Japanese.

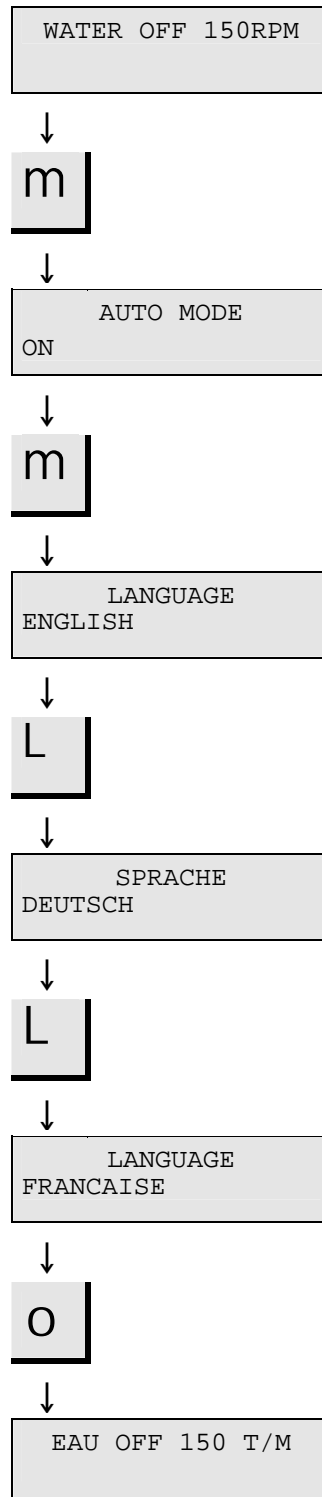
Press SELECT **L** until you reach the language you prefer.

Press STOP **O** to accept the choice. The Operation Menu will appear again in the language you have chosen.

Remember to set the same language on all connected RotoSystem modules.

Setting the Communication: see Advanced Operation Section.

Language Setting for  
RotoPol-22

Example screen

The MENU **m** key activates the setup menu.

Press **m** again. The language screen appears.

Press SELECT **L** to choose English, German, French or Japanese.

Press SELECT **L** until you reach the language you prefer.

Press STOP **O** to accept the choice. The Operation Menu will appear again in the language you have chosen.

- Repeat the above procedure for the next disc.



Remember to set the same language on all connected RotoSystem modules. Setting the Communication: see Advanced Operation Section.

## 2. Basic Operations
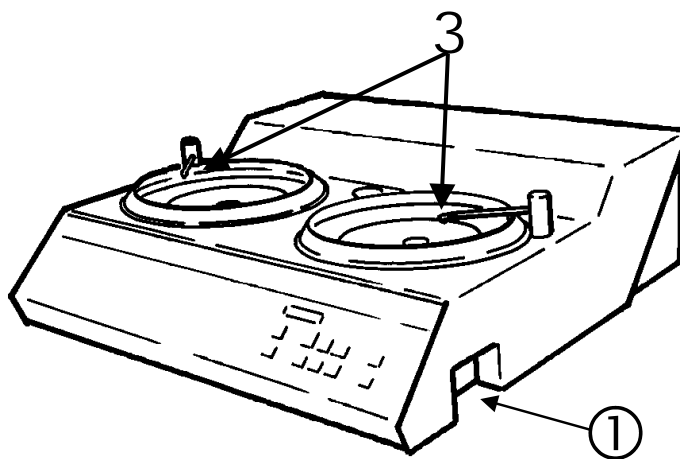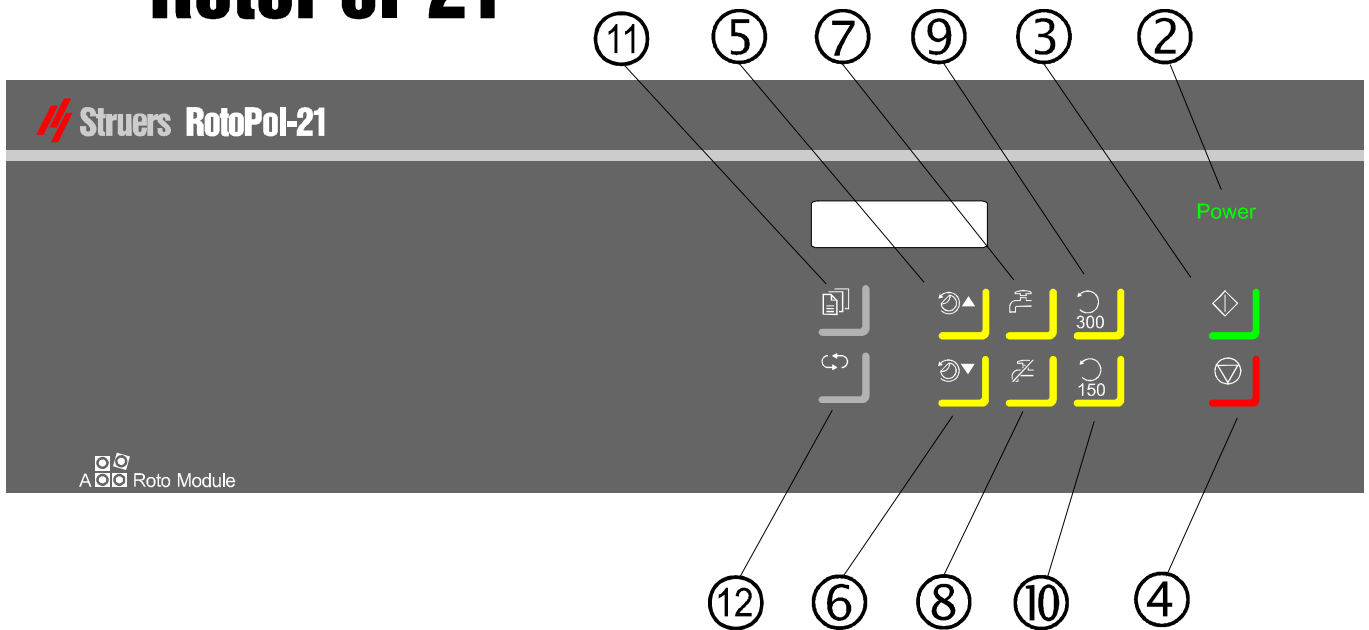
### Using the Controls

Front Panel of RotoPol-21


# RotoPol-21

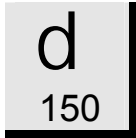*RotoPol-21/-22/-25  
Instruction Manual*


**Front Panel Controls for RotoPol-21**

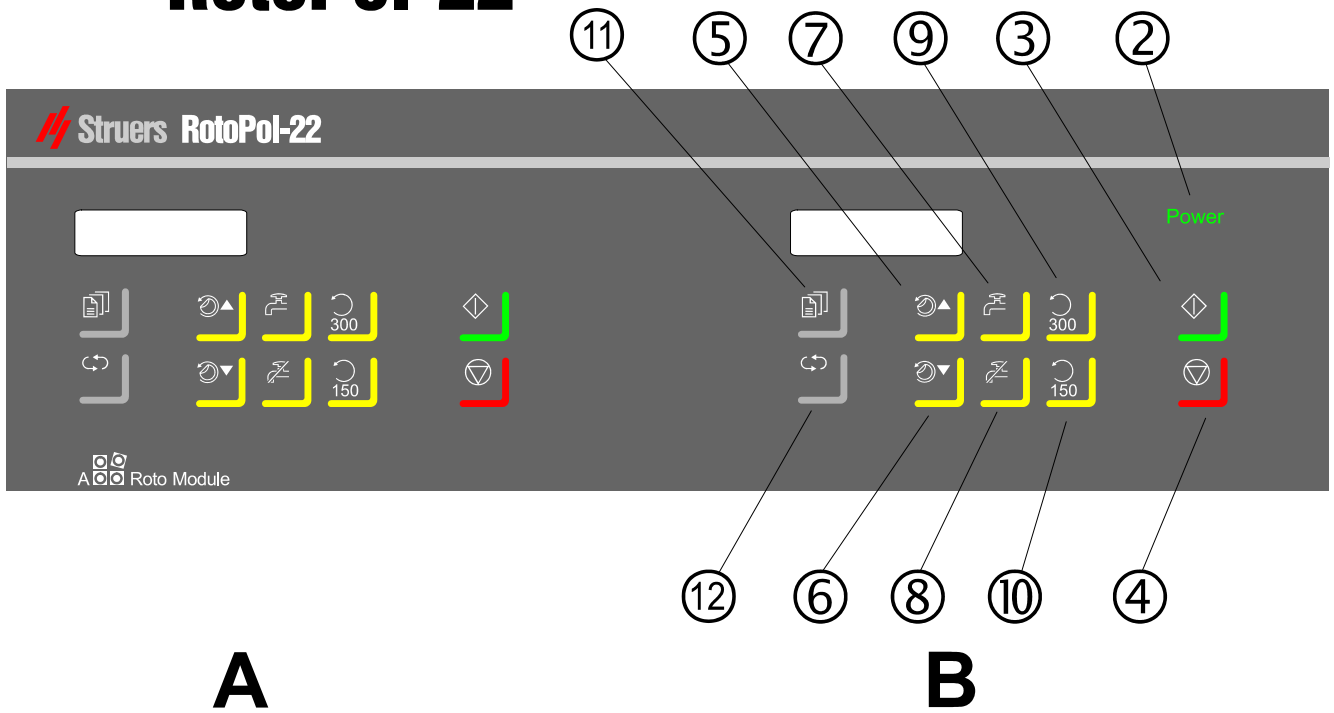*Communication is set to STANDARD mode*

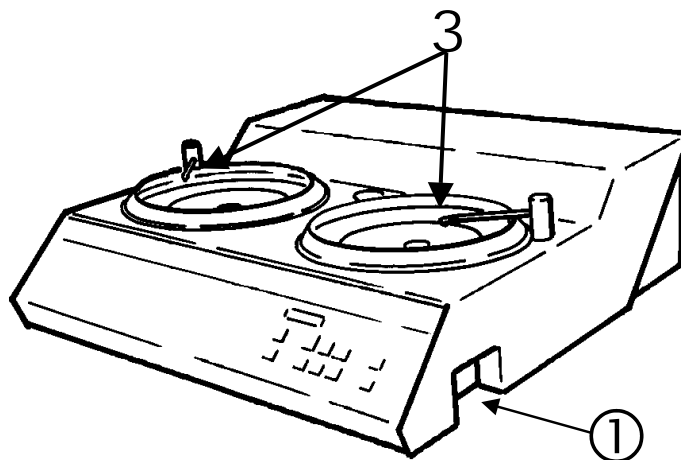
Name		Function	Name	Key	Function
① MAIN SWITCH		The main switch is located to the right of the cabinet. Turn clockwise to switch on the power.	⑧ WATER OFF		Select WATER OFF.
② POWER INDICATOR		Lights up when the main power is on.	⑨ SPEED		Select 300 RPM.
③ START		Starts the disc and any connected RotoSystem equipment set in Auto mode.	⑩ SPEED		Select 150 RPM.
④ STOP		Stops the disc and any connected RotoSystem equipment set in Auto mode. Exits the setup menu.	1 MENU		Activates the setup menu or moves down in the menu structure.
⑤ DISC		Scrolls up in the operation menu	2 SELECT		Changes values in the setup menu.
⑥ DISC		Scrolls down in the operation menu.	3 WATER TAP		Adjust the water at the tap.
⑦ WATER ON		Select WATER ON: press the key briefly. Manual Dosing: keep pressing the key. Release the key to stop the water.			

Front Panel of RotoPol-22

# RotoPol-22


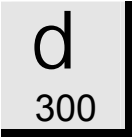
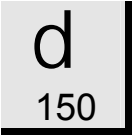
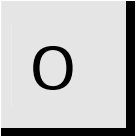
- A** Front panel controls for the disc on the left-hand side. The figure numbers are identical to the front panel controls for the right-hand disc.
- B** Front panel controls for the disc on the right-hand side.



**Front Panel Controls for RotoPol-22 (Identical for side A and B)**

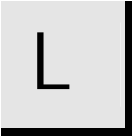*Communication is set to STANDARD mode*

Name		Function	Name	Key	Function
① MAIN SWITCH		The main switch is located to the right of the cabinet. Turn clockwise to switch on the power.	⑧ WATER OFF		Select WATER OFF.
② POWER INDICATOR		Lights up when the main power is on.	⑨ SPEED		Select 300 RPM.
③ START		Starts the disc and any connected RotoSystem equipment set in Auto mode.	⑩ SPEED		Select 150 RPM.
④ STOP		Stops the disc and any connected RotoSystem equipment set in Auto mode. Exits the setup menu.	1 MENU		Activates the setup menu or moves down in the menu structure.
⑤ DISC		Scrolls up in the operation menu	2 SELECT		Changes values in the setup menu.
⑥ DISC		Scrolls down in the operation menu.	3 WATER TAP		Adjust the water at the tap.
⑦ WATER ON		Select WATER ON: press the key briefly. Manual Dosing: keep pressing the key. Release the key to stop the water.			

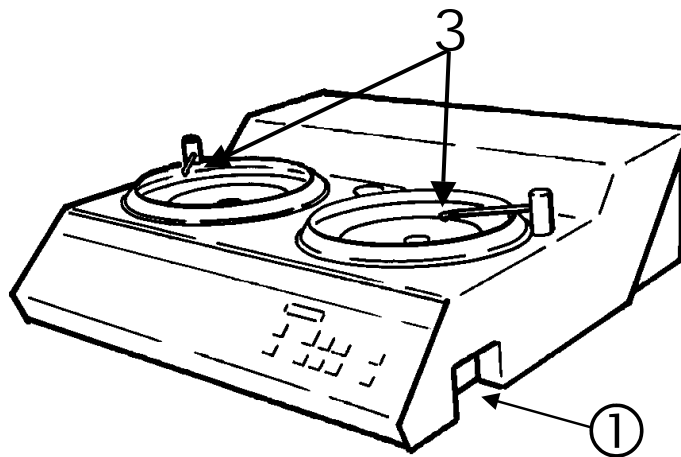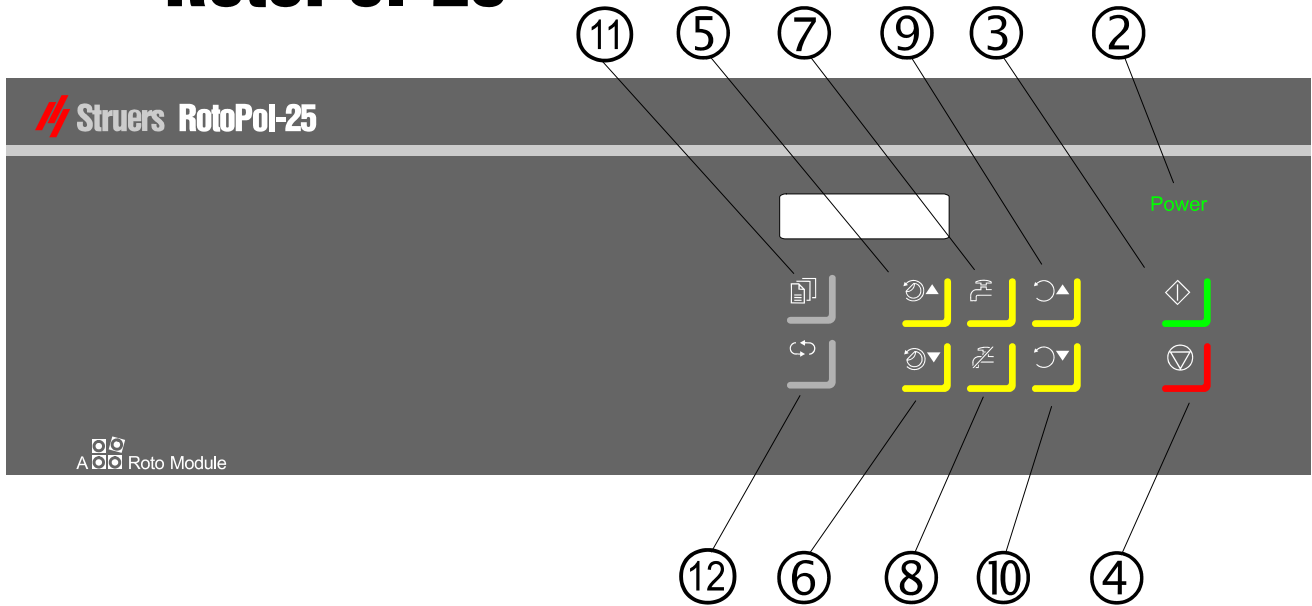
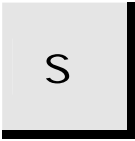
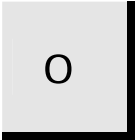
Front Panel of RotoPol-25

# RotoPol-25

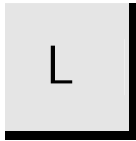**Front Panel Controls for RotoPol-25**

*Communication is set to STANDARD mode*

Name		Function	Name	Key	Function
① MAIN SWITCH		The main switch is located to the right of the cabinet. Turn clockwise to switch on the power.	⑧ WATER OFF		Select WATER OFF.
② POWER INDICATOR		Lights up when the main power is on.	⑨ SPEED		Increase the speed
③ START		Starts the disc and any connected RotoSystem equipment set in Auto mode.	⑩ SPEED		Decrease the speed
④ STOP		Stops the disc and any connected RotoSystem equipment set in Auto mode. Exits the setup menu.	1 MENU		Activates the setup menu or moves down in the menu structure.
⑤ DISC		Scrolls up in the operation menu	2 SELECT		Changes values in the setup menu.
⑥ DISC		Scrolls down in the operation menu.	3 WATER TAP		Adjust the water at the tap.
⑦ WATER ON		Select WATER ON: press the key briefly. Manual Dosing: keep pressing the key. Release the key to stop the water.			

## **General Use of Keys**

- WATER ON **W**, WATER OFF **W** and SPEED **d** can only be selected freely at User Select.
- MENU **m**, SELECT **L** and DISC **A** can only be activated when the machine has been stopped.
- SELECT **L** can only be activated when MENU **m** is selected.

## Acoustic Signals

- When pressing a key, a short beep indicates that the command has been accepted, whereas a long beep indicates that the key cannot be activated at the moment.

## **Reading the Display**

WATER ON 300RPM GRINDING
-----------------------------

- The first line on the screen shows the setting of parameters WATER and SPEED.
- The second line, if any, shows a selected preset mode (GRINDING or POLISHING).

## **Routine Operations**

### *Mounting a Disc*

- Place the disc on the turntable and turn it until the three driving pins engage with the holes in the turntable.

### *Starting the Disc*

- Press START **S** . The disc starts rotating. Selected functions are executed, and connected RotoSystem modules start operating.

### *Stopping the Disc*

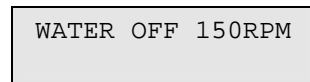
- Press STOP **O** . The disc stops rotating. Selected functions and connected RotoSystem modules stop.

**User Select on  
RotoPol-21 and -22**

*Selecting Water and Speed*

You can either select water and speed yourself or press DISC A to select one of the preset GRINDING and POLISHING modes.

*Setting to User Select*

The user Select screen has an empty line at the bottom.

- Press W or **W** to select WATER ON or WATER OFF.

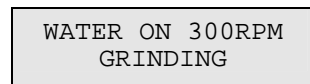
- Press SPEED **d** 150 or **d** 300 to select speed

Press DISC A N or A O to change between the settings.

*Setting to GRINDING*


The standard setting GRINDING is preset to WATER ON and 300RPM.

Press DISC A N or A O to change between the settings.

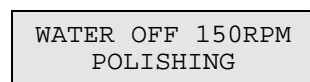*Setting to POLISHING*

The standard setting POLISHING is preset to WATER OFF and 150RPM.

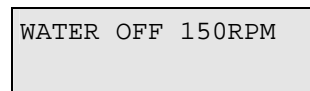
**User Select on  
 RotoPol-25**

*Selecting Water*

You can either select water yourself or press DISC A N to select one of the preset GRINDING and POLISHING modes.

RotoPol-25 is protected against unintentional cooling water supply on the surface when running in POLISHING mode.

*Setting to User Select*

The User Select screen has an empty line at the bottom.

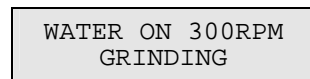
- Press W or **W** to select WATER ON or WATER OFF.
- Press SPEED d O or d N to select speed.

Press DISC A N or A O to change between the settings.

*Setting to GRINDING*

The standard setting GRINDING is preset to WATER ON.

- Press SPEED d O or d N to select speed.

Press DISC A N or A O to change between the settings.

*Setting to POLISHING*

The standard setting POLISHING is preset to WATER OFF.

- Press SPEED d O or d N to select speed.

*Settings with a  
RotoCom Memory Unit*

If RotoPol is connected to RotoCom and communication is in ROTOCOM mode, the following screen will appear:

WATER ON 300RPM  
PG-Paper #120

(Sample screen)

- The first line on the screen shows the settings of parameters for WATER and SPEED.

DISC A N and A O toggles between the different surfaces defined in RotoCom.

**STRUERS CONSUMABLES**

Struers offers a wide range of:

- Consumables for both plane and fine grinding.
- Polishing cloths, diamond and oxide abrasive materials.

*RotoPol-21/-22/-25*  
*Instruction Manual*

# Reference Guide

Table of Contents	Page
<b>1. Advanced Operations</b>	
Selection of Disc.....	22
Mounting of Specimen Mover (Optional).....	22
RotoForce-3 and -4.....	22
RotoForce-1.....	22
Mounting of Multidoser (optional).....	22
RotoCom Memory Unit (optional).....	22
Setting the Communication.....	23
Recirculation Cooling Unit (Optional).....	23
Configuration of RotoPol-22.....	24
Setting AUTO MODE when RotoForce-3 or -4 is mounted.....	24
Setting 2-DISC SYSTEM when RotoForce-1 is mounted....	24
Struers Metalog Guide™.....	25
Oxide Polishing.....	26
Overview of Menu Screens in the Setup Menu.....	27
RotoPol-21 and 25 Setup Menus.....	27
RotoPol-22 Setup Menus.....	28
<b>2. Accessories and Connected Equipment</b> .....	29
<b>3. Trouble-shooting</b> .....	30
<b>4. Maintenance</b>	
Daily Service.....	32
Weekly Service.....	32
Replacing the Cooling Water.....	32
Checking the Recirculation Cooling Unit.....	32
<b>5. Technical Data</b> .....	33

## **1. Advanced Operations**

### **Selection of Disc**

RotoPol is designed for wet grinding and polishing discs of up to 250 mm dia. However, the whole range of Struers discs can be mounted.

For automatic preparation of specimen holders mounted in specimen movers, only aluminium discs should be used. See the instructions for RotoForce-3 or RotoForce-4.

### **Mounting of Specimen Mover (Optional)**

*RotoForce-3 and -4*

RotoPol-21/22/25 can be equipped with a RotoForce-3 or RotoForce-4 specimen mover. To mount the specimen mover, see the appropriate Instruction Manual.

*RotoForce-1*

RotoPol-21 and -25 can be equipped with one RotoForce-1 specimen mover. RotoPol-22, having two independently operating discs, can be equipped with one or two RotoForce-1 specimen movers. To mount the specimen mover, see the appropriate Instruction Manual.

***IMPORTANT***

If RotoPol-21/22/25 has to be relocated,  
the specimen mover must be removed.

### **Mounting of Multidoser (optional)**

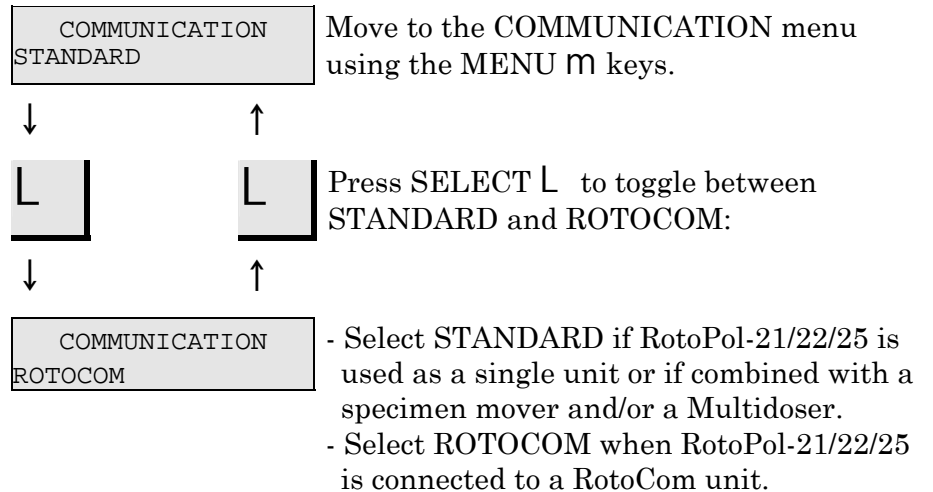
RotoPol-21/22/25, combined with a specimen mover, can be equipped with *Multidoser* for automatic dosing of suspensions and lubricants. To connect Multidoser, see the appropriate Instruction Manual.

### **RotoCom Memory Unit (optional)**

RotoPol-21/22/25, combined with RotoForce-3 or RotoForce-4 and Multidoser, can be equipped with a RotoCom memory unit. To connect RotoCom to RotoPol, see the appropriate Instruction Manual.

## Setting the Communication


Press MENU  $\mathfrak{M}$  to activate the setup menu.

- Press STOP  $\mathfrak{O}$  when the desired setting

## Recirculation Cooling Unit (Optional)
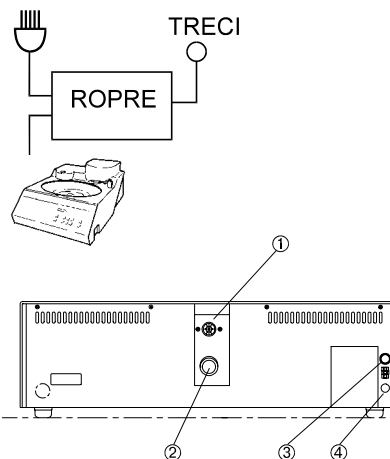
### ④ Socket for Recirculation

A Recirculation Cooling Unit (TRECI) can be connected by means of a Recirculation Cooling Unit Connector (ROPRE).

- Connect the cable from the pump in TRECI to ROPRE, according to the diagram of ROPRE.
- Connect the cable with plug from ROPRE to the socket marked Recirculation on the back of the RotoPol unit.
- Connect the cable from ROPRE to the mains.
- Check that the pump rotates in the direction indicated on the pump. If the direction is incorrect, switch two of the phases in the connection to the pump in TRECI (3-phase versions only).
- Replace the outlet tube on the pump in TRECI with the non-return valve with elbow and hexagon nipple from ROPRE. Struers recommends that you seal the threads.
- Connect the pressure hose of RotoPol to the hexagon nipple. Remember to mount the gasket.
- Select the reduction piece from ROPRE which fits the drain hose of RotoPol. Mount the piece on the outlet angle of the TRECI.
- Shorten the drain hose if necessary. Fix the drain hose on the reduction piece: strip away a short length of the steel reinforcement in the drain hose and mount a hose clamp on the stripped end.
- Mount the outlet angle on TRECI.

### **REMEMBER**

Avoid sharp bends in the drain hose and make sure that the hose slopes throughout its length.

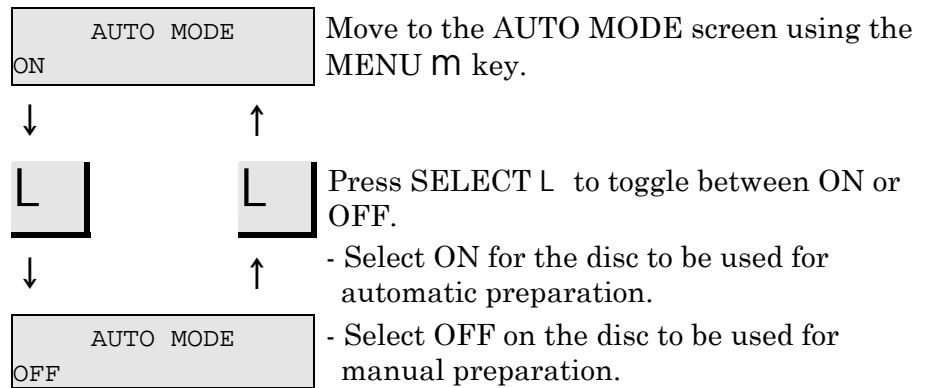
## Configuration of RotoPol-22

### *Setting AUTO MODE when RotoForce-3 or -4 is mounted*

As RotoPol-22 contains two individually controlled discs, there are some additional options.

One disc can be used for automatic preparation and the second disc can be used for manual preparation.

- Press MENU  to activate the setup menu.

- Press STOP  when the desired setting appears.


If both discs are to be used with the RotoForce sample movers, they should both be set to AUTO MODE ON.


### *Setting 2-DISC SYSTEM when RotoForce-1 is mounted*

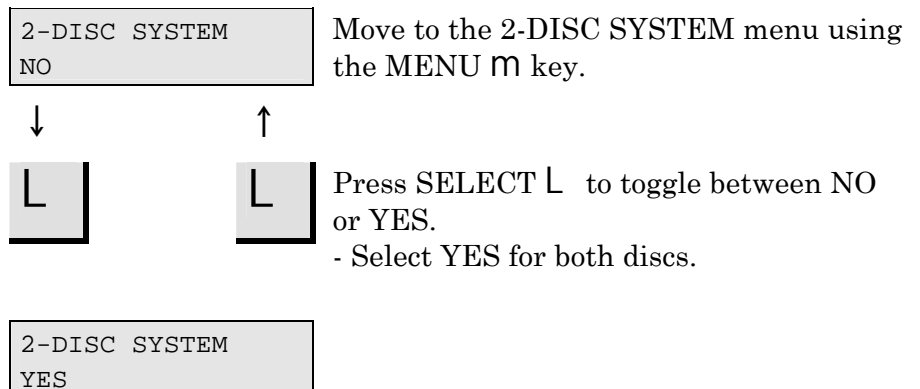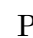1 or 2 RotoForce-1 sample movers can be mounted on RotoPol-22.

To start only RotoForce-1 with the respective disc, set 2-DISC SYSTEM to YES.

- Press MENU  to activate the setup menu.

- Press STOP  when the desired setting appears.

## **Struers Metalog Guide™**

RotoPol-21/22/25 is designed for both grinding and polishing. In Struers Metalog Guide™ you will find a detailed description of grinding/polishing for automated mechanical specimen preparation.

## **Metalog Guide**

Struers Metalog Guide™ offers preparation methods for the most common materials, based on a simple analysis of two key properties: hardness and ductility. Finding the right method, including choice of consumables, is easy. Always consult Struers Metalog Guide™ for the correct preparation method for the actual specimens.

Struers Metalog Guide™ contains 6 useful chapters:

- **Metalogram**: a quick and safe guide to the right preparation method.
- **Metalog Methods**: a complete catalogue of preparation methods, based on Struers' extensive materialographic experience, and employing Struers' range of consumables.
- **Preparation Philosophy**: the basics of modern specimen preparation, seen from a professional point of view.
- **Metalog Process**: the materialographic preparation process from start to finish, logically explained.
- **Metalog Master**: a combined trouble-shooting guide and supply of in-depth information on the processes of mechanical preparation, including an expert system for the solving of preparation problems.
- **Metalog Code**: quick access to the relevant consumables for the chosen preparation methods.

### **Struers Metalog Guide™**

A complete guide to materialographic specimen preparation. Contact your local dealer for a free copy of Metalog Guide™.

## **Oxide Polishing**

For polishing with Oxide Polishing Suspension: at the end of the polishing process, select WATER ON W. Adjust the water tap if necessary. Let the water flow until the oxide polishing suspension has been cleaned away (about 10 seconds).

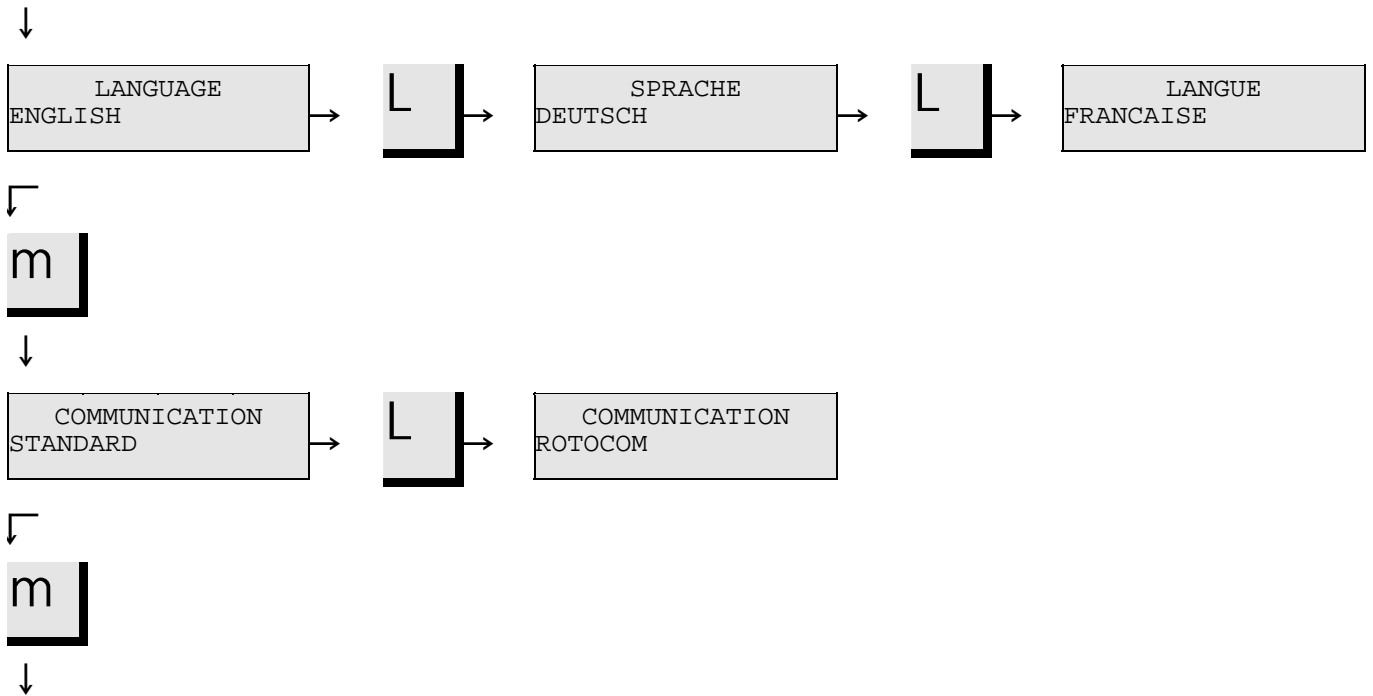
***IMPORTANT***

If a Recirculation Cooling Unit is connected, special precautions must be taken:

- the water coming from the built-in water tap cannot be used to wash the oxide polishing suspension away. Use clean water.
- The outflow must not be conducted to the Recirculation Cooling Unit. Direct the outlet to the drain or collect the outflow in a bucket.

**Overview of Menu Screens  
in the Setup Menu**


*RotoPol-21 and 25 Setup Menus*

The settings in the left-hand column of the screens show the default values.

*RotoPol-21/-22/-25  
Instruction Manual*

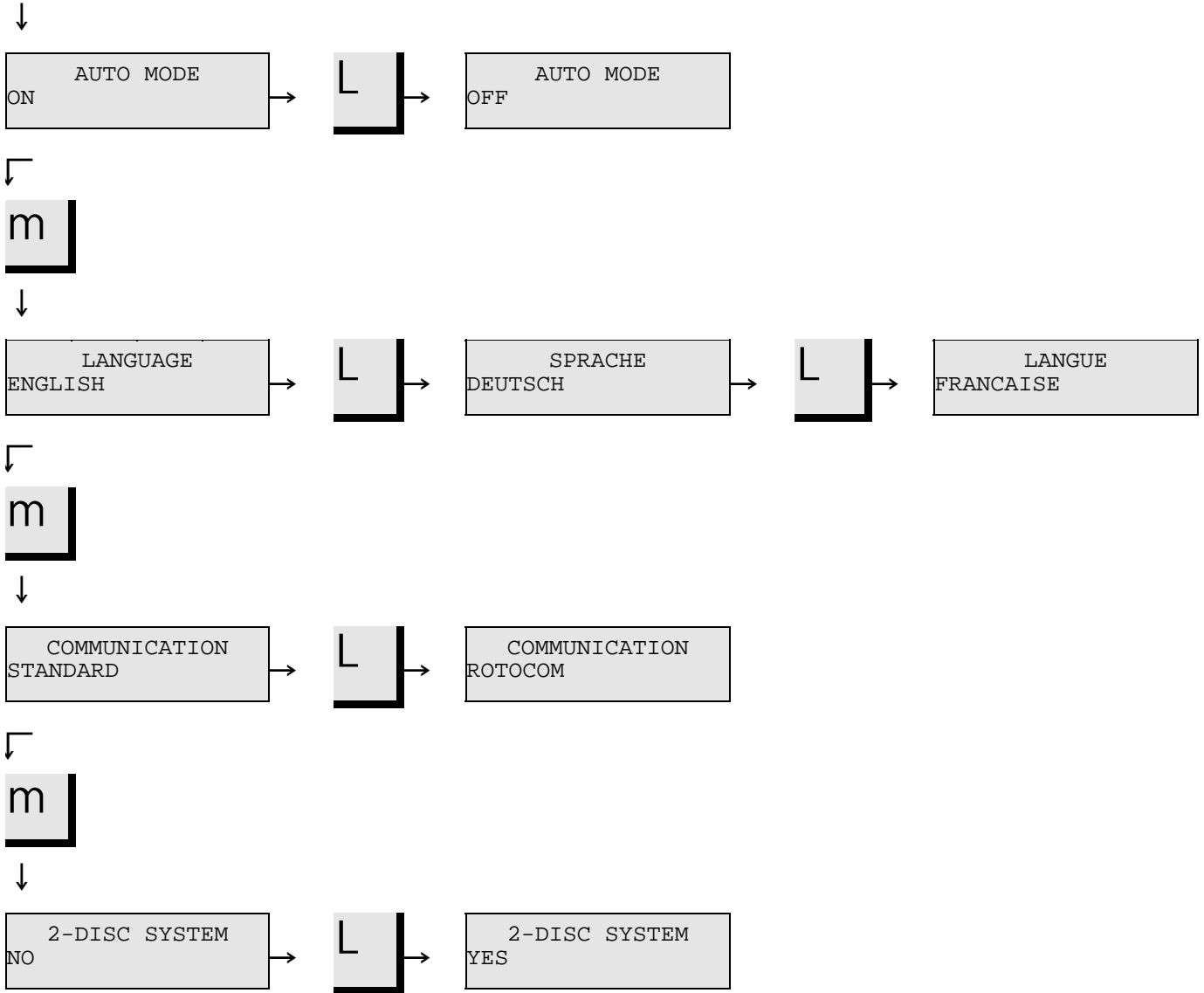*RotoPol-22 Setup Menus*

The settings in the left-hand column of the screens show the default values.

## **2. Accessories and Connected Equipment**

<b>Specification</b>	<b>Code</b>
<b>Accessories</b>	
<i>Lid</i>	
Lid for RotoPol-21/-22/-25	ROLID
<i>Wet Grinding Discs with Ring</i>	
250 mm dia., aluminium	ROTIF
<i>Polishing Discs</i>	
250 mm dia., disc for magnetic fixation	DEMIF
250 mm dia., aluminium	DEDIF
<b>Connected Equipment</b>	
<i>RotoForce-4, specimen mover</i>	ROFO4
<i>RotoForce-3, specimen mover</i>	ROFO3
<i>RotoForce-1, specimen mover</i>	ROFO1
<i>Multidoser (6 Bottles)</i> for automatic dosing of 6 liquids (1-5 suspensions, 0-4 lubricants and 1 OP-Suspension)	MULDO
<i>Multidoser Base Unit (3 Bottles)</i> for automatic dosing of 3 liquids (0-3 suspensions, 0-3 lubricants)	MULBA
<i>RotoCom, memory unit</i>	ROCOM
<i>Recirculation Cooling Unit</i>	TRECI
<i>Recirculation Cooling Unit Connector</i> for connection of Recirculation Cooling Unit (TRECI) to grinding/polishing machines	ROPRE

### 3. Trouble-shooting

Error	Explanation	Action
<b>Error Message</b>		
OVERLOAD	The motor has been overloaded for a period and the thermo switch has stopped the motor. The motor stops.	Wait until the motor has cooled down. When the display returns to the previous message, the machine can be re-started.
<b>Acoustic Signal</b>		
Long beep.	The selected function is protected. It cannot be activated.	Check the setting.
<b>Machine Problems</b>		
Noise when the machine starts or the machine will not pull.	The belt is not tight enough.	Call a Struers service technician. The belt must be tightened.
Disc rotates in the wrong direction (not for RotoPol-25).	Wrong connection to the power supply.	Switch two of the phases.
Motor buzzes (not for RotoPol-25).	One phase lacking.	Check all fuses of local power supply.
Functions are not carried out. RotoCom and/or Multidoser do not operate.	Fuse at the rear of RotoPol blown.	Replace the fuse.
For RotoPol-25 only: The machine does not operate.	One of the fuses on the right-hand side of the cabinet blown.	Replace the fuse.
Water not draining away.	Drain hose squeezed.	Straighten the hose.
	Drain hose clogged.	Clean the hose.
	Drain hose does not slope downward.	Adjust the hose to an even slope.
Water dripping underneath the machine.	The water does not drain away.	See above.
Cooling water stops.	Wrong software setting.	Check software setting.
	Water tap on mains closed.	Open for water.
	Built-in water tap closed.	Open for water.
	Built-in water tap blocked.	Clean water tap.
	Filter at the water inlet blocked.	Clean filter.
Cooling water drips after stop.	Defect in the solenoid valve.	Call a Struers service technician. The solenoid valve must be replaced.
The disc vibrates.	Dirt on the underside of the disc or on the turntable.	Clean the contact face between the disc and turntable.

*RotoPol-21/-22/-25*  
*Instruction Manual*

JAPANESE TEXTS	All RotoPol modules have the possibility of showing the screen texts in Japanese. When choosing a European language in the LANGUAGE menu, you might accidentally press M MENU at the Japanese option.	To return to English, German or French texts, just follow this procedure:  <i>RotoPol -21 and -25:</i> Press O STOP. Press M MENU. Press L SELECT until you reach the language you prefer. Press O STOP.  <i>RotoPol -22:</i> Press O STOP. Press M MENU two times. Press L SELECT until you reach the language you prefer. Press O STOP.
----------------	---	---

## 4. Maintenance

### Daily Service

Clean all accessible surfaces with a moist cloth.

### Weekly Service

Remove the grinding/polishing disc and remove all dirt from the tray and the drain tube.

### Replacing the Cooling Water

*Only relevant when a Recirculation Cooling Unit is connected:*  
Replace the cooling water in the Recirculation Cooling Unit at least once a month.

- Place a Disposable Plastic Container (TREPO) in the tank and fold it over the edge.
- Fill the tank with 30 l of water. NB! Too high a level of coolant in the tank could damage the pump.
- Add Struers Additive according to the instructions.
- Put the lid on the tank.
- Place the pump in the lid

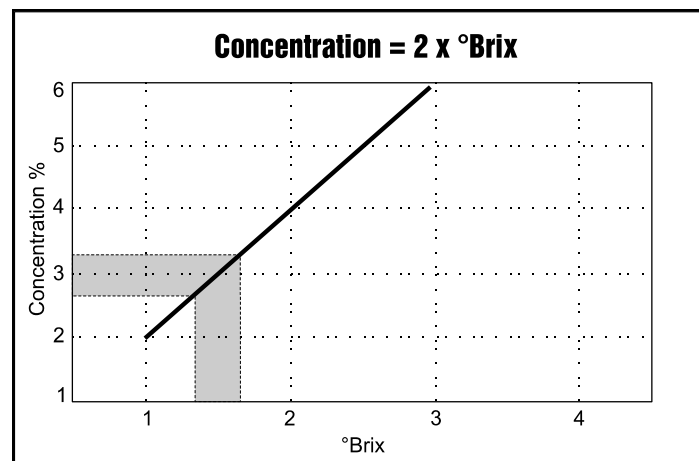
#### IMPORTANT

Always maintain the correct concentration of Struers Additive in the cooling water (percentage stated on the container of the Additive). Remember to add Struers Additive each time you refill with water.

### Checking the Recirculation Cooling Unit

The cooling unit should be checked for cooling water every week and refilled when the flushing pump cannot reach the cooling water. Remember to add Struers Additive for Cooling Fluid: One part of Additive for 33 parts of water.

The concentration of additive should always lie between 2.7 and 3.3%; the pH should be around 9. Replace the cooling water if pH is  $\leq 8.0$ . To check the concentration of additive, use a refractometer (0-10 Brix, see diagram.). Concentration =  $2 \times$  Brix. Add Struers Additive for Cooling Fluid if the concentration is lower than 2.7%. Add water if the concentration is higher than 3.3%.

## 5. Technical Data

Subject		Specifications		
		RotoPol-21	RotoPol-22	RotoPol-25
Disc	Diameter	200-250mm / 8-10"	200-250mm / 8-10"	200-250mm / 8-10"
	Speed	150/300rpm	150/300rpm	40-600rpm in steps of 10
	Rotational direction	Anti-clockwise	Anti-clockwise	Anti-clockwise
	No. of motors	1	2	1
	Motor	200W / 0.27HP at 150 rpm 400W / 0.54HP at 300 rpm	2 x 200W / 0.27HP at 150 rpm 2 x 400W / 0.54HP at 300 rpm	550W / 0.75HP at 600 rpm
	Torque at disc	Continuous 12.7Nm / 9.4lb-ft Max. > 18.0Nm / 13.3lb-ft	Continuous 12.7Nm / 9.4lb-ft Max. > 18.0Nm / 13.3lb-ft	Continuous 7.3Nm / 5.4lb-ft Max. > 14.6Nm / 10.8lb-ft
Connected specimen movers	RotoForce-1	1	2	1
	RotoForce-3 and -4	1	1	1
Start/stop	RotoForce-1 and Water recirculation	Yes	Yes	Yes
Automatic stop	RotoForce-3/4 and Multidoser	Yes	Yes	Yes
Data exchange	RotoCom	Yes	Yes	Yes
Software and electronics	LC display	2x16 characters	2x16 characters	2x16 characters
	Controls	Touch pads	Touch pads	Touch pads
	Memory	EPROM/RAM/NV-RAM	EPROM/RAM/NV-RAM	EPROM/RAM/NV-RAM
Fittings	No. of water taps	2	2	2
	Splash rings	2	2	2
Safety standards	General	89/392/EEC & 91/386/EEC & 93/44/EEC	EN292-1, EN292-2, EN60204-1 (IEC204-1)	
	EMC	89/336/EEC & 92/231/EEC	EN50081-1, EN50082-1	FCC part 15, subpart B
Noise level	At idle running, at a distance 1.0 m / 39.4" from the machine	52 dB (A)	54 dB (A)	58 dB (A)
Surrounding temperature		5-40°C / 41-104°F	5-40°C / 41-104°F	5-40°C / 41-104°F
Humidity	Non condensing	0-95%RH	0-95%RH	0-95%RH

*RotoPol-21/-22/-25*  
*Instruction Manual*

Subject		Specifications		
		RotoPol-21	RotoPol-22	RotoPol-25
Supply	Power	3 (3L+PE) phases	3 (3L+PE) phases	1 (N+L+ PE) phase
	Power consumption	660W / 0.90HP	1260W / 1.7HP	600W / 0.8HP
	Voltage / frequency	Max continuous load		
	3 x 200V / 50Hz	2.2A	4.0A	-
	3 x 200-210V / 60Hz	2.0A	3.6A	-
	3 x 220-230V / 50Hz	2.3A	4.2A	-
	3 x 220-240V / 60Hz	2.1A	3.8A	-
	3 x 380-415V / 50Hz	1.3A	2.4A	-
	3 x 460-480A / 60Hz	1.1A	2.0A	-
	1 x 100V / 50-60Hz	-	-	8.0A
	1 x 110-120V / 50-60Hz	-	-	8.0A
	1 x 220-240V / 50-60Hz	-	-	4.0A
	Pressure for tap water	1-10 bar / 14.5-145 psi	1-10 bar / 14.5-145 psi	1-10 bar / 14.5-145 psi
	Water inlet	½" or ¾"	½" or ¾"	½" or ¾"
Water outlet	ø32 mm / 1 1/4"	ø32 mm / 1 1/4"	ø32 mm / 1 1/4"	
Dimensions and weight	Width	700mm / 27.6"	700mm / 27.6"	700mm / 27.6"
	Depth	700mm / 27.6"	700mm / 27.6"	700mm / 27.6"
	Height	215mm / 8.5"	215mm / 8.5"	215mm / 8.5"
	Weight	37.5kg / 82.7lbs	47kg / 103.6lbs	39kg / 86.0lbs

# Quick Reference Guide for RotoPol-21 and -22

## Mounting a Disc

- Place the disc on the turntable and turn it until the three driving pins engage with the holes in the turntable.

## Starting the Disc

- Press **START S** . The disc starts rotating. Selected functions are executed and connected RotoSystem modules start operating.

## Stopping the Disc

- Press **STOP O** . The disc stops rotating. Selected functions and connected RotoSystem modules stop.

## Grinding/Polishing

- Mount the required surface on a disc.
- Mount the disc on the turntable.
- Select the desired functions, if RotoCom is not connected.
- Press **START S** .

## Settings

### Setting to User Select

WATER OFF 150RPM

The User Select screen has an empty line at the bottom.

- Press **W** or **W** to select WATER ON or WATER OFF.

- Press **SPEED d 150** or **d 300** to select speed.

Press **DISC A N** or **A O** to change between the settings.

↓                      ↑  
r                      A O

### Setting to GRINDING

WATER ON 300RPM  
GRINDING

The standard setting GRINDING is preset to WATER ON and 300RPM.

Press **DISC A N** or **A O** to change between the settings.

↓                      ↑  
A N                      A O

### Setting to POLISHING

WATER OFF 150RPM  
POLISHING

The standard setting POLISHING is preset to WATER OFF and 150RPM.

# Quick Reference Guide for RotoPol-25

## Mounting a Disc

- Place the disc on the turntable and turn it until the three driving pins engage with the holes in the turntable.

## Starting the Disc

- Press **START S** . The disc starts rotating. Selected functions are executed and connected RotoSystem modules start operating.

## Stopping the Disc

- Press **STOP O** . The disc stops rotating. Selected functions and connected RotoSystem modules stop.

## Grinding/Polishing

- Mount the required surface on a disc.
- Mount the disc on the turntable.
- Select the desired functions, if RotoCom is not connected.
- Press **START S** .

## Settings

### Setting to User Select

WATER OFF 150RPM

The User Select screen has an empty line at the bottom.

- Press **W** or **W** to select WATER ON or WATER OFF.

- Press **SPEED d 150** or **d 300** to select speed.

Press **DISC A N** or **A O** to change between the settings.

↓                      ↑  
**A N**                      **A O**

### Setting to GRINDING

WATER ON 300RPM  
 GRINDING

The standard setting GRINDING is preset to WATER ON

- Press **SPEED d O** or **d N** to select speed.

Press **DISC A N** or **A O** to change between the settings.

↓                      ↑  
**A N**                      **A O**

### Setting to POLISHING

WATER OFF 150RPM  
 POLISHING

The standard setting POLISHING is preset to WATER OFF.

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*RotoPol-21/-22/-25*  
*Spare Parts and Diagrams*

---

**Always state *Serial No* and *Voltage/frequency*  
if you have technical questions or when ordering spare parts.**

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Struers assumes no responsibility for errors in the manual text/illustrations. The information in this manual is subject to changes without notice. The manual may mention accessories or parts not included in the present version of the equipment.

The contents of this manual is the property of Struers. Reproduction of any part of this manual without the written permission of Struers is not allowed.

All rights reserved. © Struers 2007.

**Struers A/S**  
Pederstrupvej 84  
DK-2750 Ballerup  
Denmark  
Telephone +45 44 600 800  
Fax +45 44 600 801

---

# Spare Parts and Diagrams

<b>Table of contents</b>	<b>Drawing</b>
<b>RotoPol-21</b>	
<i>Drawings</i>	
Bottom, complete, RotoPol-21 .....	15150021E
Water Tap, Complete .....	15200050B
<i>Diagrams</i>	
Wiring Diagram .....	15153100A
<b>RotoPol-22</b>	
<i>Drawings</i>	
Bottom, complete, RotoPol-22 .....	15160022E
Water Tap, Complete .....	15200050B
<i>Diagrams</i>	
Wiring Diagram (2 pages) .....	15163100B
<b>RotoPol-25</b>	
<i>Drawings</i>	
Bottom, complete, RotoPol-25 .....	15170023G
Electrical Drawer, Complete, RotoPol-25.....	15170043B
Water Tap. Complete .....	15200050B
<i>Diagrams</i>	
Wiring Diagram .....	15173100C

Some of the drawings may contain position numbers  
not used in connection with this manual.

*RotoPol-21/-22/-25*  
*Spare Parts and Diagrams*

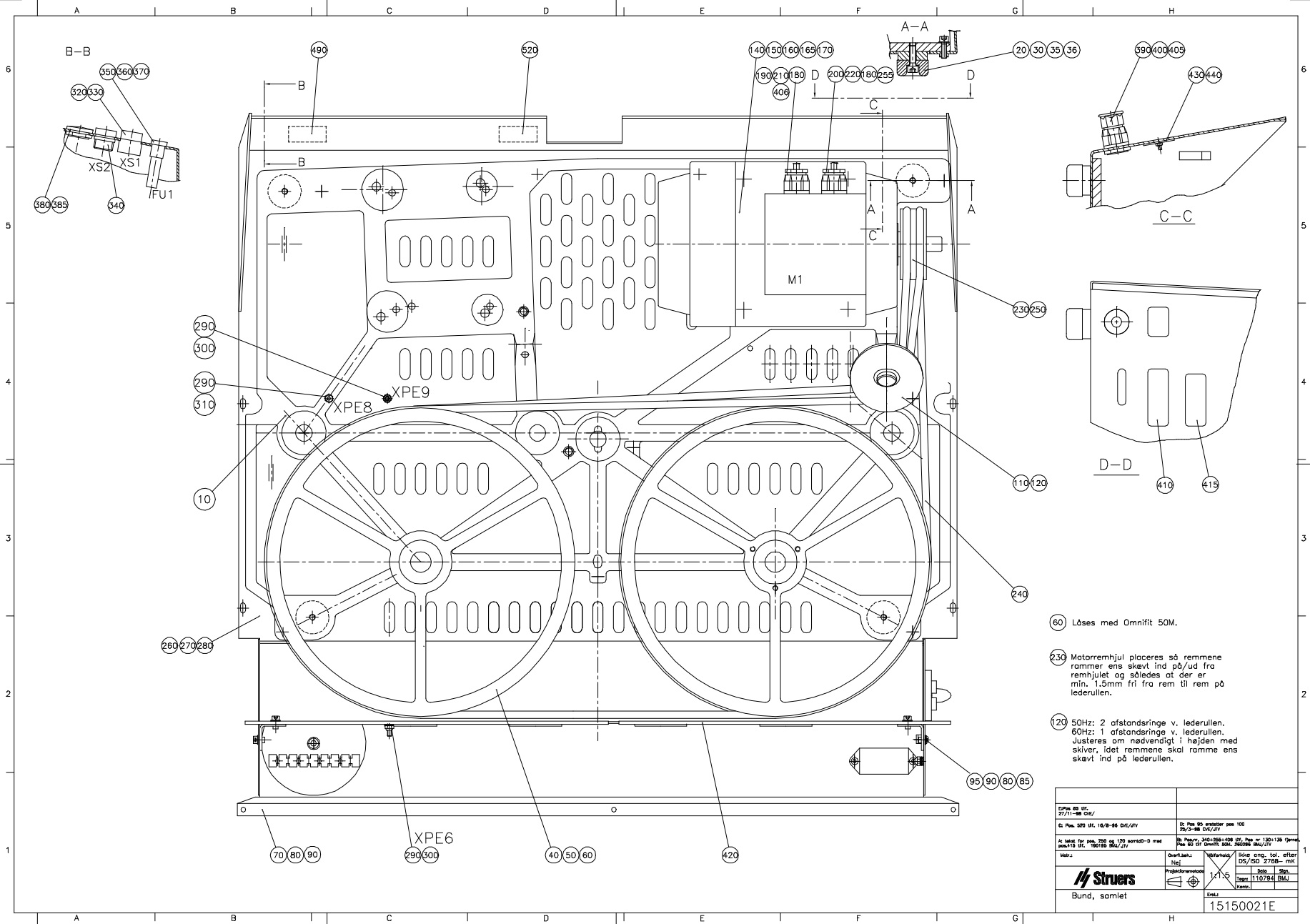
<b>Drawing</b>	<b>Pos.</b>	<b>Spare Part</b>	<b>Part No.</b>
Various Parts	070	<b>RotoPol-21</b>	
		Splash ring	14910120
		Pressure hose 2x3/4"	2NU93020
15150021		<b>Bottom, complete, RotoPol-21</b>	
	240	Driving belt, 50Hz (2 pcs.)	14950126
	240	Driving belt, 60Hz (2 pcs.)	14950125
	360	Fuse holder (bayonet)	2XS19816
	370	2.50A T fuse glass tube (10 pcs.)	RFU14000
15200050		<b>Water Tap, Complete</b>	15200050
		Gaskets for water tap	R4910071
	090	Water tube and tap, complete	15202901

*RotoPol-21/-22/-25*  
*Spare Parts and Diagrams*

<b>Drawing</b>	<b>Pos.</b>	<b>Spare Part</b>	<b>Part No.</b>
Various Parts	070	<b>RotoPol-22</b>	
		Splash ring	14910120
		Pressure hose 2x3/4"	2NU93020
<b>15150022</b>		<b>Bottom, complete, RotoPol-22</b>	
	270	Driving belt, 50Hz (2 pcs.)	14960085
	270	Driving belt, 60Hz (2 pcs.)	14960086
	400	Fuse holder (bayonet)	2XS19816
	410	2.50A T fuse glass tube (10 pcs.)	RFU14000
<b>15200050</b>		<b>Water Tap, Complete</b>	15200050
		Gaskets for water tap	R4910071
	090	Water tube and tap, complete	15202901

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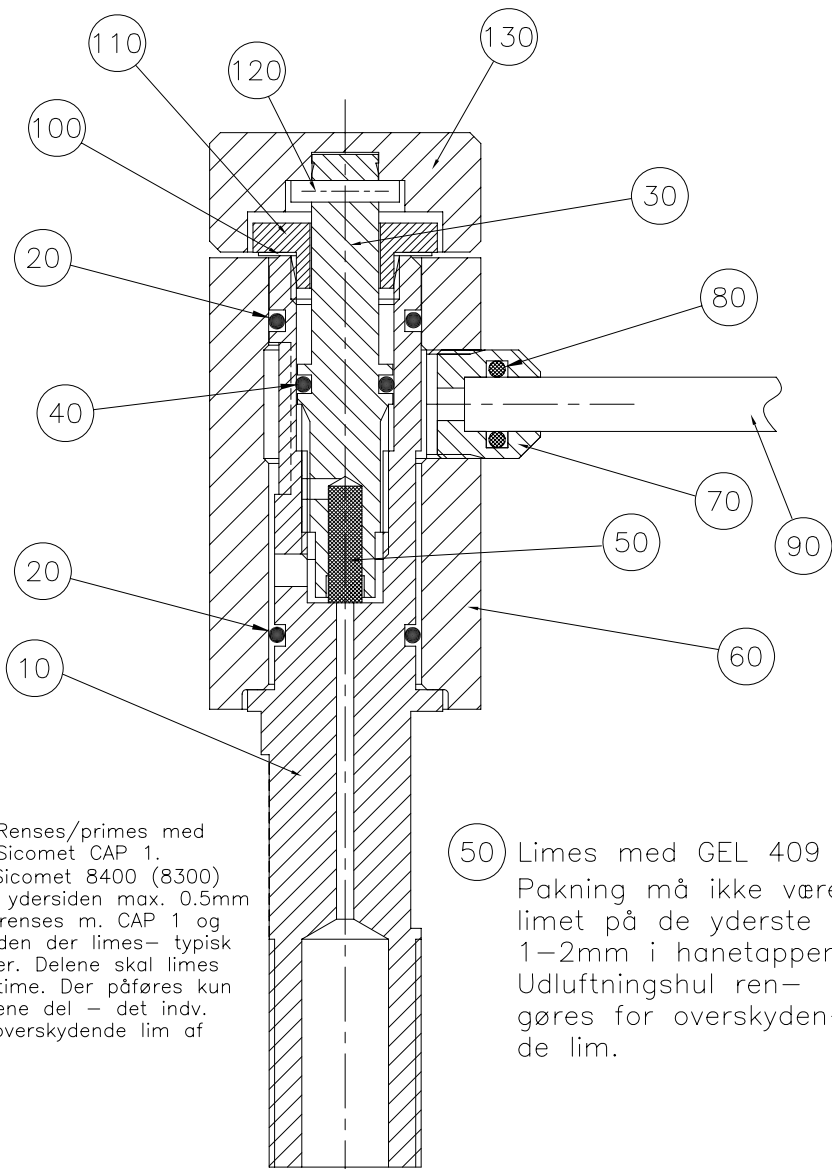
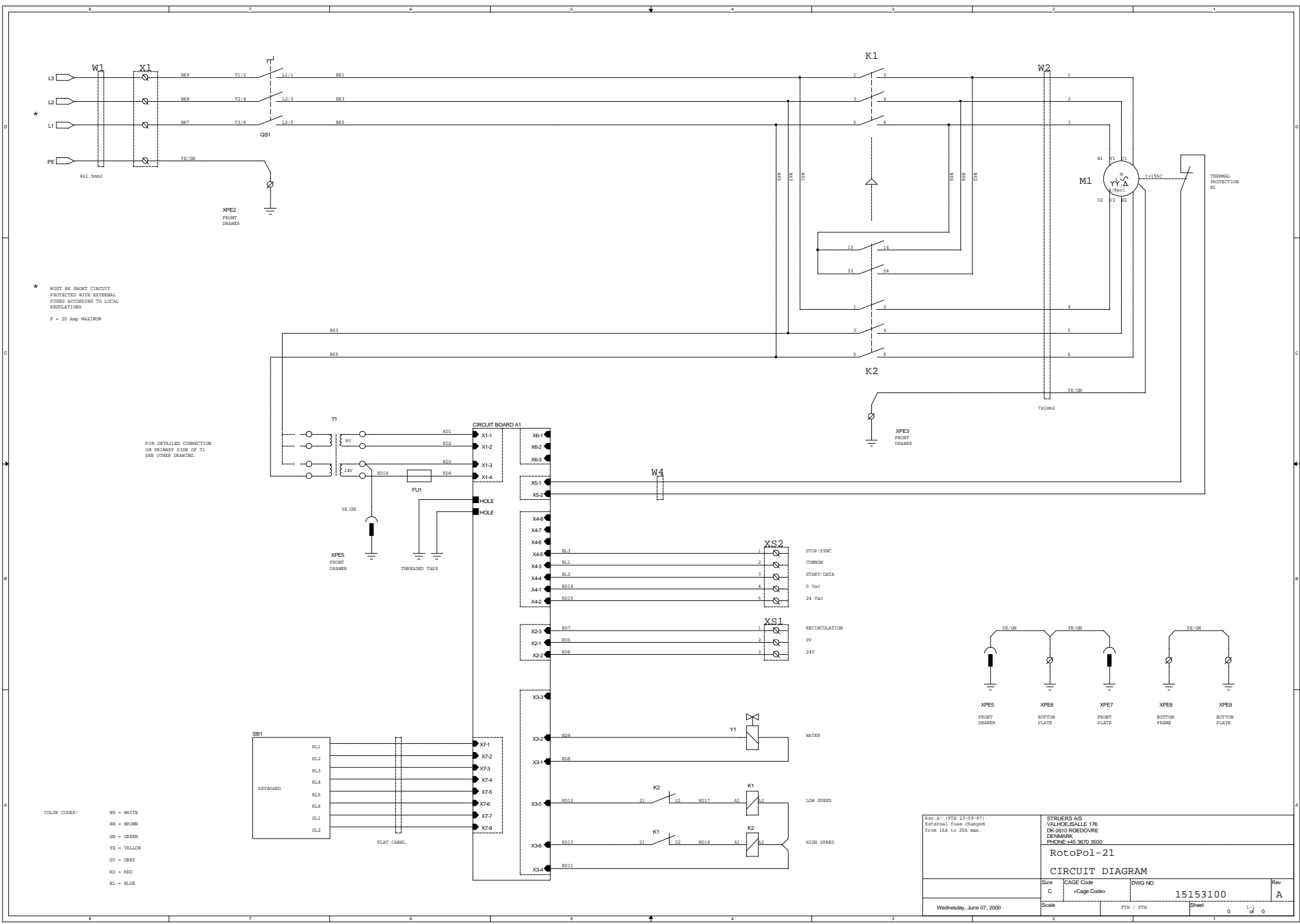
\* MUST BE SHORT CIRCUIT PROTECTED WITH EXTERNAL FUSES ACCORDING TO LOCAL REGULATIONS  
F = 20 Amp MAXIMUM


FOR DETAILED CONNECTION ON PRIMARY SIDE OF T1 SEE OTHER DRAWING.

COLOR CODES:  
WH = WHITE  
BN = BROWN  
GN = GREEN  
YE = YELLOW  
GY = GREY  
RD = RED  
BL = BLUE

Rev. A1 (FTH 23-99-97) External fuse changed from 16A to 20A max.		STRAUERS AS VALHOEJUSALLE 176 DK-2610 ROEDOVRE DENMARK PHONE: +45 3670 3500	
<b>RotoPol-21</b>			
<b>CIRCUIT DIAGRAM</b>			
Size C	CAGE Code <Cage Code>	DWG NO 15153100	Rev A
Wednesday, June 07, 2000		Scale FTH / FTH	Sheet 0 13 of 0

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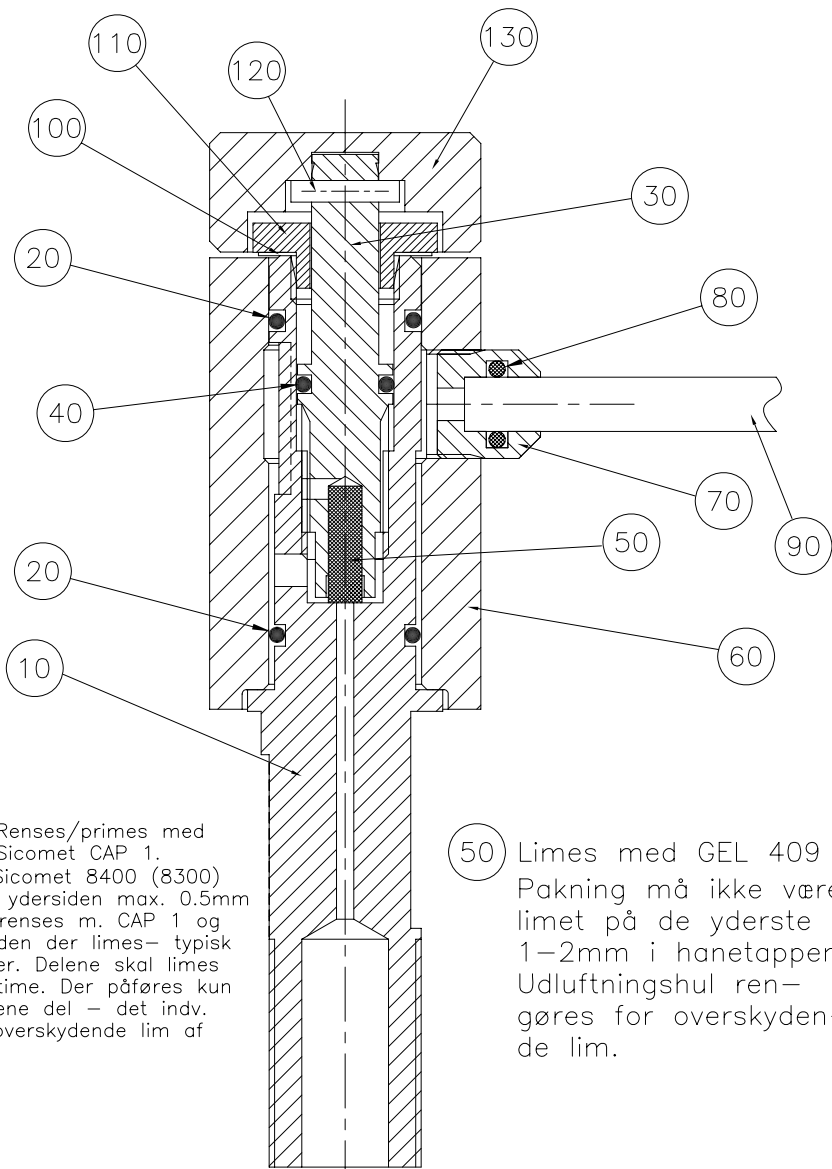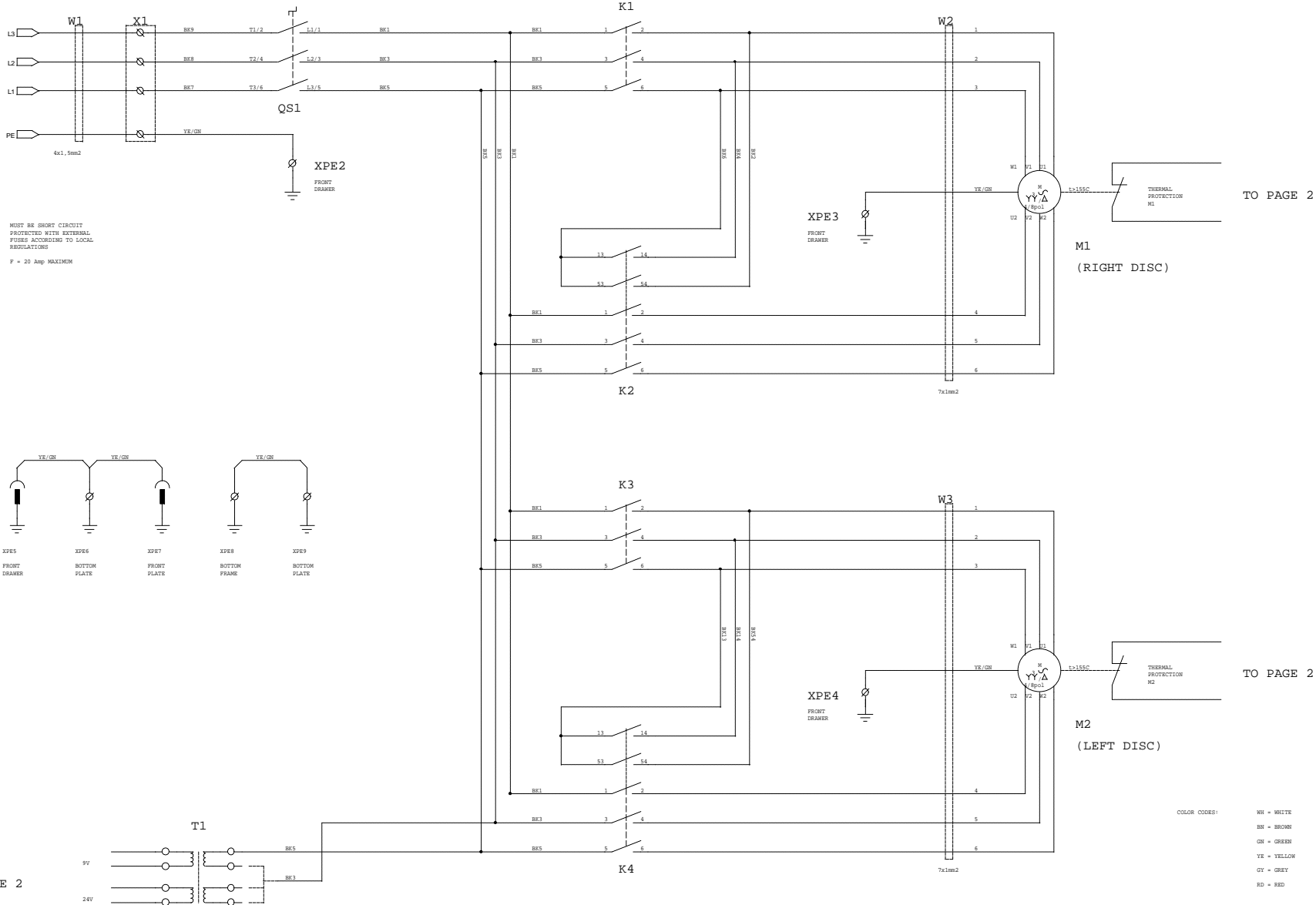
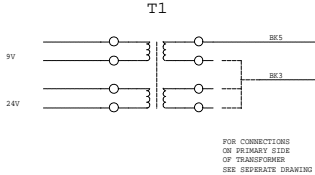
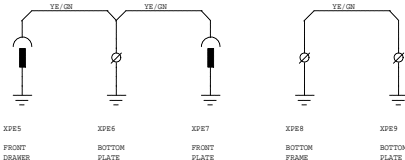
\* MUST BE SHORT CIRCUIT PROTECTED WITH EXTERNAL FUSES ACCORDING TO LOCAL REGULATIONS  
 F = 20 Amp MAXIMUM

FOR CONNECTIONS ON PRIMARY SIDE OF TRANSFORMER SEE SEPARATE DRAWING


COLOR CODES:  
 WH = WHITE  
 BR = BROWN  
 GN = GREEN  
 YE = YELLOW  
 GR = GREY  
 RD = RED

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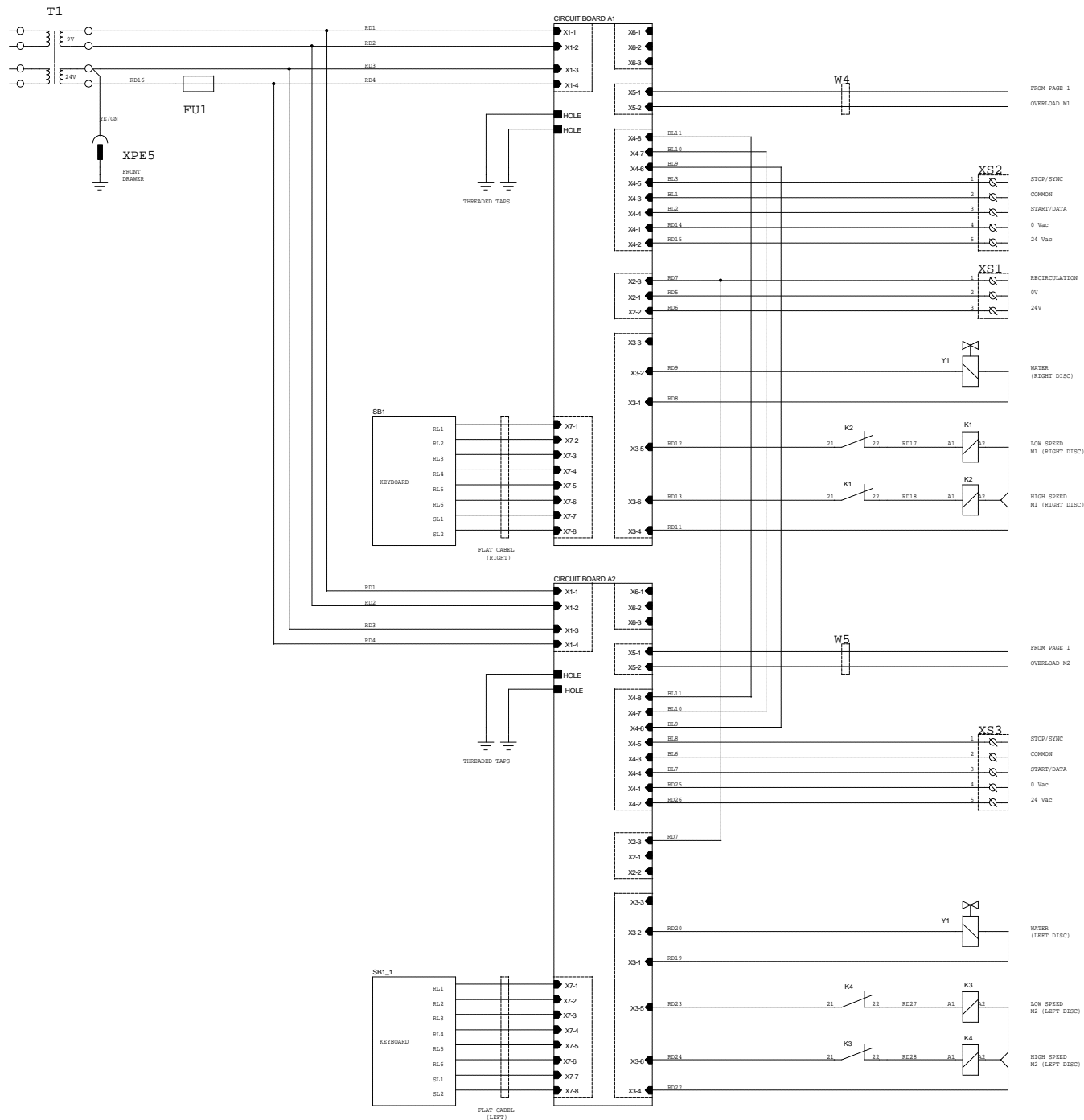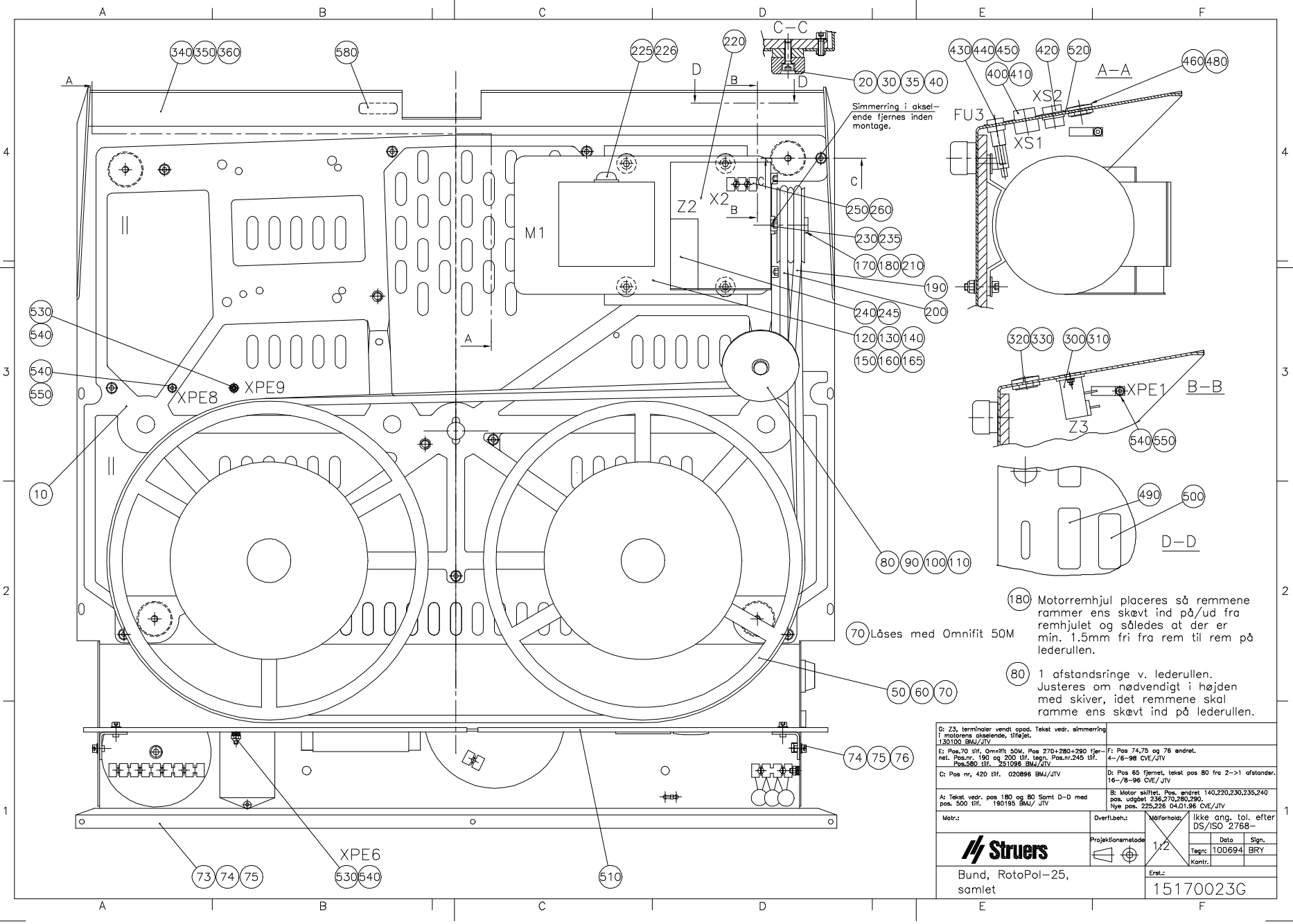
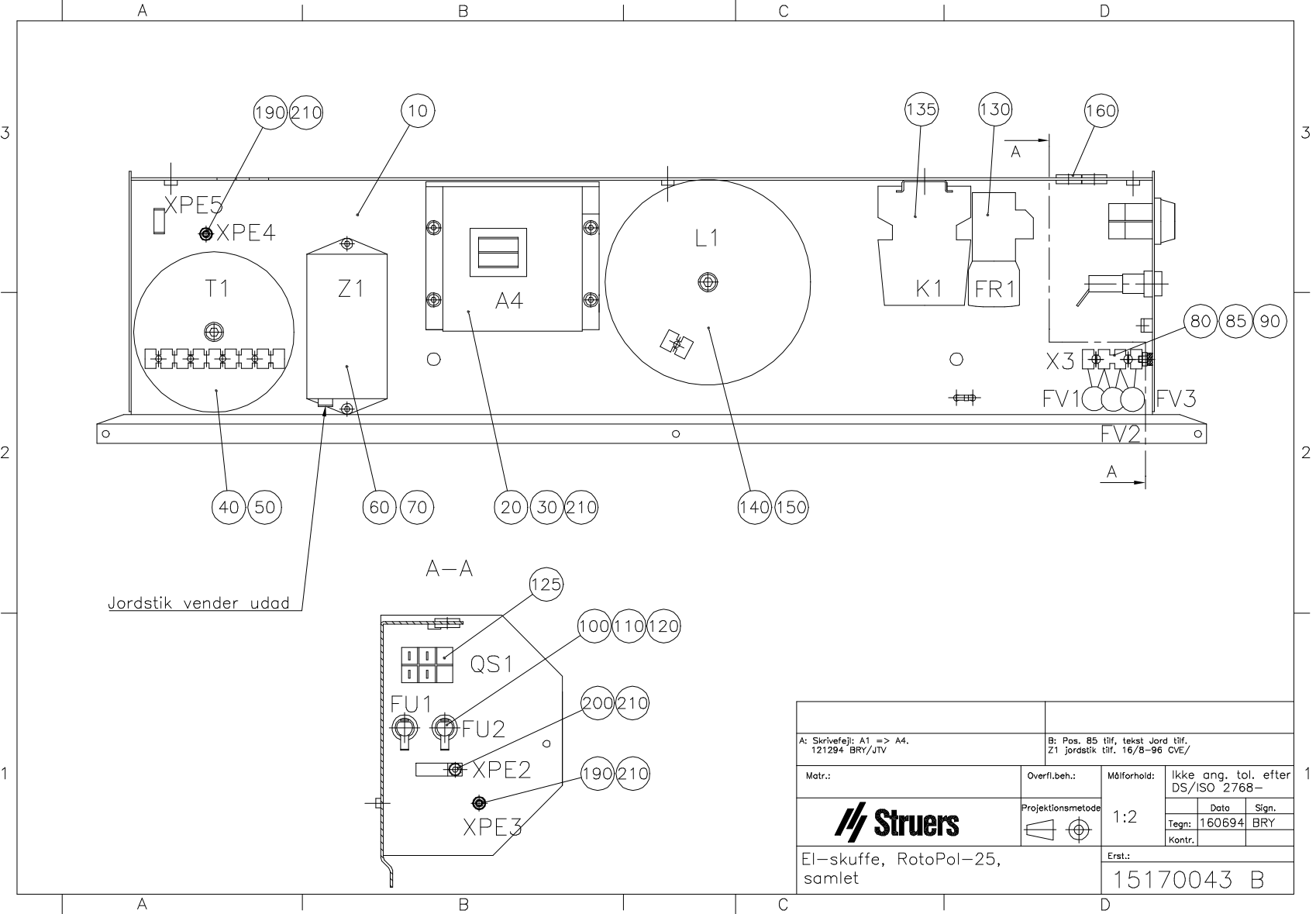
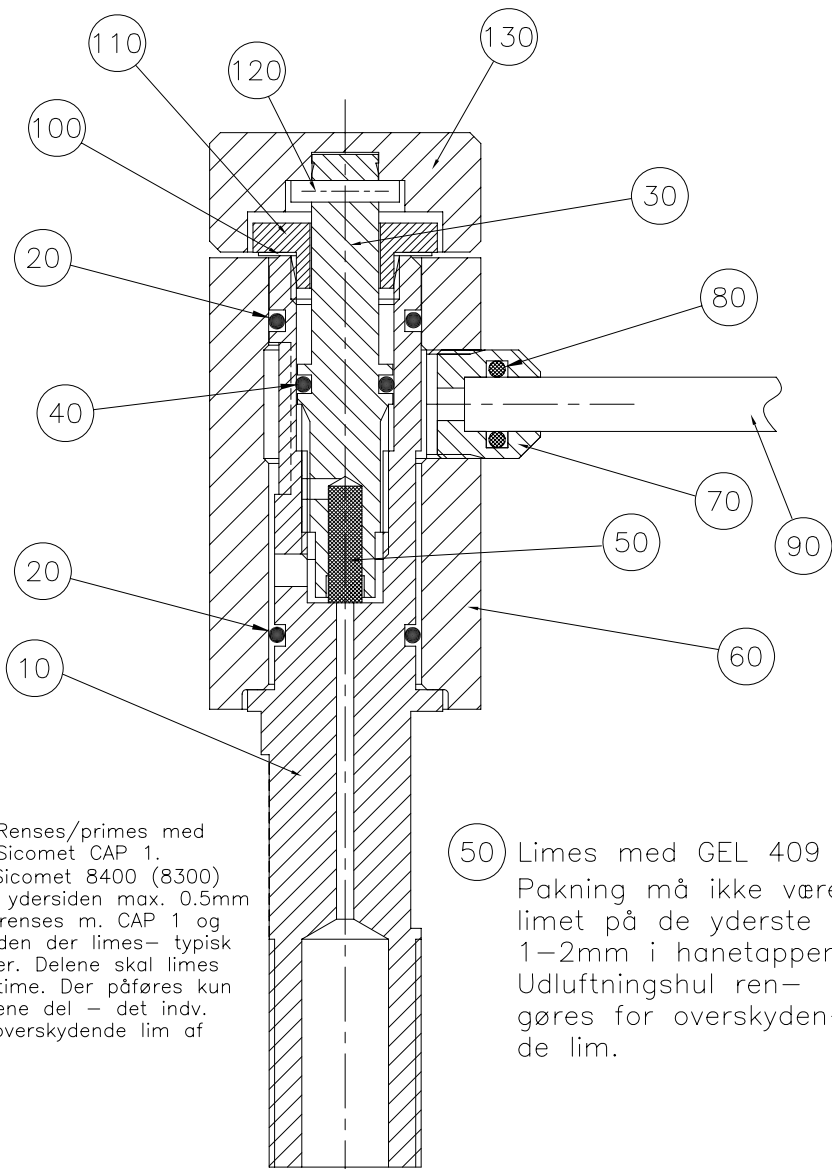
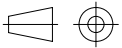
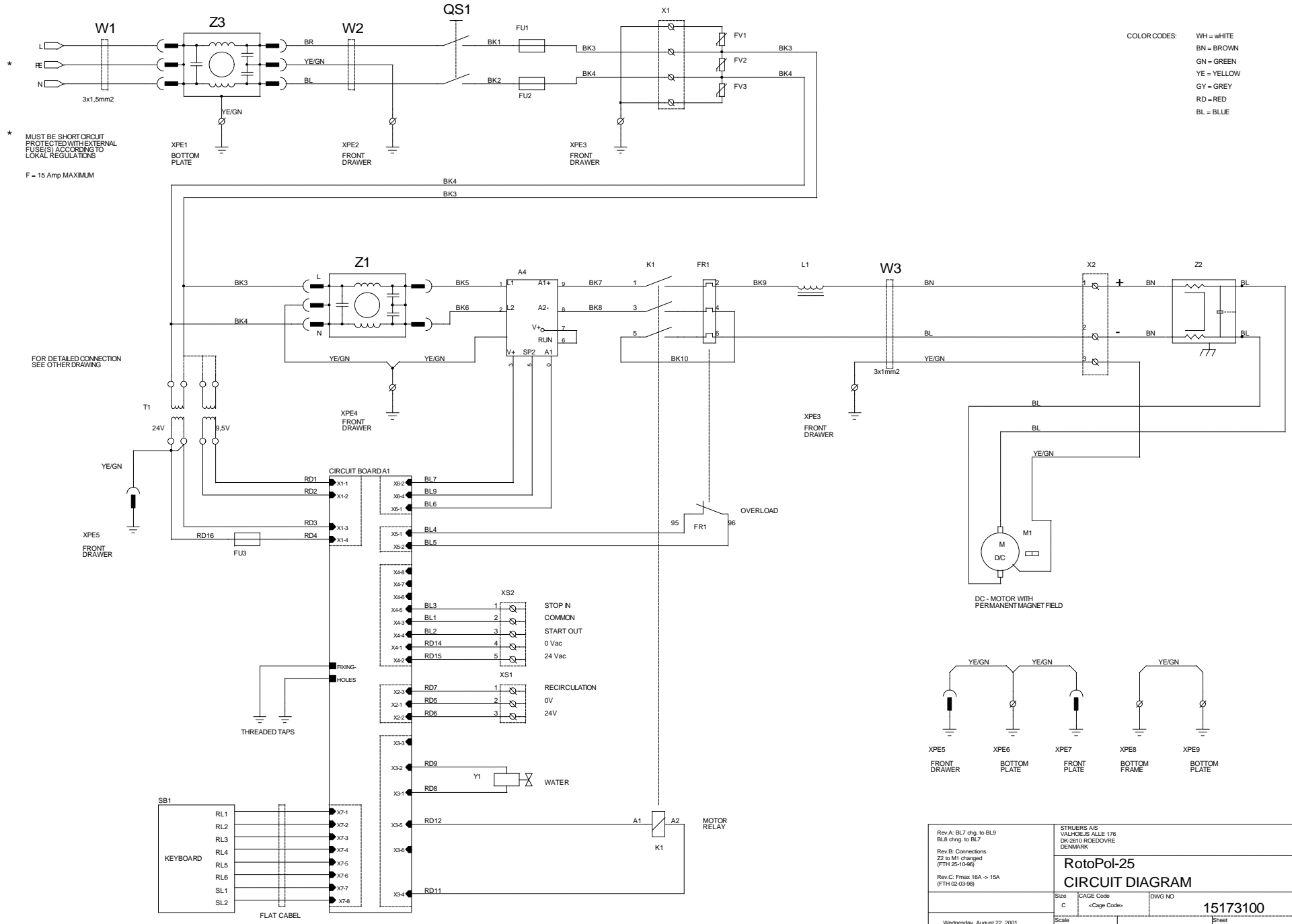
\* MUST BE SHORTCIRCUIT PROTECTED WITH EXTERNAL FUSE(S) ACCORDING TO LOCAL REGULATIONS  
 F = 15 Amp MAXIMUM

FOR DETAILED CONNECTION SEE OTHER DRAWINGS

Rev.A: BL7 chg. to BL9 BL8 chng. to BL7 Rev.B: Connections Z2 to K1 changed (FTH 25-10-96) Rev.C: Fmax 16A -> 15A (FTH 02-03-96)		STRUERS AS VALHØJES ALLE 176 DK-2610 ROEDOVRE DENMARK	
<b>RotoPol-25</b> <b>CIRCUIT DIAGRAM</b>			
Size C	CAGE Code <Cage Code>	DWG NO 15173100	Rev C
Wednesday, August 22, 2001		Scale	Sheet 1 of 1

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Herewith declares that

<i>Product Name:</i>	RotoPol-31, -35
<i>Type No:</i>	492, 528
<i>Machine Type:</i>	Grinding and polishing machine

is in conformity with the provisions of the following directives:

**Safety of Machinery** 89/392/EEC and 91/368/EEC and 93/44/EEC according to the following standard(s):  
EN292-1:1991, EN292-2:1991, EN60204-1:1992.

**EMC-Directive** 89/336/EEC and 92/31/EEC according to the following standard(s):  
EN50081-1:1991, EN50082-1:1992.

**Low Voltage Directive** 73/23/EEC and 93/68/EEC according to the following standard(s):  
EN60204-1:1992.

**Supplementary Information** The equipment complies with the American standards:  
FCC part 15, subpart B and UL508.

The above has been declared according to the global method, module A

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Klaus Kisbøll, Entwicklungsleiter, Struers A/S
**Fabricant** Struers A/S
Klaus Kisbøll, Chef du service développement, Struers A/S