LaboPol-21

LaboPol-25

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LaboPol-21/ -25

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English

Declaration of Conformity

**Manufacturer,
responsible for
Technical File**

Struers A/S
Pederstrupvej 84
DK-2750 Ballerup, Denmark
Telephone +45 44 600 800

Herewith declares that

<i>Product Name:</i>	LaboPol-21
<i>Type No:</i>	529
<i>Machine Type:</i>	Grinding and polishing machine

is in conformity with the provisions of the following directives:

Safety of Machinery 2006/42/EC according to the following standard(s):
EN ISO 12100:2010, EN 60204-1:2006/AC:2010.

EMC-Directive 2014/30/EU according to the following standard(s):
EN 61000-6-1:2007, EN 61000-6-3:2007/A1:2011.

RoHS 2011/65/EU according to the following standard(s):
EN 50581:2012.

Supplementary Information The equipment complies with the American standards:
UL508.

The above has been declared according to the global method, module A

Date: 23.02.2016

Christian Skjold Heyde,
Vice President, R & D and Production, Struers ApS

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English

Declaration of Conformity

**Manufacturer,
responsible for
Technical File**

Struers A/S
Pederstrupvej 84
DK-2750 Ballerup, Denmark
Telephone +45 44 600 800

Herewith declares that

<i>Product Name:</i>	LaboPol-25
<i>Type No:</i>	586
<i>Machine Type:</i>	Grinding and polishing machine

is in conformity with the provisions of the following directives:

Safety of Machinery

2006/42/EC according to the following standard(s):
EN ISO 12100:2011, EN ISO 13849-1:2008/AC:2009, EN ISO 13849-2:2014,
EN 60204-1:2006/AC:2010, EN ISO 13857:2008, EN 349 + A1:2010, EN 1037 + A1:2010.

EMC-Directive

2004/108/EC according to the following standard(s):
EN 61000-6-1:2007, EN 61000-6-3:2007/A1:2011.

RoHS

2011/65/EU according to the following standard(s):
EN 50581:2012.

**Supplementary
Information**

The equipment complies with the American standards:
UL508.

The above has been declared according to the global method, module A

Date: 13.02.2015

Christian Skjold Heyde,
Vice President, R & D and Production, Struers A/S

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Some of the drawings may contain position numbers
not used in connection with this manual.

**Always state *Serial No* and *Voltage/frequency*
if you have technical questions or when ordering spare parts.**

The following restrictions should be observed, as violation of the restrictions may cause cancellation of Struers legal obligations:

Instruction Manuals: Struers Instruction Manual may only be used in connection with Struers equipment covered by the Instruction Manual.

Service Manuals: Struers Service Manual may only be used by a trained technician authorised by Struers. The Service Manual may only be used in connection with Struers equipment covered by the Service Manual.

Struers assumes no responsibility for errors in the manual text/illustrations. The information in this manual is subject to changes without notice. The manual may mention accessories or parts not included in the present version of the equipment. The contents of this manual is the property of Struers. Reproduction of any part of this manual without the written permission of Struers is not allowed.

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
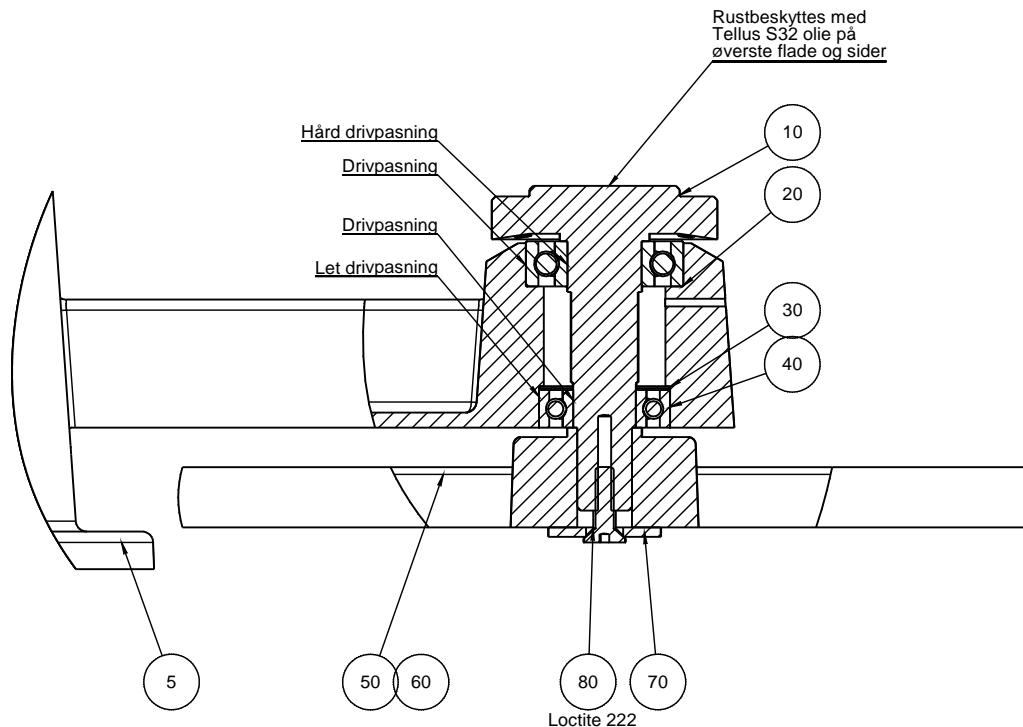
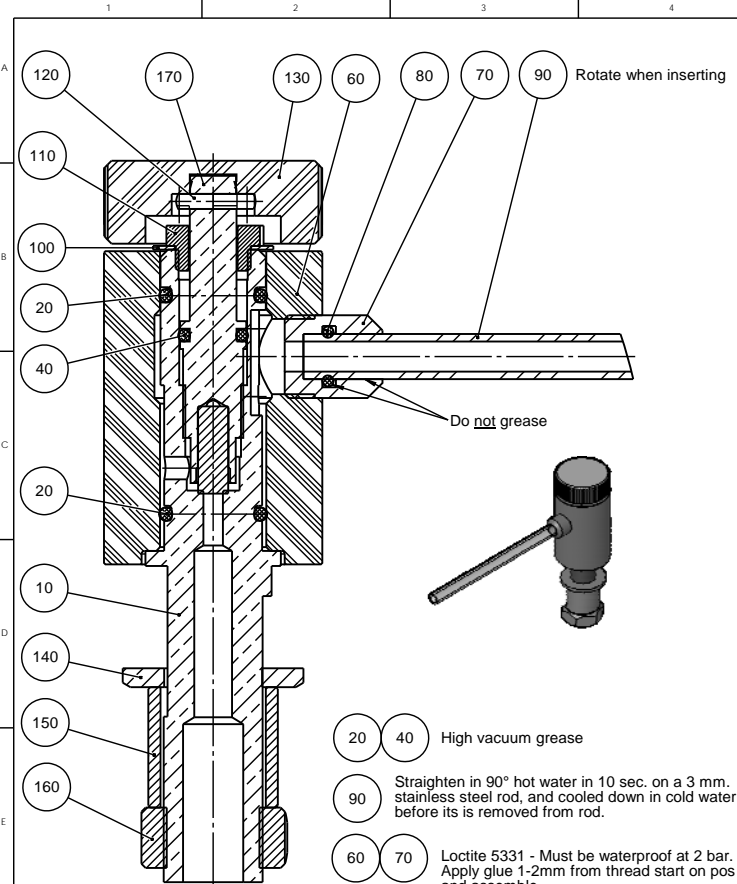
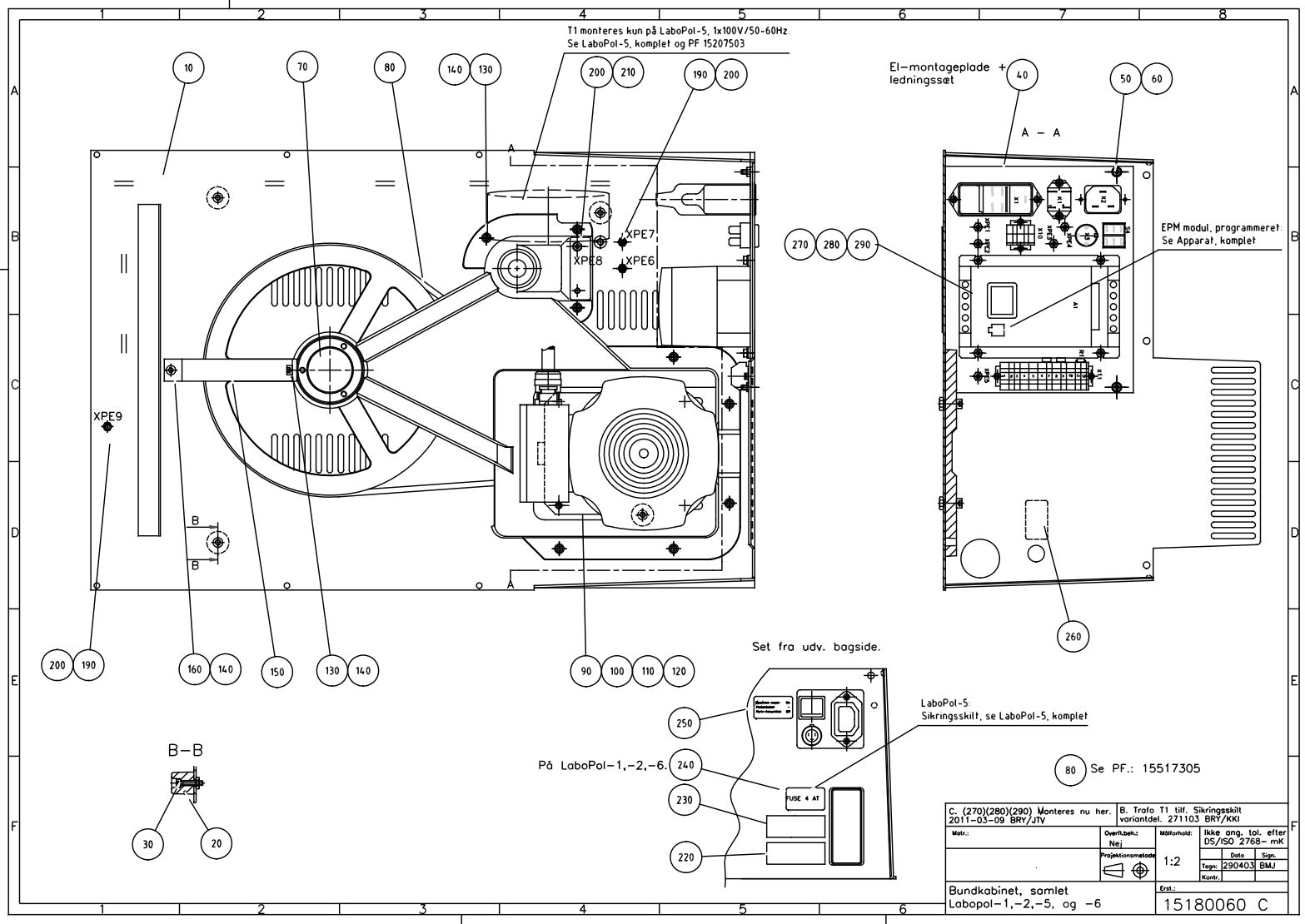
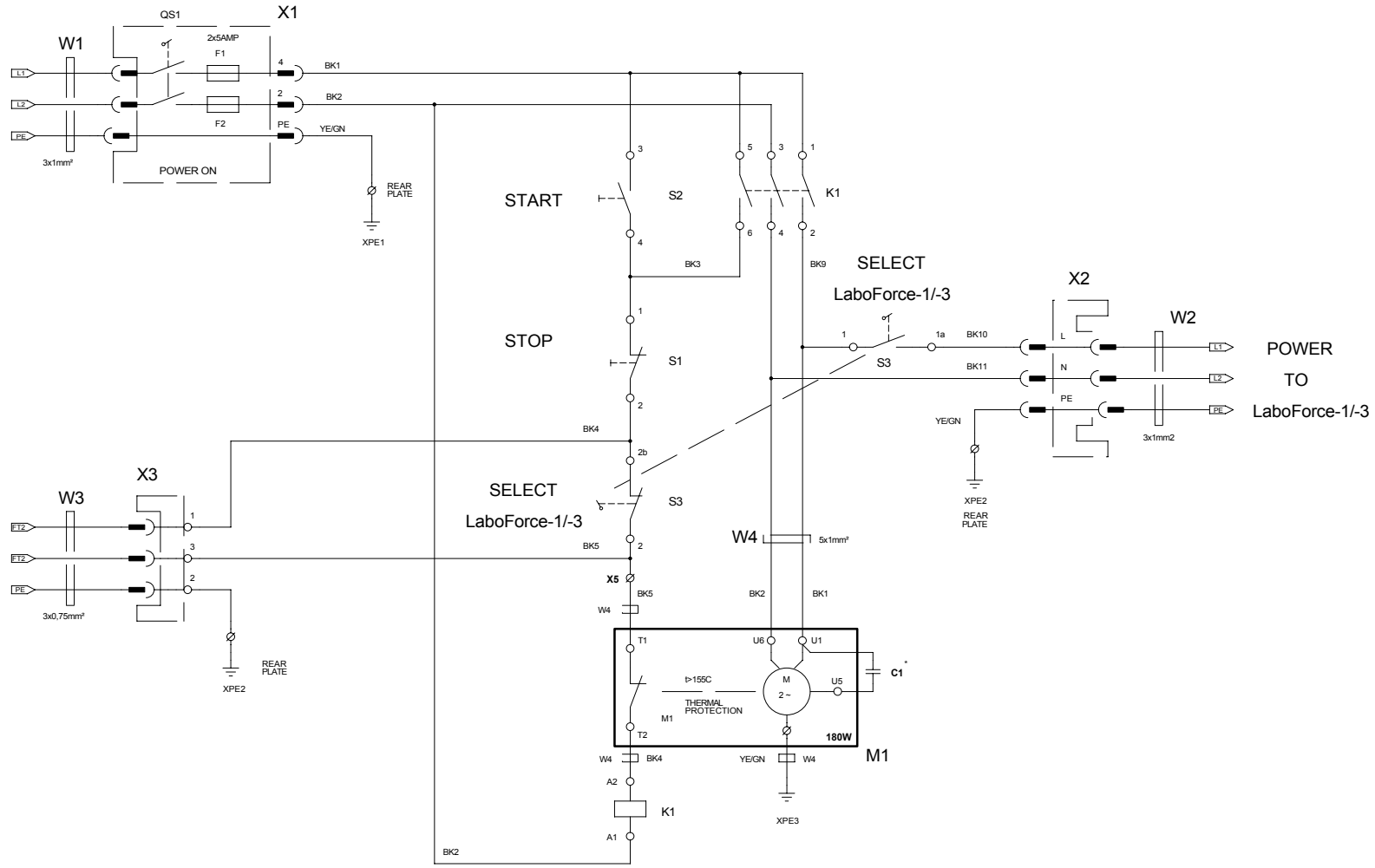
MUST BE SHORT CIRCUIT PROTECTED WITH EXTERNAL FUSES ACCORDING TO LOCAL REGULATIONS.
F = 16Amp MAXIMUM

FROM LaboForce-3
THERMAL PROTECTION M2
Comment:
Use Dummy for LaboForce-1:


ALL WIRINGS 0,75mm², EXCEPT OTHERWISE MARKED.

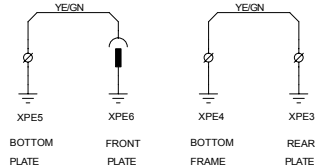- COLOR CODES:
- BK = BLACK
 - BN = BROWN
 - RD = RED
 - YE = YELLOW
 - GN = GREEN
 - BL = BLUE
 - GY = GREY
 - WH = WHITE



MAINS SUPPLY:		C1
100V/50Hz.		50µF
100V/60Hz.		40µF
110-120V/60Hz.		30µF
220-240V/50Hz.		10µF
220-240V/60Hz.		8µF


STRUERS A/S VALBOESALLE 178 DK-2610 ROEDOVRE DENMARK PHONE: +45 3670 3500		CIRCUIT DIAGRAM		LaboPol-21	
Thursday, September 13, 2001	Size A2	CAGE CODE	DWG NO	15293100	Rev
Monday, December 11, 2000	SCALE	DEM/DEM	Sheet	1	of 1

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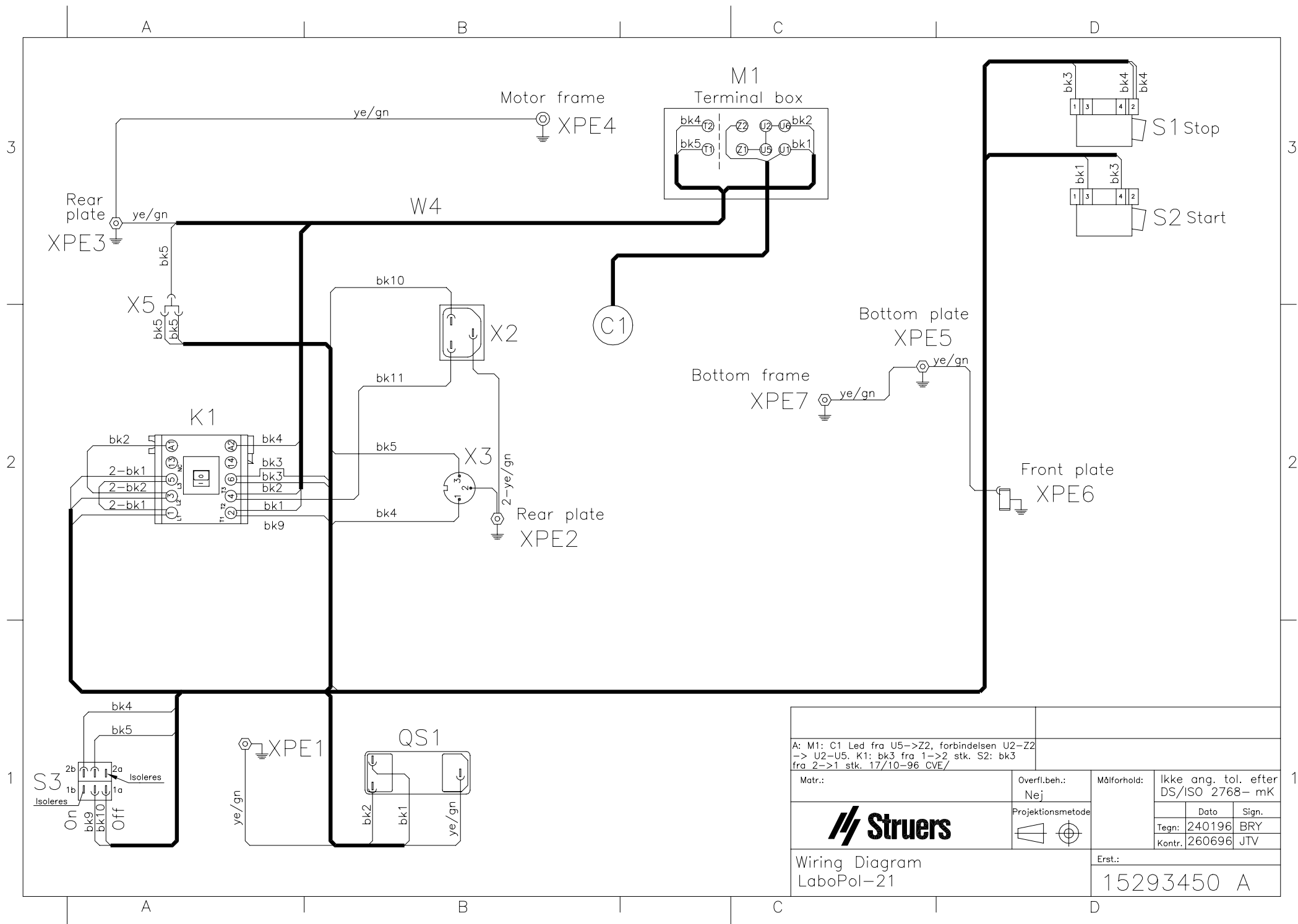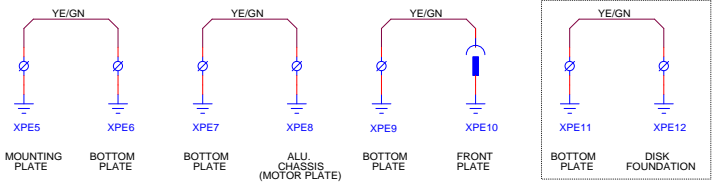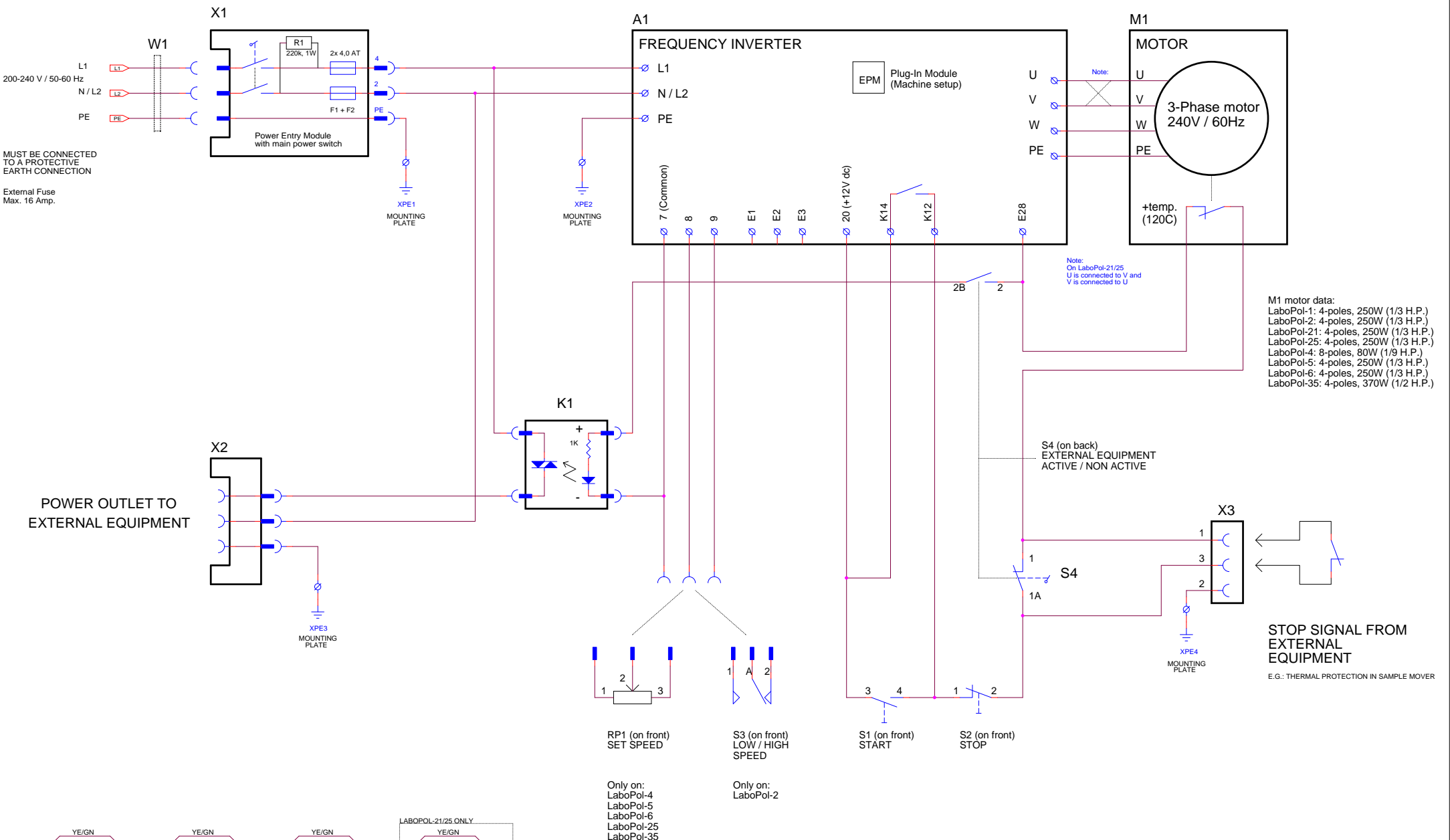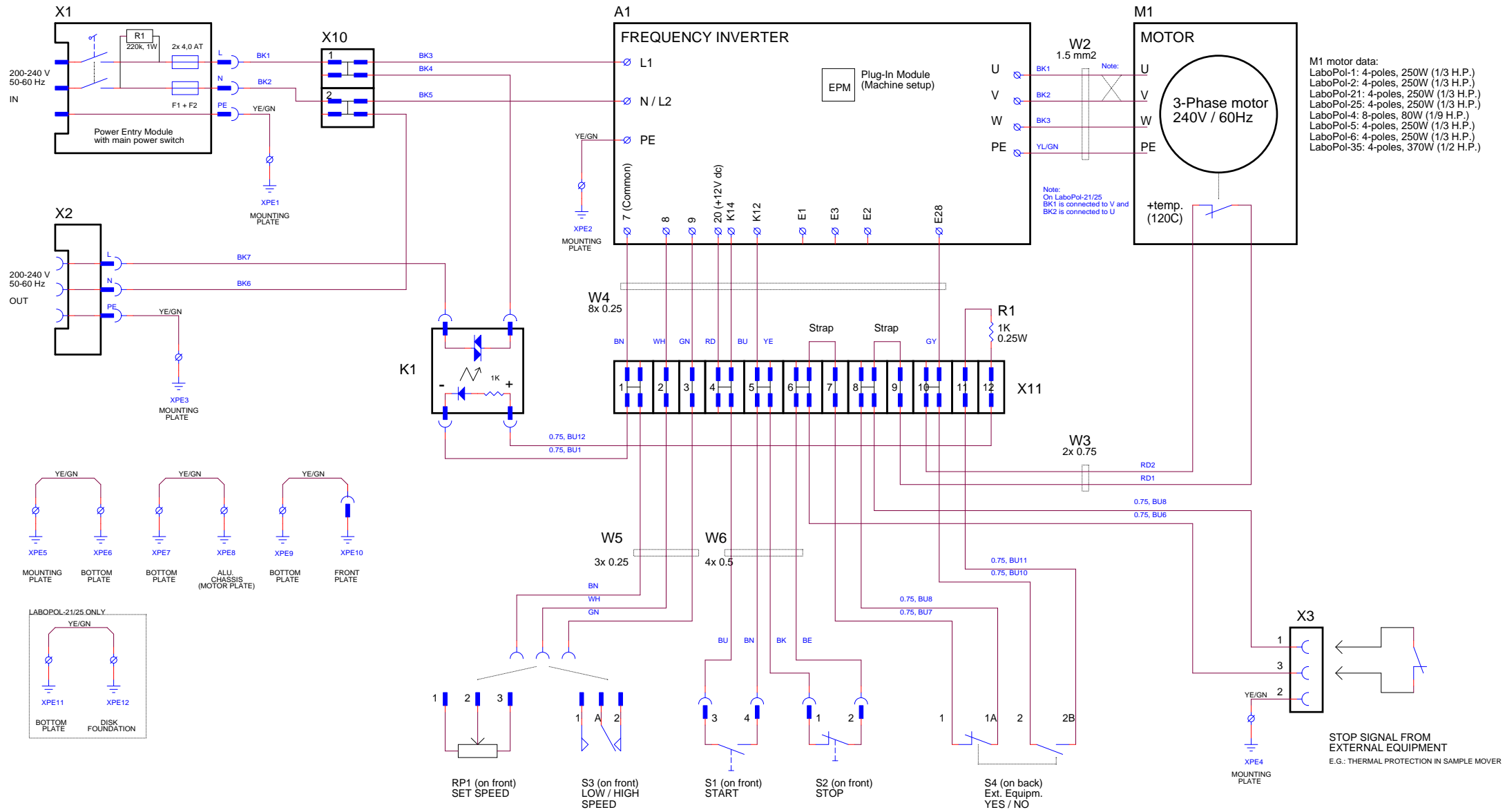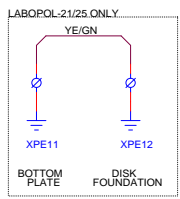
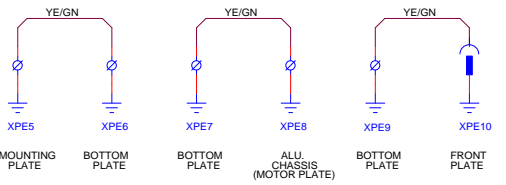
ALL WIRINGS AWG-16, EXCEPT OTHERWISE MARKED.

COLOR CODES:
 BK = BLACK
 BN = BROWN
 RD = RED
 OG = ORANGE
 YE = YELLOW
 GN = GREEN
 BU = BLUE
 VT = VIOLETT
 GY = GREY
 WH = WHITE
 BE = BEIGE

Only on:
 LaboPol-4
 LaboPol-5
 LaboPol-6
 LaboPol-25
 LaboPol-35

Only on:
 LaboPol-2

Rev.E: SPE 2012.05.29
 W2 1 mm2->1.5 mm2

Rev.B: FTH 10-07-2003 VALHOEJS ALLE 176 DK-2610 ROEDOVRE DENMARK PHONE: +45 3670 3500	STRUERS A/S VALHOEJS ALLE 176 DK-2610 ROEDOVRE DENMARK PHONE: +45 3670 3500	
Rev.C: FTH 15-02-2008 R1 resistor added in Power Switch X1	LABOPOL-SERIES, WIRING DIAGRAM	
Rev.D: FTH 28-08-2009 LaboPol-35 added	Size A2	CAGE Code <Cage Code>
Tuesday, May 29, 2012	Scale	DWG NO 15863450
		Rev E
		Sheet 1 of 1

Pederstrupvej 84
DK-2750 Ballerup
Denmark

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All rights reserved. © Struers 2015