# LaboPress-3

## Instruction Manual

Manual No.: 15087003

Date of Release 11.01.2008

*LaboPress-3*  
*Instruction Manual*

<b>Table of Contents</b>	<b>Page</b>
User's Guide .....	1
Reference Guide.....	13
Quick Reference Card .....	24

---

Always state *Serial No* and *Voltage/frequency* if you have technical questions or when ordering spare parts. You will find the Serial No. and Voltage on the type plate of the machine itself. We may also need the *Date* and *Article No* of the manual. This information is found on the front cover.

The following restrictions should be observed, as violation of the restrictions may cause cancellation of Struers legal obligations:

**Instruction Manuals:** Struers Instruction Manual may only be used in connection with Struers equipment covered by the Instruction Manual.

**Service Manuals:** Struers Service Manual may only be used by a trained technician authorised by Struers. The Service Manual may only be used in connection with Struers equipment covered by the Service Manual.

Struers assumes no responsibility for errors in the manual text/illustrations. The information in this manual is subject to changes without notice. The manual may mention accessories or parts not included in the present version of the equipment.

The contents of this manual is the property of Struers. Reproduction of any part of this manual without the written permission of Struers is not allowed.

All rights reserved. © Struers 2008.

**Struers A/S**  
Pederstrupvej 84  
DK-2750 Ballerup  
Denmark  
Telephone +45 44 600 800  
Fax +45 44 600 801

---

## **LaboPress-3** **Safety Precaution Sheet**

### **To be read carefully before use**

1. The operator should be fully aware of the use of the machine according to the Instruction Manual.
2. The machine must be placed in a well ventilated room on a working table with adequate height for convenient operation.
3. Be sure that the actual voltage corresponds to the voltage stated on the back of the machine and on the heating/cooling unit. The machine must be earthed.
4. Be sure that the water connections are mounted correctly and without leaks. The main water supply should be turned on when the machine is in use.  
Struers recommend that the mains water supply is shut off or disconnected if the machine is to be left unattended
5. Be sure that the outlet hose is safely attached to the water outlet system.
6. Be sure that the mounting unit is correctly assembled on the press before starting the process.
7. Be sure that the top closure with upper ram is either correctly mounted on the mounting cylinder or completely removed from the mounting cylinder before starting the press.
8. Do not operate the mounting press with a higher force/pressure than recommended for the actual cylinder diameter and resin in Struers Application Guide for Hot Mounting.
9. Following a heating cycle, ensure the mounting cylinder is cooled for a minimum of two minutes before opening.
10. Do not operate the machine whilst assembling or disassembling the mounting unit.

---

The equipment is designed for use with consumables supplied by Struers. If subjected to misuse, improper installation, alteration, neglect, accident or improper repair, Struers will accept no responsibility for damage(s) to the user or the equipment.

Dismantling of any part of the equipment, during service or repair, should always be performed by a qualified technician (electromechanical, electronic, mechanical, pneumatic, etc.).

---

# User's Guide

## Table of Contents

	Page
<b>1. Getting Started</b>	
Checking the Contents .....	2
LaboPress-3.....	2
Mounting Unit .....	2
Unpacking LaboPress-3.....	2
Placing LaboPress-3 .....	2
Transport Screw Cap.....	2
Getting Acquainted with LaboPress-3.....	3
Noise Level.....	4
Supplying Power .....	4
Supplying Water .....	4
Water Inlet.....	4
Water Outlet.....	4
Assembling the Mounting Unit.....	5
Removing the Cover .....	5
Installing the Lower Ram .....	5
Installing the Mounting Unit.....	5
Installing the Cover.....	6
Installing the Dust Protection Ring .....	6
Disassembling the Mounting Unit.....	6
Dropping the Lower Ram .....	6
Removing the Dust Protection Ring .....	6
Removing the Cover .....	6
Removing the Mounting Unit .....	7
Removing the Lower Ram .....	7
Changing the Mounting Unit.....	7
<b>2. Basic Operations</b>	
Using the Controls.....	8
Front Panel Controls of LaboPress-3 .....	8
Acoustic Signals.....	8
Main Switch .....	8
Placing the Specimen .....	10
Pouring Resin over the Specimen.....	10
Placing Two Specimens .....	10
Installing the Top Closure.....	11
Starting the Mounting Process .....	12
Preheating with Low Pressure .....	12
Stopping the Mounting Process .....	12
Removing the Top Closure .....	12

## 1. Getting Started

### Checking the Contents

*LaboPress-3*

In the packing box you should find the following parts:

- 1 LaboPress-3 machine
- 1 Pressure hose
- 1 Filter gasket
- 1 Gasket
- 1 Reduction ring with gasket
- 2 Measuring spoons for mounting resin
- 1 Funnel
- 1 Air filter
- 1 Set of Instruction manuals

*Mounting Unit*

- 1 Mounting unit
- 1 Top closure with upper ram
- 1 Lower ram
- 1 Dust protection ring
- 1 Piston pin
- 1 Mould release agent (FASTI)
- 1 Scraper (PROAN)

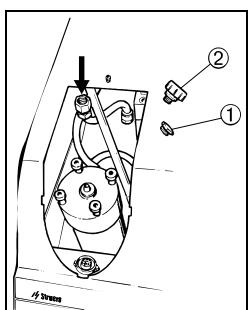
### Unpacking LaboPress-3

LaboPress-3 is detached from the bottom of the packing case by removing the four screws from below.

### Placing LaboPress-3

LaboPress-3 should be placed on a steady table with an adequate working height. The machine must be placed close to the power supply, water mains and water outlet facilities. If water recirculation is used, there must be room under the table for the Recirculation Cooling Unit (TRECI).

### Transport Screw Cap


Carry out the following procedure before using the machine for the first time. Exchange the transport screw cap ① with the enclosed air filter ②, to equalise the pressure in the hydraulic system.

The transport screw cap ① is situated underneath the cover for the mounting unit.

*LaboPress-3*  
*Instruction Manual*

**Getting Acquainted with  
LaboPress-3**

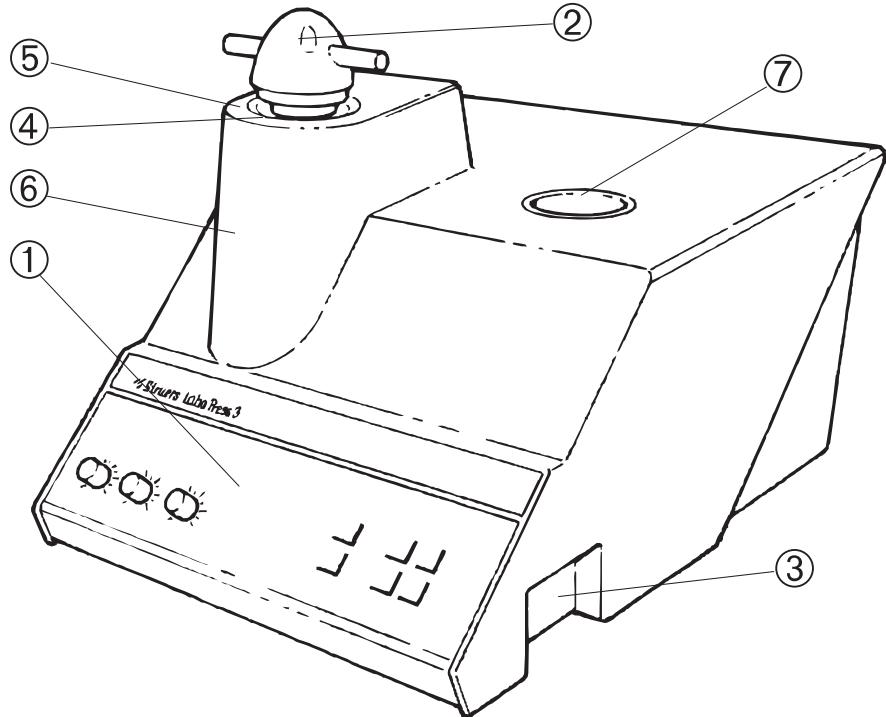Take a moment to familiarise yourself with the location and names of the LaboPress-3 components.

- ① Front panel
- ② Top closure
- ③ Main switch
- ④ Mounting unit
- ⑤ Dust protection ring
- ⑥ Cover for mounting unit
- ⑦ Top closure holder

## Noise Level

The noise level of the machine is 60 dB (A) measured when the pump is running, at a distance of 1.0 m/39.4" from the machine.

## Supplying Power

LaboPress-3 is factory mounted with an electric cable. Mount a plug on the cable:

Brown: phase  
Blue: neutral  
Yellow/green: earth

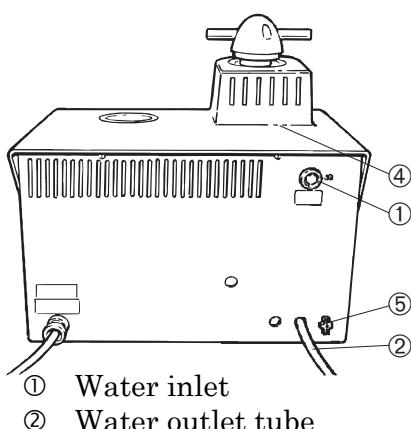
***IMPORTANT***

Check that the mains voltage corresponds to the voltage stated on the type plate on the machine. Do not operate the machine before the transit screw cap has been exchanged with the enclosed air filter.

## Supplying Water

Water for cooling may be supplied from the water mains or a Recirculation Cooling Unit.

### Water Inlet

Mount the pressure hose onto the water inlet tube (pos.1) on the back of the machine:

- Insert the filter gasket in the coupling nut with the flat side against the pressure hose.
- Tighten the coupling nut completely.

Mount the other end of the pressure hose on the water mains tap for cold water:

- Mount the reduction ring with gasket on the water mains tap, if necessary.
- Introduce the gasket and tighten the coupling nut completely.

***IMPORTANT***

Only connect to cold water

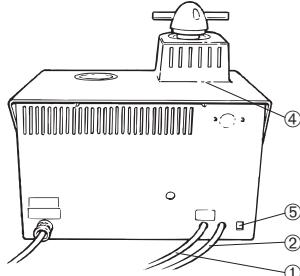
### Water Outlet

- Lead the water outlet tube (pos. 2) to the drain and be absolutely sure to place the hose with a steady downward slope and without any obstructions.
- Attach the water outlet tube to the water outlet system.

## Assembling the Mounting Unit

### Removing the Cover

- Press the catch (pos. 4).
- Press gently on both sides of the cover to release it, and lift it up.

④ Catch


### Installing the Lower Ram

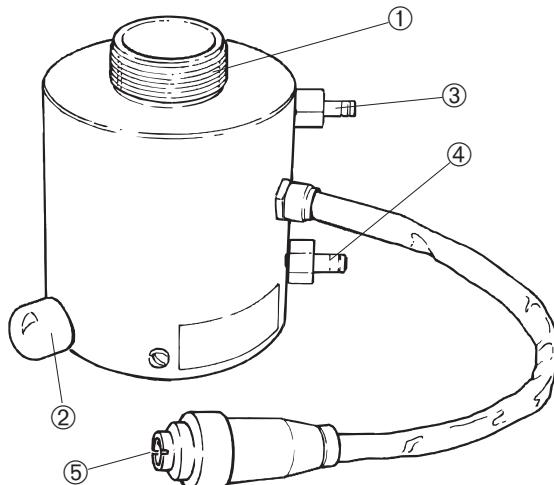- Place the lower ram on the top of the piston rod.
- Turn the lower ram so that the holes in the axle journal on top of the rod and ram line up, and insert the piston pin.
- Make sure that the ends of the pin do not protrude.

**IMPORTANT**

The axle journal is part of the machine's safety mechanism to protect against damage to the machine. Please see: Maintenance. If it breaks, it must only be replaced with a Struers replacement part or the safety mechanism may not function.


### Installing the Mounting Unit

- ① Mounting cylinder
- ② Fixation screw
- ③ Upper quick coupling for cooling water
- ④ Lower quick coupling for cooling water
- ⑤ Plug


**IMPORTANT**

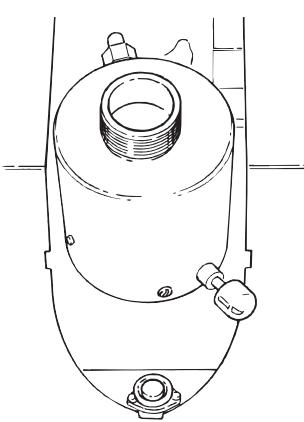Do not operate the machine whilst installing the mounting unit

- Unscrew the fixation screw about 10 mm.
- Place the mounting cylinder over the lower ram with the fixation screw in the position shown.
- Turn the mounting unit in a clockwise direction until it stops.
- Tighten the fixation screw completely.
- Mount the tube with the straight quick coupling on the lower quick coupling (position 4) of the mounting unit and push to connect. Make sure that the ring comes to stop in the extreme end of the quick coupling.
- Mount the tube with the elbow quick coupling on the upper quick coupling of the mounting unit (pos. 3) and push to connect. Make sure that the ring comes to stop in the extreme end of the quick coupling.
- Mount the plug (pos. 5) in the socket. Tighten the coupling nut.

#### *Installing the Cover*

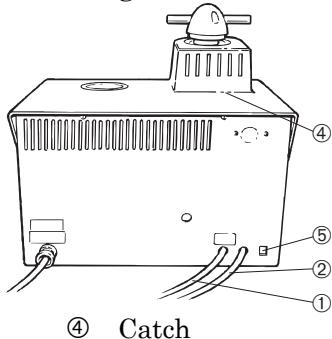- Insert the barb into the front of the opening for the cover.
- Insert the barbs on both sides of the cover into the cabinet. Press the sides of the cabinet gently.
- Press the rear end of the cover in so that the catch engages.

#### *Installing the Dust Protection Ring*

Place the dust protection ring around the mounting cylinder. The concave side should face upwards.

### **Disassembling the Mounting Unit**

#### *Dropping the Lower Ram*

- Switch on the main power of the machine.
- Press the key RAM DOWN ▼ to lower the ram to its lower limit.

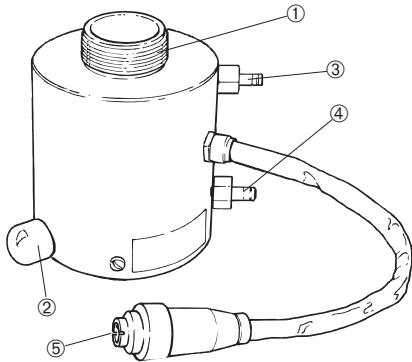
#### *Removing the Dust Protection Ring*

Lift out the dust protection ring.

#### *Removing the Cover*

- Press the catch (position 4 above).
- Press gently on both sides of the cover, to release it, and lift it up.

## Removing the Mounting Unit

**IMPORTANT**

Do not operate the machine whilst removing the mounting unit

- ① Mounting cylinder
  - ② Fixation screw
  - ③ Upper quick coupling for cooling water
  - ④ Lower quick coupling for cooling water
  - ⑤ Plug
- Disconnect the plug (pos. 5) in the socket. Loosen the coupling nut and pull.
  - Disconnect the tube with elbow quick coupling on the upper quick coupling on the mounting unit (pos. 3). Pull the ring. Wait 5 seconds to allow the water to flow out of the cooling coil.
  - Disconnect the tube with straight quick coupling on the lower quick coupling on the mounting unit (pos. 4). Pull the ring.
  - Unscrew the fixation screw about 10 mm.
  - Turn the mounting unit in an anti-clockwise direction until it stops.
  - Lift the mounting unit.

## Removing the Lower Ram

- Push the piston pin out of the lower ram and lift it off.

**IMPORTANT**

The axle journal is part of the machine's safety mechanism to protect against damage to the machine. Please see: Maintenance. If it breaks, it must only be replaced with a Struers replacement part or the safety mechanism may not function.

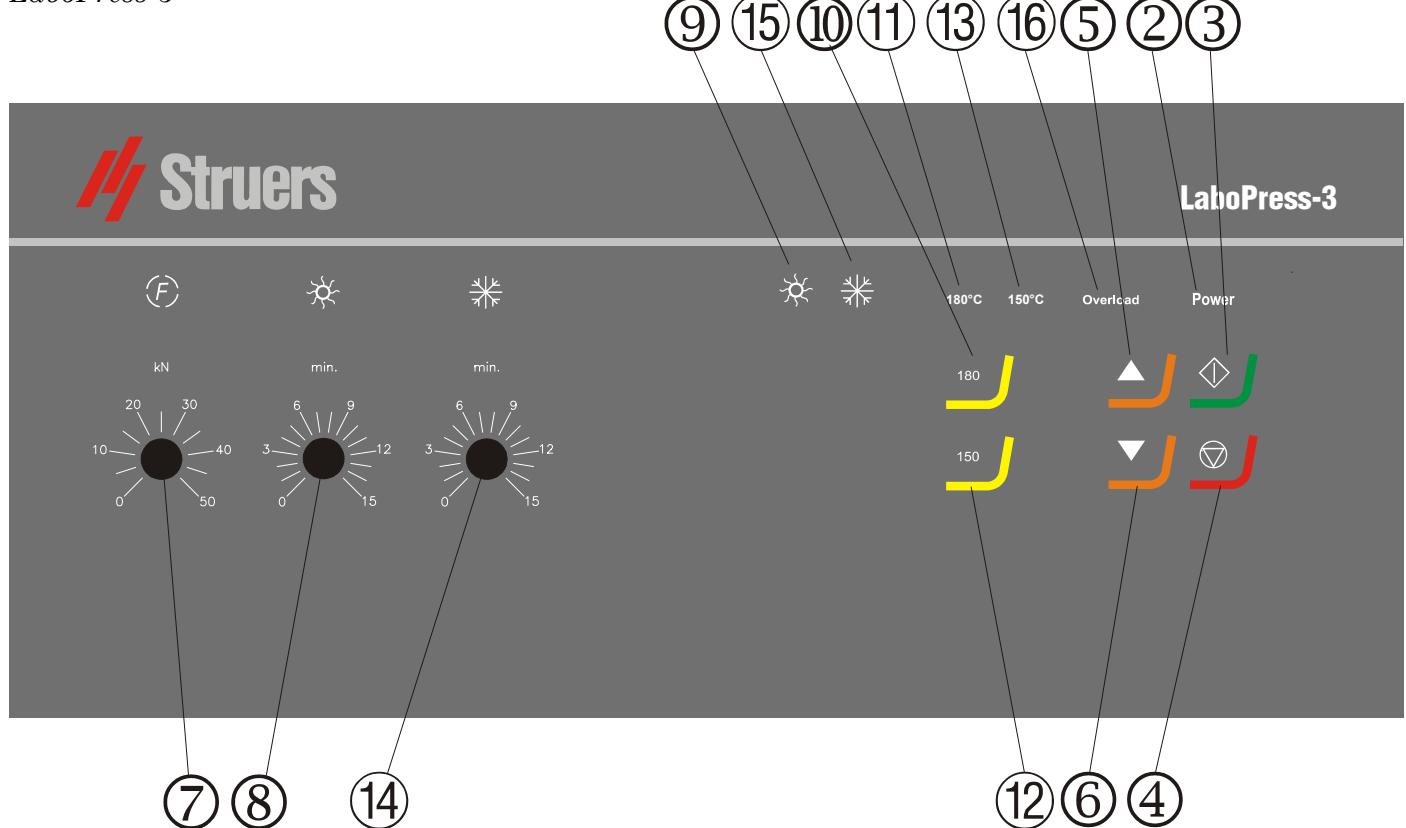
## Changing the Mounting Unit

Follow the instructions for Removing the Mounting Unit and Installing the Mounting Unit.

## 2. Basic Operations

### Using the Controls


Front Panel Controls of  
LaboPress-3

### Acoustic Signals

- ⌚ *Short Beep:* when a key is pressed, a short beep indicates that the command has been accepted.
- ⌚ *Long Beep:* a long beep indicates that the key is inactive at that moment.
- ⌚ *Three long beeps:* indicate that the mounting process is finished.
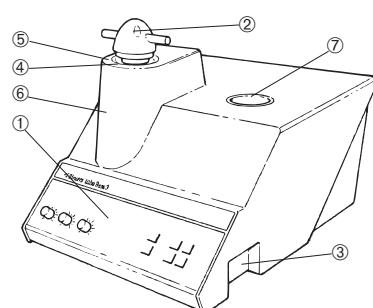
### Main Switch

- ① Front panel
- ③ Main switch

*LaboPress-3*  
*Instruction Manual*

Pos. No.	Key/Knob	Function	Pos. No.	Key/Knob	Function
① MAIN SWITCH		Turns the main power of the machine on/off. The main switch is located to the right of the cabinet.	⑨ HEATING INDICATOR		Lights up when heating is on.
② POWER		Lights up when the main power is on.	⑩ HEATING		Selects a heating temperature of 180°C.
③ START		Starts the automatic mounting process.	⑪ HEATING INDICATOR		Lights up when the heating temperature is set to 180°C.
④ STOP		Stops the mounting process, followed by automatic release of pressure. Stops the movement of the lower ram.	⑫ HEATING		Selects a heating temperature of 150°C.
⑤ RAM UP		Starts the upward movement of lower ram. When started: stops the upward movement of the lower ram. The ram automatically stops when the upper limit is reached.	⑬ HEATING INDICATOR		Lights up when the heating temperature is set to 150°C.
⑥ RAM DOWN		Starts the downward movement of lower ram. When started: Stops the downward movement of the lower ram. The ram automatically stops when the lower limit is reached.	⑭ COOLING		Adjustable knob for setting the cooling time.
⑦ FORCE		Adjustable knob for setting the force (pressure).	⑮ COOLING INDICATOR		Lights up when cooling is on.
⑧ HEAT		Adjustable knob for setting the heating time.	⑯ OVERLOAD		Lights up if the motor has been overloaded.

## **Placing the Specimen**

- Press and hold RAM UP ▲ to raise the lower ram to its upper limit.
- Apply "*Mould Release Agent*" to the surface of the lower ram.
- Place the specimen on the ram. The specimen must be clean, dry and free from grease. The distance between the specimen and the cylinder wall must be minimum 3 mm to avoid cracks in the resin.

## *Pouring Resin over the Specimen*

- Press and hold RAM DOWN ▼ to lower the ram to its lower limit.
- Fill a suitable amount of resin into the cylinder by means of the enclosed funnel.

**Always** make sure there is sufficient resin to cover the sample after compression. Please note that the volume of the resin reduces when the granulate becomes compressed. If insufficient resin is used, the rams may come in contact with the sample, and the cylinder may be damaged.

### ***Important***

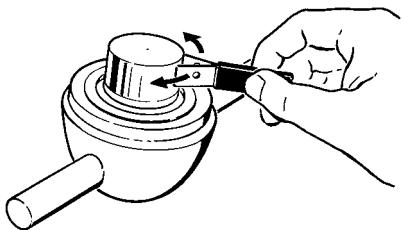
The "*Mould Release Agent*" must always be applied to the mounting rams as a thin layer to prevent the resins sticking to the surface. By means of Struers AntiStick a thin layer of stearate powder can easily be dabbed on the rams.

## **Placing Two Specimens**

- Follow the instructions in Placing the Specimen.
- Apply "*Mould Release Agent*" to all surfaces of the intermediate ram.
- Place the intermediate ram on top of the resin.
- Place the second specimen on the intermediate ram and fill the cylinder with an adequate quantity of resin.

Again, make sure you use adequate resin!

### Installing the Top Closure

- Remove resin dust from the upper part of the mounting cylinder.
- Clean the cylindrical surface of the upper ram. Cured resin can easily be removed without damage to the surface of the ram using the scraper supplied.
- Apply "*Mould Release Agent*" to all accessible surfaces of the upper ram.
- Place the top closure with the upper ram on the mounting cylinder.
- Press the top closure straight down, turning it counter-clockwise until you hear a click.
- Press the top closure down and turn it clockwise until its lower limit.
- Turn the top closure a quarter of a turn back.

***Important***

If the ram does not fit in the cylinder then check ram and cylinder for cured resin. The tolerance between the cylinder and the ram is very fine and even small amounts of resin from previous mounting may cause problems.

### **Starting the Mounting Process**

- Set the force.
- Set the heating time.
- Set the heating temperature (180/150°C).
- Set the cooling time.
- Press START , and the process will run automatically.
- The mounting process can be followed on the heating and cooling indicators.

### *Preheating with Low Pressure*

Set the force to a low value to begin with. Adjust the force after the desired period of preheating.

### **Stopping the Mounting Process**

- The machine automatically stops and relieves the pressure when the cooling time has elapsed. The machine can be stopped at any time during the mounting process by pressing STOP .

*If you have stopped the machine during the mounting process:*

Cool the mounting cylinder in minimum 2 min before opening, after a heating period. Please note that the mount possibly will be destroyed.

### **Removing the Top Closure**

When the mounting process is finished:

- Turn the top closure counter-clockwise until released from the threads.
- Press RAM UP  to raise the lower ram to its upper limit.
- Place the top closure in the top closure holder.

# Reference Guide

## Table of Contents

## Page

### **1. Advanced Operations**

Installation of Recirculation Cooling Unit (Optional) .....	14
Socket for Recirculation .....	14

### **2. Struers Metalog Guide™ .....**

15

### **3. Application Guide for Hot Mounting .....**

15

### **4. Accessories .....**

16

### **5. Consumables .....**

17

### **6. Trouble-Shooting .....**

18

### **7. Maintenance**

Daily Service .....	20
Monthly Service .....	20
Cleaning under the Lower Ram.....	20
Cleaning the Rams .....	20
Removing the Upper Ram .....	21
Lubricating the Threads for the Top Closure.....	21
Decalcifying the Cooling Coil in the Mounting Unit.....	21
Replacing the Cooling Water.....	22
Checking the Recirculating Cooling Unit.....	22

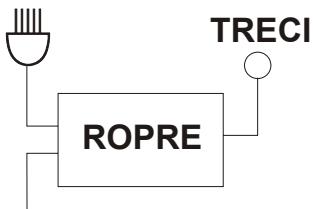
### **8. Technical Data .....**

23

## 1. Advanced Operations


### Installation of Recirculation Cooling Unit (Optional)

A Recirculation Cooling Unit (TRECI) can be connected by means of a Recirculation Cooling Unit Connector (ROPRE).

LaboPress-3/  
ProntoPress-10/ProntoPress-20

#### Socket for Recirculation

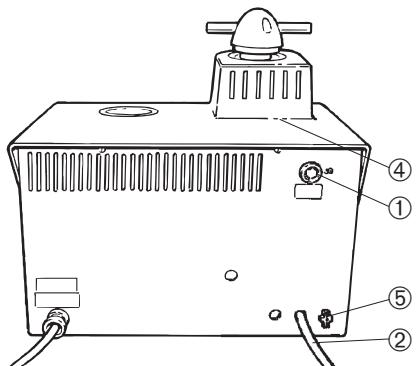- ② Outlet hose
- ⑤ Socket for recirculation

- Connect the cable with plug from ROPRE to the socket (pos. 5) on the back of the mounting press.
- Connect the cable from ROPRE to the mains.
- Check that the pump rotates in the direction indicated on the pump. If the direction is incorrect, switch two of the phases in the connection to the pump in TRECI (3-phase versions only).
- Replace the outlet tube on the pump in TRECI with the non-return valve with elbow and hexagon nipple from ROPRE. Struers recommends that you seal the threads.
- Connect the pressure hose of the mounting press to the hexagon nipple. Remember to mount the gasket.
- Lead the outlet hose (pos. 2) to the inlet on TRECI. Be absolutely sure to place the hose with a steady downward slope and without obstructions.
- Attach the water outlet to the inlet of TRECI.

## **2. Struers Metalog Guide™**

Struers Metalog Guide™ offers preparation methods for the most common materials, based on a simple analysis of two key properties: hardness and ductility. Finding the right method, including choice of consumables is easy. Always consult Struers Metalog Guide™ for the correct preparation method for the actual specimens.

Struers Metalog Guide™ contains 6 useful chapters:

A quick and safe guide to the right preparation method.

*Metalogram*

*Metalog Method:*

A complete catalogue of preparation methods, based on Struers' extensive materialographic experience, and employing Struers' range of consumables.

*Preparation Philosophy*

The basics of modern specimen preparation, seen from a professional point of view.

*Metalog Process*

The materialographic preparation process from start to finish, logically explained.

*Metalog Master*

A combined trouble-shooting guide and supply of in-depth information on the processes of mechanical preparation, including an expert system for the solving of preparation problems.

*Metalog Code*

Quick access to the relevant consumables for the chosen preparation methods.

### **Struers Metalog Guide™**

A complete guide to materialographic specimen preparation. Contact your local dealer for a free copy of Metalog Guide™.

## **3. Application Guide for Hot Mounting**

Useful mounting data and hints can be found in the Struers [Application Guide for Hot Mounting](#) or visit the Struers website on <http://www.struers.com> and see the section on **Knowledge**

## 4. Accessories

Type	Description	Code
<b>Mounting Unit</b> consisting of cylinder, upper and lower ram, heating/cooling unit and top closure	25 mm dia. 30 mm dia. 1 ¼" dia. 1 ½" dia. 40 mm dia. 50 mm dia.	CYNMI CYNTE CYNQU CYNHA CYNIF CYNAM
<b>Intermediate Ram</b> for making 2 simultaneous mounts	25 mm dia. 30 mm dia. 1 ¼" dia. 1 ½" dia. 40 mm dia. 50 mm dia.	CYLDO CYLDO CYLDO CYLDO CYLDO CYLDO
<b>Recirculation Cooling Unit</b>		TRECI
<b>Recirculation Cooling Unit Connector</b> for connection of Recirculation Cooling Unit (TRECI) to LaboPress-3.		ROPRE

## 5. Consumables

Resin	Application	Specific property	Material	Contents	Code
<b>ConduFast</b> (Resin 1)	Electropolishing	Electrically conductive Very low shrinkage	Acrylic resin with iron powder filler Thermoplastic, Grey	1 kg	RESFE
<b>SpeciFast</b> (Resin 3)	Glass clear mounts Porous samples Surface electrical insulator for ConduFast	Transparent Medium shrinkage	Acrylic Thermoplastic	1.5 kg 7.5 kg 25 kg	RESTH RESYV RESFA
<b>PolyFast</b>	Edge retention	Very low shrinkage Medium removal rate	Phenolic resin with carbon filler Thermosetting, Black	1 kg 7.5 kg	FAPSA FAPME
	Examination in scanning electron microscope	Electrically conductive Low emission in the vacuum chamber.			
<b>IsoFast</b> (Resin 4)	Edge retention Planeness	Low shrinkage Low removal rate	Dialylphthalate with glass fibre filler Thermosetting, Green	1 kg 7.5 kg	RESDI RESTY
<b>DuroFast</b> (Resin 5)	Edge retention Planeness	Very low shrinkage Good adhesion Very low removal rate	Epoxy with mineral filler Thermosetting, Black	1 kg 7.5 kg	RESIF RESEN
<b>MultiFast</b>	Routine examination, Backup resin	Low shrinkage Medium removal rate	Phenolic resin with wood flour filler Thermosetting, Black	2.5 kg 7.5 kg 25 kg	RESRU RESUK RESLA
<b>MultiFast Green</b>	Routine examination Colour marking	Low shrinkage Medium removal rate	Phenolic resin with wood flour filler Thermosetting, Green	2.5 kg 7.5 kg 25 kg	FAGSA FAGME FAGLA
<b>MultiFast Brown</b>	Routine examination Colour marking	Low shrinkage Medium removal rate	Phenolic resin with wood flour filler Thermosetting, Brown	2.5 kg 7.5 kg 25 kg	FABSA FABME FABLA

Product	Application	Specific property	Material	Contents	Code
<b>Pre-Mount</b>	Serial mounting of uncomplicated shapes	Easy to handle Low shrinkage Medium removal rate	Phenolic resin with wood flour filler In tablet form Thermosetting, Black	<b>25 mm/ 1" dia.</b> approx. 1100 pcs. 10 kg	RESUN
				<b>30 mm/ 1.1/4" dia.</b> approx. 760 pcs. 10 kg	RESGI
				<b>40 mm/ 1 1/2" dia.</b> approx. 460 pcs. 10 kg	RESAM

Product	Application	Material	Contents	Code
<b>Mount Release Wax</b>	To avoid sticking of the mounts to the rams	Wax	24 g	PREWA
<b>AntiStick</b>	To avoid sticking of the mounts to the rams	Very fine stearate powder	8 g	FASTI

## 6. Trouble-Shooting

Display/error	Cause	Action
<b>Error Message</b>		
Overload	The motor has been overloaded.	Let the motor cool down and start the mounting press again. If this does not help: Call a Struers service technician.
<b>Acoustic Signals</b>		
Long beep.	The command can not be accepted.	Check User's Guide, Section 2.
<b>Machine Problems</b>		
Process will not start.	Heating time and cooling time are set at zero.	Set the correct times.
Insufficient compression.	Incorrect setting of force/pressure.	Set the correct parameter.
	Defect in the pressure system.	Call a Struers service technician.
Insufficient heating.	Incorrect setting of time or temperature for heating.	Set the correct parameters.
	Defect in the heating system.	Call a Struers service technician.
Insufficient cooling.	Incorrect setting of time for cooling.	Set the correct parameter.
	Mains tap supplying cooling water is either closed or not opened sufficiently. Filter at the water inlet is blocked. The water is too hot.	Open the tap. Clean the filter.
	Insufficient water within the Recirculation Cooling Unit. The water is too hot.	Fill to the correct water level. See the section: Maintenance.
	Defect in the cooling system.	Call a Struers service technician.
Cooling water drips underneath the machine.	The quick coupling is not mounted correctly.	Remove the cover from the mounting unit and check the quick coupling connections.
The top closure will not start threading on the mounting cylinder.	The top closure is not mounted correctly.	Press the top closure straight down, turning it counter-clockwise until you hear a click. Turn the top closure clockwise.
	The upper ram is too hot.	Let the top closure and the upper ram cool down. Reduce the mounting temperature.
	Cured resin inside the mounting cylinder.	Clean the mounting cylinder with a brass wire brush.
	Cured resin on the cylindrical surface of the upper ram.	Clean the ram with the scraper supplied.
	The top closure has been dropped on the floor/table, causing a bulge in the edge of the ram.	Replace the upper ram.

*LaboPress-3*  
*Instruction Manual*

Error	Cause	Action
	Damage to the threaded connection in the top closure or to the upper ram.	Remove the upper ram from the top closure (see instructions in chapter Maintenance). Try both the following procedures to identify the problem: <ul style="list-style-type: none"> <li>▪ Try to mount the top closure, without the upper ram. If this is not possible, call a Struers service technician.</li> <li>▪ Try to insert the disconnected upper ram into the mounting cylinder. If this is not possible, call a Struers service technician.</li> </ul>
The top closure cannot be screwed down completely.	Dirt in the threads on the top closure and the mounting cylinder.	Clean the threads. Use only a dry lubricant powder.
	The disc for thermal insulation, placed on the top of the upper ram, has a larger diameter than the upper ram.	Call a Struers service technician.
The top closure can not be loosened.	The top closure has not been turned $\frac{1}{4}$ turn back before starting the process.  Cured resin on the cylindrical surface of the top ram.  Dirt in the threads of the top closure.	To release the top closure: <ul style="list-style-type: none"> <li>- Move the lower ram up and down several times.</li> </ul> If this does not help: <ul style="list-style-type: none"> <li>- Put heating on for 1 min. and try again.</li> </ul> If this does not help: <ul style="list-style-type: none"> <li>- Complete a mounting process without force, but with heating and cooling time of about 15 min.</li> </ul> If this does not help: <ul style="list-style-type: none"> <li>- Remove the two handles on the top closure.</li> <li>- Remove the plastic cover from the top closure (See Maintenance).</li> <li>- Loosen the top closure with a fork spanner.</li> </ul>

## 7. Maintenance

### Daily Service

- Clean all accessible surfaces with a moist cloth.

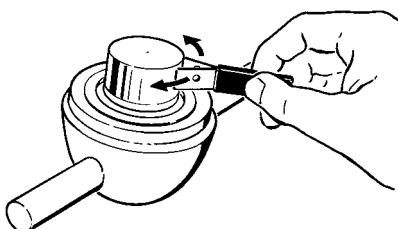
### Monthly Service

#### *Cleaning under the Lower Ram*

When mounting, some old cured resin will fall down from the lower ram and built up underneath. The cured resin must be cleaned away, to avoid it harming the machine. The axle journal securing the lower ram has been designed to break in this case, thus preventing serious damage to the machine.

- Remove the mounting unit. See this section.
- Remove the resin under the lower ram with a cloth or a soft brush.
- Install the mounting unit. See this section.

### Cleaning the Rams

Before each mounting, check that there is no resin left on the flat surface of the rams from previous operations. Clean the cylindrical surface of the upper ram and optional intermediate ram as this may cause problems with travel in the cylinder and mounting the top closure. Cured resin can easily be removed without any damage to the surface of the rams using the scraper supplied.


If the rams have been scratched on the polished surfaces, disassemble and polish them with 6 or 3 µm diamonds on a polishing cloth.

#### ***IMPORTANT***

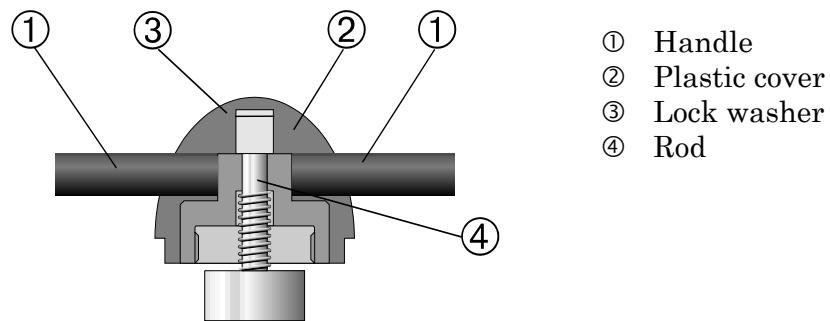Do not polish the surface on the sides of the rams. If the sides of the rams are heavily scratched, they must be scrapped.

## Removing the Upper Ram

- Take the cover plugs out of the ends of the handles and unscrew them with an allen key. (pos. 1).
- Remove the plastic cover (pos.2).
- Lift the upper ram and remove the lock washer (pos.3).
- Pull out the upper ram. Do not remove the rod (pos. 4) from the upper ram unless absolutely necessary.
- If it is necessary to remove the rod, hold the upper ram in a vice or a similar device. The ram surface must be protected with plastic or soft metal.

## Lubricating the Threads for the Top Closure

In case there is any friction in the threads, they should be lubricated with a dry lubricant such as MoS<sub>2</sub> powder or graphite.

### ***IMPORTANT***

Please observe that the threads in the joint between the top closure and the mounting cylinder must not be lubricated with oil or grease. As the mounting cylinder operates at high temperatures, only a dry lubricant, resistant to high temperatures, can be used.

## Decalcifying the Cooling Coil in the Mounting Unit

When using cooling water from the mains tap in areas with a high chalk or mineral content, deposits can build up in the cooling coil. This may reduce the cooling effect, so once a year the cooling coil should be decalcified.

- Remove the mounting unit. (See this section.)
- Flush the coil with a mild decalcifying fluid, such as that used for coffee machines.
- Flush the cooling coil with clean water.
- Re-install the mounting unit.

## Replacing the Cooling Water

*Only relevant when a Recirculation Cooling Unit is connected:*

- Replace the cooling water in the Recirculation Cooling Unit at least once a month.
- Place a Disposable Plastic Container (TREPO) in the tank and fold it over the edge.
- Fill the tank with 30 l of water. NB! Too high a level of coolant in the tank might damage the pump.
- Add Struers Additive according to the instructions.
- Put the lid on the tank.
- Place the pump in the lid

## Checking the Recirculation Cooling Unit

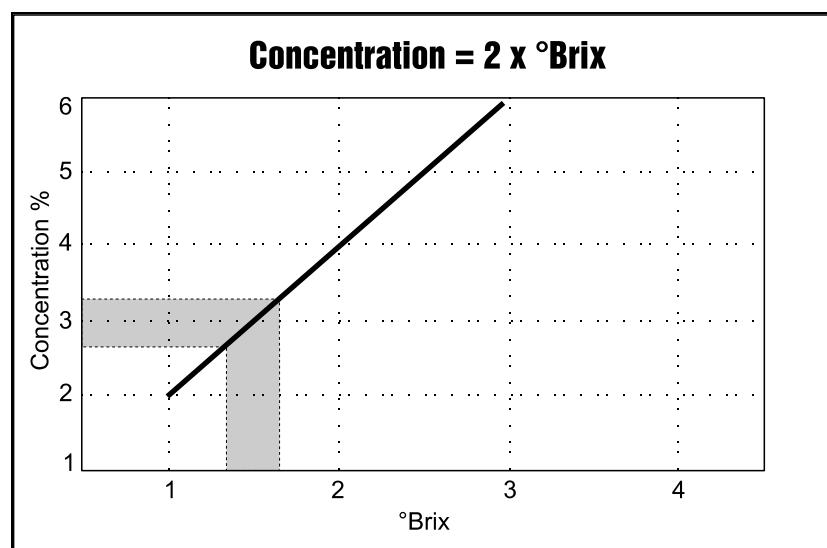
### ***IMPORTANT***

Always maintain the correct concentration of Struers Additive in the cooling water (percentage stated on the container of the Additive). Remember to add Struers Additive each time you refill with water.

The cooling unit should be checked for cooling water every week and refilled when the flushing pump cannot reach the cooling water. Remember to add Struers Additive for Cooling Fluid: One part of Additive for 33 parts of water.

The concentration of additive should always lie between 2.7 and 3.3%; the pH should be around 9. Replace the cooling water if pH is  $\leq 8.0$ .

To check the concentration of additive, use a refractometer (0-10 Brix, see diagram.). Concentration =  $2 \times$  Brix. Add Struers Additive for Cooling Fluid if the concentration is lower than 2.7%. Add water if the concentration is higher than 3.3%.

## 8. Technical Data

<b>Subject</b>	<b>Specifications</b>		
	<b>Metric/International</b>	<b>US</b>	
Mounting Units (Optional)	Diameter	25, 30, 40, 50 mm	1½" 1½"
Compression	Force on piston rod	Variable between 0 - 50 kN	Variable between 0 - 12000 Lbf
Heating (with pressure on)	Temperature	150/180 °C	302/356°F
	Time	Variable between 0 and 15 min	
Cooling (with pressure on)	Time	Variable between 0 and 15 min	
Water Supply	Tap Water or Recirculation Cooling Unit		
	Pressure for tap water	1 - 10 bar	14.5 - 145 psi
	Inlet	½" or ¾"	½" or ¾"
Supply Voltage	Power phases	1	
	Power consumption:		
	Max. Total Mounting Unit	1500W 750W	
	Voltage/frequency:	Max. continuous load	
	100-120V / 50Hz 100-120V / 60Hz 220-240V / 50Hz 220-240V / 60Hz	10A (100V) - 12A (120V) 10A (100V) - 12A (120V) 4.5A (220V)- 5.0A (240V) 4.5A (220V)- 5.0A (240V)	
Dimensions and Weight	Width	420 mm	
	Depth	580 mm	
	Height (Mounting Unit and Top Closure installed)	420 mm	
	Weight	38 kg	

# Quick Reference Card

## Placing the Specimen

- Press and hold RAM UP ▲ to raise the lower ram to its upper limit.
- Apply “*Mould Release Agent*” to the surface of the lower ram.
- Place the specimen on the ram.

## Pouring Resin over the Specimen

- Press and hold RAM DOWN ▼ to lower the ram to its lower limit.
- Fill a suitable amount of resin into the cylinder through the funnel.

## Installing the Top Closure

- Remove resin dust from the upper part of the mounting cylinder.
- Clean the cylindrical surface of the upper ram. Cured resin can easily be removed without damage to the surface of the ram using the scraper supplied.
- Apply “*Mould Release Agent*” to all accessible surfaces of the upper ram.
- Place the top closure with the upper ram on the mounting cylinder.
- Press the top closure straight down, turning it counter-clockwise until you hear a click.
- Press the top closure down and turn it clockwise until its lower limit.
- Turn the top closure a quarter of a turn back.

## Starting the Mounting Process

- Set the force.
- Set the heating time.
- Set the heating temperature (180/150°C).
- Set the cooling time.
- Press START ◇, and the process will run automatically.

## Stopping the Mounting Process

- The machine automatically stops and relieves the pressure when the cooling time has elapsed.

## Removing the Top Closure

When the mounting process is finished:

- Turn the top closure counter-clockwise until released from the thread.
- Press RAM UP ▲ to raise the lower ram to its upper limit.
- Place the top closure in the top closure holder.

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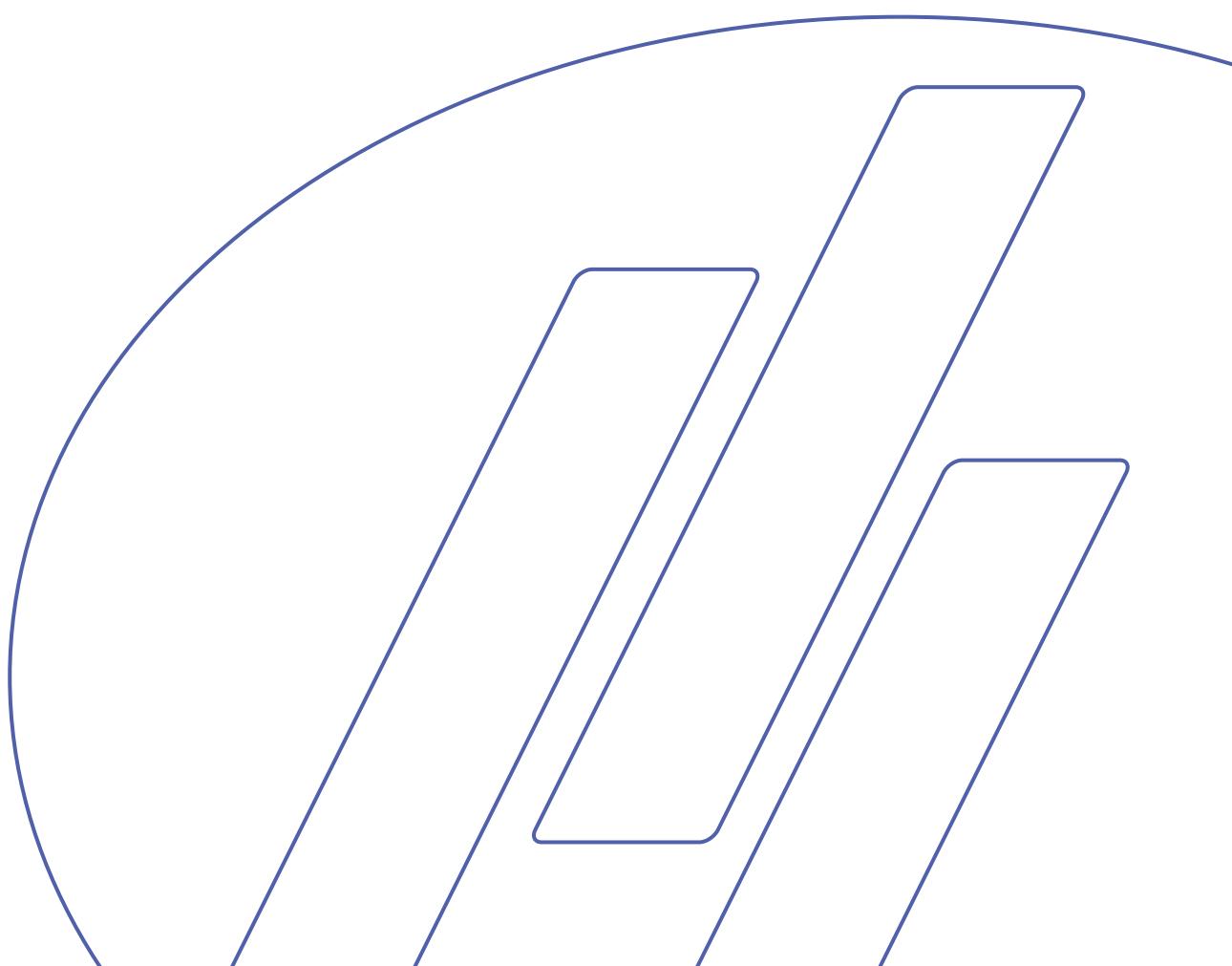
*LaboPress-1/-3  
Spare Parts and Diagrams*

---

**Always state *Serial No* and *Voltage/frequency*  
if you have technical questions or when ordering spare parts.**

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***Please note***

Some of the drawings may contain position numbers  
not used in connection with this manual.

*LaboPress-1/-3*  
*Spare Parts and Diagrams*

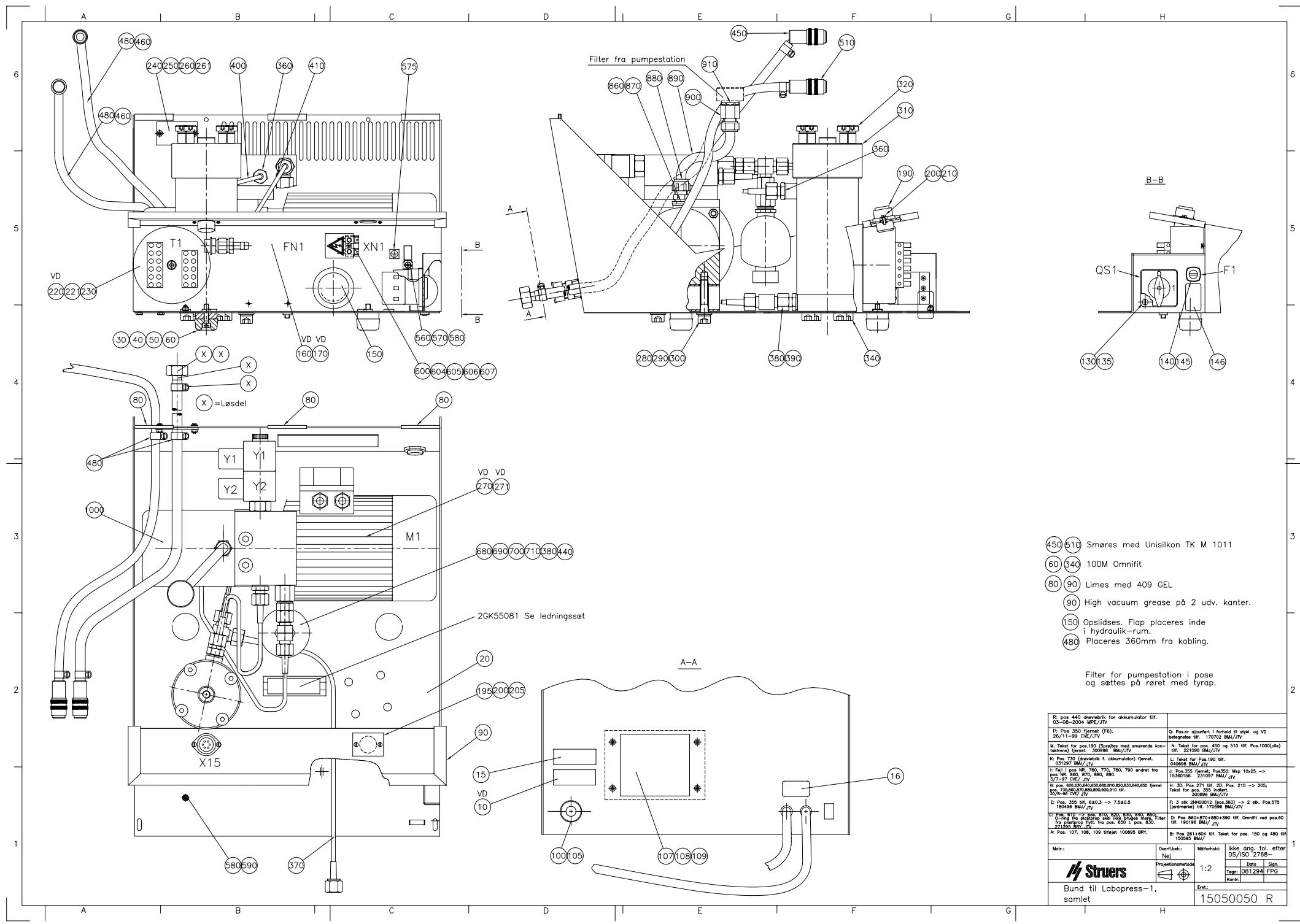
Drawing	Pos.	Spare Part	Cat No.
<b>15050050</b>		<b>Bottom, complete</b>	
	30+40	Rubber foot with screw and spacer	R4910140
	145	Fuse 3.15A T, glass tube	359MP109A
	350	Axle journal	15360228
	450	Coupling 45°	2NH11445
	510	Straight coupling with hose nipple Air Filter (lid for hydraulic oil)	2NH10014 2YF90005
<b>15050001</b>		<b>LaboPress-1, cabinet, complete</b>	
	50	Top cabinet 505	15050115
	70	Top closure holder	15050385
		<b>Various parts</b>	
		Funnel ø80/15	2GP38508
		Measuring spoon	50300000
		Allan Key 5	2GR00050
		Pipe Wrench 13mm	2GR00006
		Timer clock	62000025
		<b>LaboPress-3 Only</b>	
		Pressure water hose	2NU93020
		Gasket with filter ¾"	2IX20410
		Gasket ø11/ø24x1.5	13590359
		Reducing piece ½" ¾"	2NG30013
<b>15360036</b>		<b>Cylinder Unit ø25</b>	
	155	Cover plate ø25	15368005
	190	Fixation screw	15050040
		Lower ram, ø25	15360200
		Piston pin, for ø25 ram	15360215
		Dust protection ring	15360313
		Scraper (PROAN)	05366901

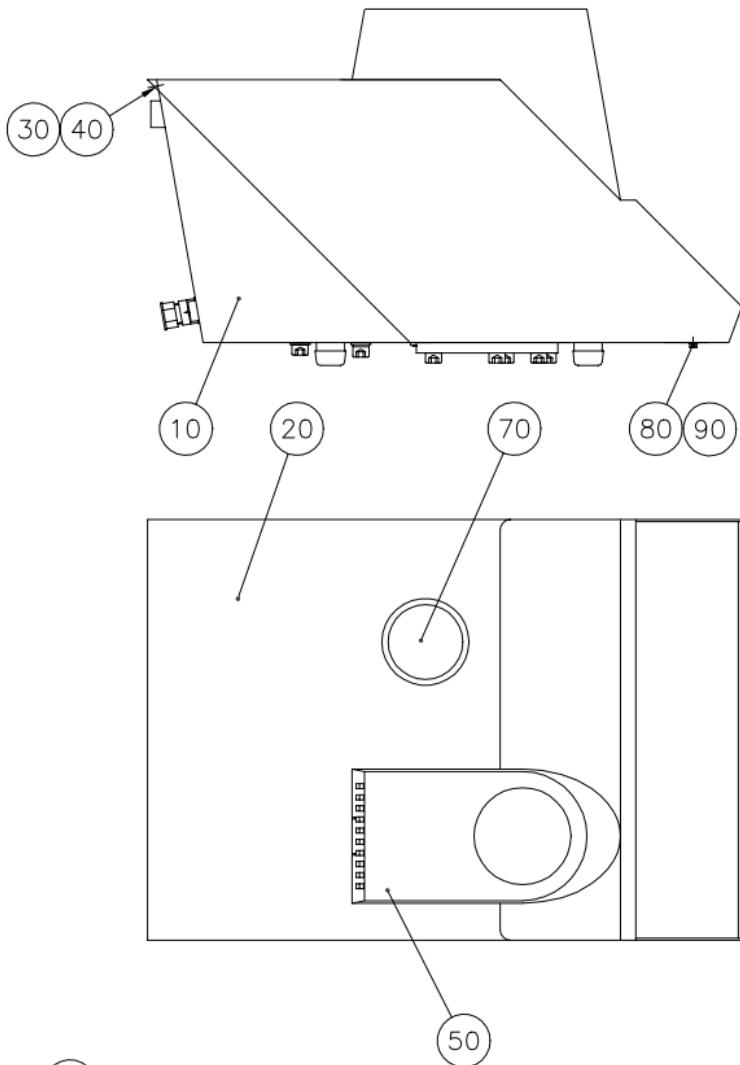
*LaboPress-1/-3*  
*Spare Parts and Diagrams*

Drawing	Pos.	Spare Part	Cat No.
<b>15360026</b>		<b>Top closure ø25</b>	
	30	Top closure ø25, complete	15360020
	30	Insulation, disc for ø25 top closure	15360300
	40	Pressure spring ø12.5x1.0	2GF10125
	50	Top part of ø25 top closure	15360170
	60	Lock washer	15050241
	70	Insulation shell for ø25 top closure	15360180
	80	Handle for top closure	15360027
<b>15360036</b>		<b>Cylinder unit ø30</b>	
	155	Cover plate ø30/1¼"	15368006
		Lower ram ø30 for ø7	15360201
		Piston pin, ø7	15360216
<b>15360026</b>		<b>Top Closure ø30</b>	
	30	Top Closure ø30, complete	15360021
	30	Insulation disc for ø30/ø1¼" top closure	15360301
	50	Top part of ø30/ø1¼" top closure	15360171
	70	Insulation shell for ø30/ø1¼" top closure	15360181
<b>15360036</b>		<b>Cylinder unit ø 1 ¼"</b>	
		Lower ram, ø7	15360202
		Piston pin, ø7	15360216
<b>15360026</b>		<b>Top closure ø 1 ¼"</b>	
		Top closure ø1¼", complete	15360022
<b>15360036</b>		<b>Cylinder unit ø 1 ½"</b>	
	155	Lower ram ø 1 ½"	15360203
	155	Cover plate ø40/1½"	15368007
		Piston pin, ø7	15360217
<b>15360026</b>		<b>Top closure ø 1 ½"</b>	
	30	Top closure ø1½", complete	15360023
	30	Insulation disc for ø40/ø1½"/ø50 top closure	15360302
	50	Top part of ø40/ø1½" top closure	15360172
	70	Insulation shell for ø40/ø1½" top closure	15360182

*LaboPress-1/-3*  
*Spare Parts and Diagrams*


Drawing	Pos.	Spare Part	Cat No.
<b>15360036</b>		<b>Cylinder unit ø40</b>	
		Lower ram ø40	15360204
		Piston pin, ø7	15360217
<b>15360026</b>		<b>Top closure ø40</b>	
		Top closure ø40, complete	15350024
<b>15360036</b>		<b>Cylinder unit ø50</b>	
	155	Cover plate ø50	15368008
		Lower ram ø50	15360205
		Piston pin, ø7	15360218
<b>15360026</b>		<b>Top closure ø50</b>	
		Top closure ø50, complete	15360025
	50	Top part of ø50 top closure	15360173
	70	Insulation shell for ø50 top closure	15360183

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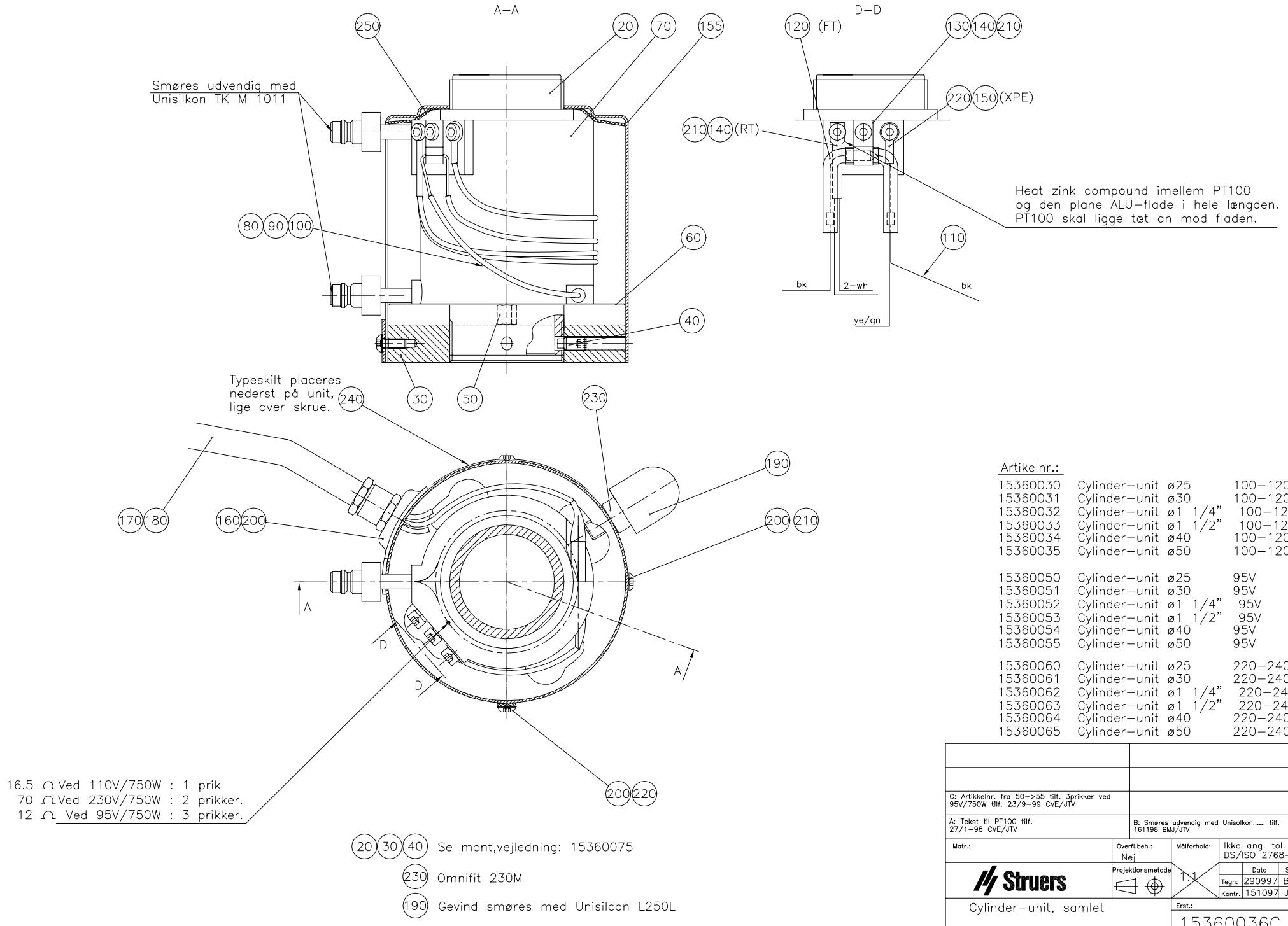
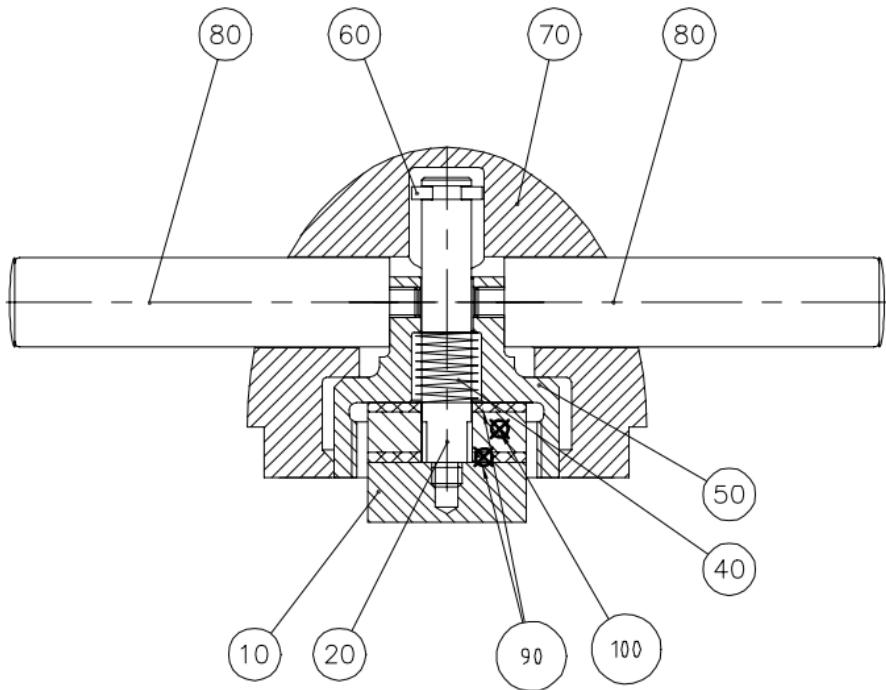
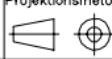
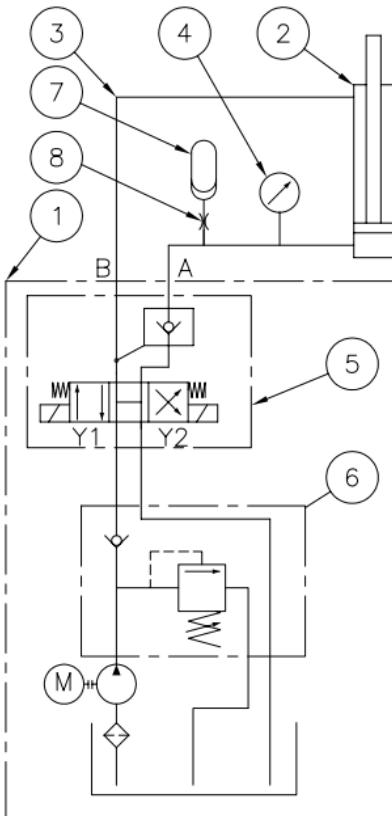
1. Hydraulic Powerpack
2. Double-acting Hydraulic Cylinder
3. ø5 Polyamid Tube
4. Pressure Gauge
5. Valve Manifold
6. Relief Valve
7. Accumulator
8. Throttle Plug

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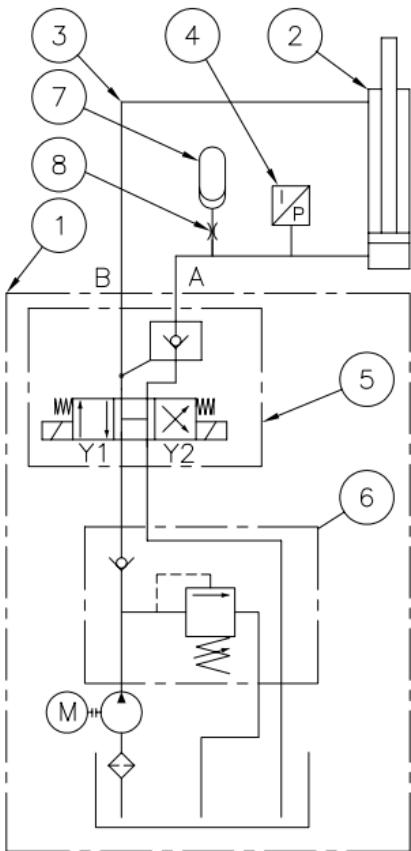
1. Hydraulic Powerpack
2. Double-acting Hydraulic Cylinder
3. ø5 Polyamid Tube
4. Pressure Transmitter
5. Valve Manifold
6. Relief Valve
7. Accumulator
8. Throttle Plug

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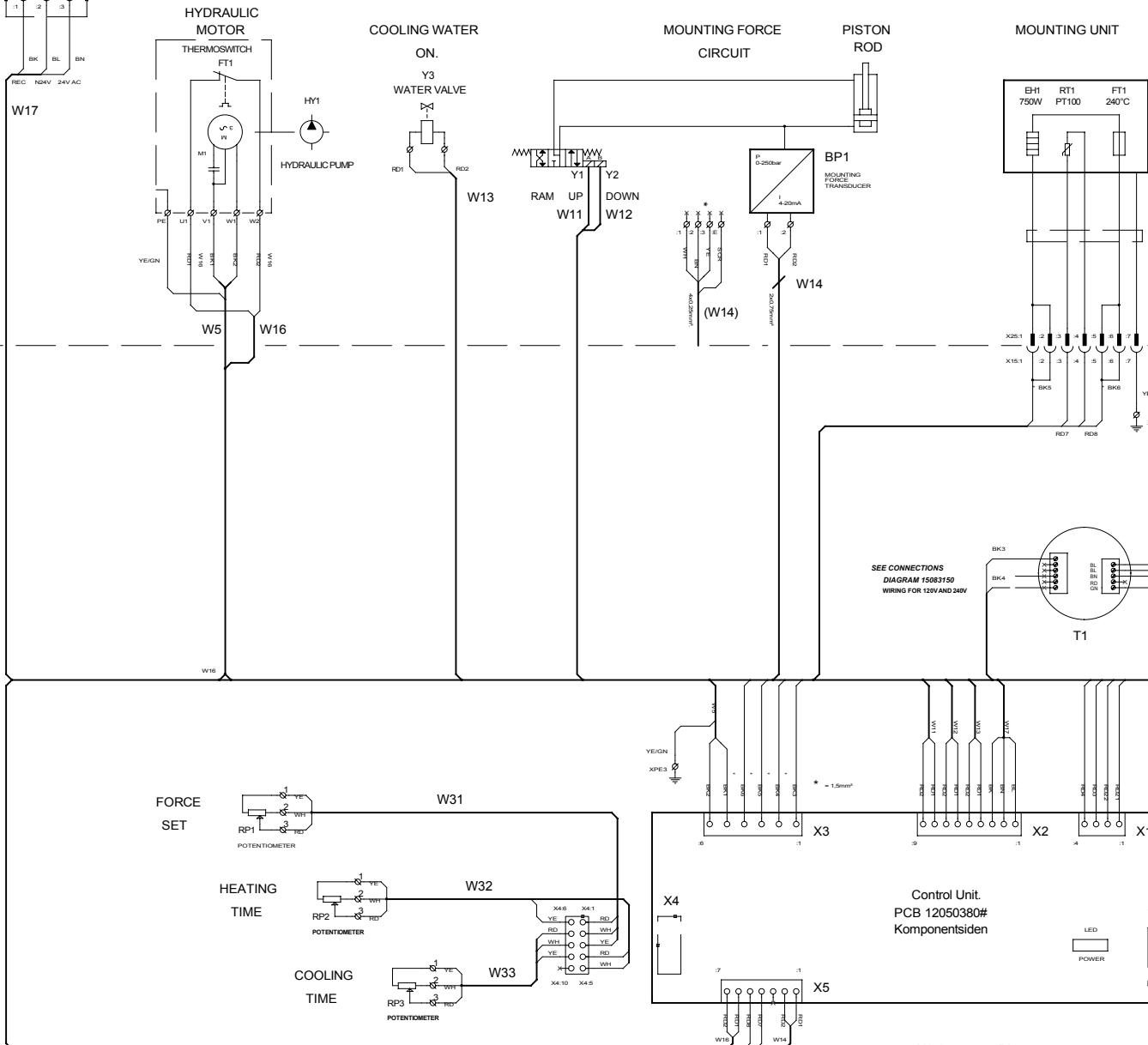Herewith declares that

<i>Product Name:</i>	LaboPress-3
<i>Type No.:</i>	508
<i>Machine Type:</i>	Mounting Apparatus

is in conformity with the provisions of the following directives:

**Safety of Machinery** 98/37/EEC according to the following standard(s):  
EN ISO 12100-1:2003, EN ISO 12100-2:2003, EN60204-1:1997.

**EMC-Directive** 2004/108/EEC according to the following standard(s):  
EN61000-6-1:2002, EN61000-6-3:2002.

**Low Voltage Directive** 2006/95/EEC according to the following standard(s):  
EN60204-1:1997.

**Supplementary Information** The equipment complies with the American standards:  
FCC part 15, subpart B and UL508.

**The above has been declared according to the global method, module A**

Date: 04.12.2007

Christian Skjold Heyde,  
Vice President, R & D and Production, Struers A/S

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Christian Skjold Heyde,  
Vice- President, R & D et Production, Struers A/S

# LaboPress-3

**Fully automatic  
mounting press for  
hot mounting of  
all materials.  
Equipped with low  
energy consuming  
mounting unit**

## Easy operation

### Setting of mounting parameters

Once the mounting parameters like: heating time, force and cooling time are set, insert the sample and resin, put on the top closure and press the start button. Your mounts will be finished in extremely short time.

## Energy saving

### New highly effective mounting units

Innovative design has merged the heating/cooling coils and the mounting cylinder into one compact and light-weight unit. This allows for a very rapid heat transfer and thus for a very fast heating and cooling. The power consumption of the heating unit could be reduced to 750 W while the efficiency compared to conventional presses could be increased by about 100 %. This means that only half the energy is sufficient to heat up the mounting unit in a far shorter time as a conventional press could do.

### Two heating levels

It is possible to set the temperature to two different levels. For mounting of heat

sensitive samples it can be set to 150 °C whereas all normal samples should be mounted at 180 °C to produce fast results.

## Different mount sizes - easy exchange

### Up to 50 mm mounts

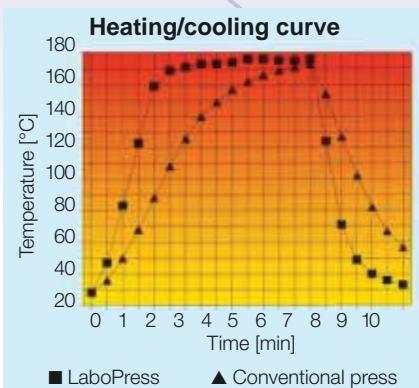
Due to the use of quick couplings for the water supply and a plug for the electrical connection, the complete mounting unit is exchanged in seconds. Six different cylinder diameters ranging from 25-50 mm and 1 1/4" - 1 1/2" are available making LaboPress-3 suitable for all possible applications.

### Intermediate ram

By using an optional intermediate ram, the capacity of the press can be doubled as two mounts can be produced at the same time.

### New top closure

The new top closure, specially developed for the new series of presses, combines the advantages of both bayonet and threaded closure.

Equipped with a triple external thread, the new closure will always "catch" like a bayonet and the threads will force the upper ram down into the cylinder, highly facilitating the work for the operator. As the threads are external, contamination with resin is highly unlikely.

## Electro-hydraulic pump

### Low energy consumption - Low noise level

The electro-hydraulic unit has an advanced motor control to stop the motor whenever the pre-set pressure is reached. This results in low energy consumption and a very low noise level.

**Struers A/S**  
Pederstrupvej 84  
DK-2750 Ballerup, Denmark  
Phone +45 44 600 800  
Fax +45 44 600 801  
struers@struers.dk  
www.struers.com

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## Technical Data

LaboPress-3		
<b>Mounting Unit</b>	Mounting dia. interchangeable	25, 30, 40, 50 mm and 1½", 1½"
	Dual mounts	Optional intermediate ram
<b>Mounting Parameters</b>	Force range, kN	0-50 Continuously adjustable
	Heating temperature, °C	150 / 180
	Heating time, min	0-15
	Cooling time, min	0-15
<b>Electronics</b>	Controls	Touch pads, knobs
	Process control for heating	Time Temperature Force
	Process control for cooling	Time Force (same as for heating)
	Automatic release of pressure	Yes
	Read out during mounting process heating temperature	Actual proces step Preset heating time, cooling time,
<b>Safety Standards</b> conforming to CE and FCC	General	89/392/EEC & 91/386/EEC & 93/44/EEC EN292-1, EN292-2, EN60204-1 (IEC 204-1)
	EMC	89/336/EEC & 92/231/EEC EN50081-2, EN50082-2 FCC eil 15, Abschnitt B
<b>Supply</b>	Electrical, no of phases	1
	Power consumption, W	
	Max total	1500
	Mounting unit	750
	Water	Tap water Min 1 bar, max 10 bar
	Recirculating cooling unit connected with ROPRE	Automatic start/stop of TRECI
<b>Dimensions and Weight</b>	Width x depth x height, mm (Mounting Unit with Top Closure installed)	420 x 580 x 420
	Weight, kg	38

## Specification

### LaboPress-3

Automatic, electro-hydraulic mounting press with pre-setting of force, heating time and time for water cooling. Mounting units in a range from 25-50 mm dia. are accommodated. The mounting unit is ordered separately

### Mounting Unit for ProntoPress-20, ProntoPress-10, LaboPress-3 and LaboPress-1

Consisting of cylinder, upper and lower ram, heating/cooling unit and top closure.

Cylinder diameter 25 mm	CYNMI
Cylinder diameter 30 mm	CYNTE
Cylinder diameter 1½"	CYNQU
Cylinder diameter 1½"	CYNHA
Cylinder diameter 40 mm	CYNIF
Cylinder diameter 50 mm	CYNAM

### Intermediate Ram

For making 2 mounts in ProntoPress-20, ProntoPress-10, LaboPress-3, LaboPress-1, PrimoPress, ProntoPress-2, PredoPress and Prestopress-3.  
Please state cylinder diameter

Struers' products are subject to constant product development. Therefore, we reserve the right to introduce changes in our products without notice.

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**Struers Ltd.**
**Struers Inc.**