

# Metallographic preparation of powder metallurgy parts

## Application Notes

Next to casting, mechanical forming and machining, powder metallurgy (P/M) technology is an important method of manufacturing metal parts. Undesirable characteristics of ingot-based metals can be greatly reduced, and desired properties of metals which would normally not alloy easily can be achieved by combining different metal powders or mixtures of metal and non metal powders.

The process of making powders, compacting them into useful shapes and then sintering them is costly, but the finished parts have some specific advantages over wrought or cast parts.

The main advantages are:



- the possibility to make fine grained homogenous structures
- the ability to form complicated shapes with close dimensional tolerances
- and the ability to produce parts with a superior surface finish.

Costly machining processes are thus reduced or eliminated and consequently there is less scrap loss compared to other forming methods. It is therefore most economical to use powder metal-

lurgy for the high volume production of small, intricately shaped, and/or very precise parts such as gears and links.

In addition, the process offers the potential to produce a wide variety of alloys with different material properties such as high temperature toughness and hardness. High speed cutting tool bits from sintered tungsten carbide powder are an example of the variety of different properties which can be achieved with the powder metallurgical process.

As the density of the compacted and sintered part influences its key properties of strength, ductility and hardness, a specific porosity is critical. For process control, metallography is used to check porosity, non-metallic inclusions and cross-contamination. In research and failure analysis, metallography is a major tool used to develop new products and improve manufacturing processes. In addition to chemical analysis, quality control also includes physical methods for checking density, dimensional changes, flow rate etc.

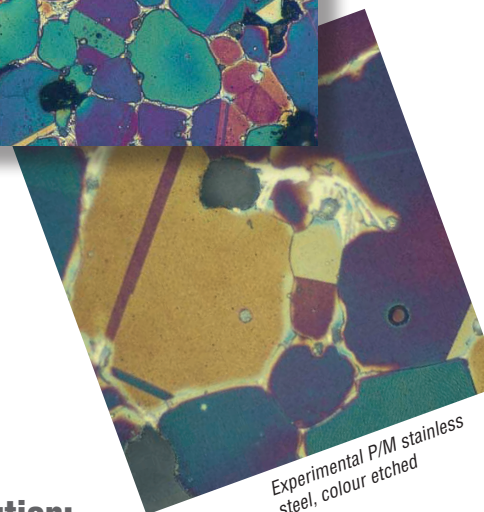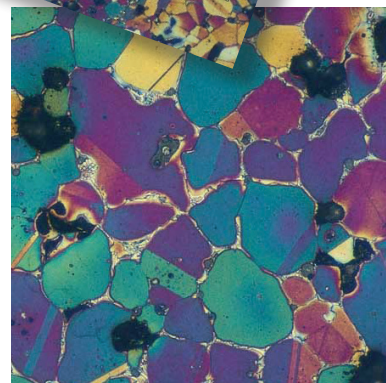

### Difficulties during metallographic preparation

#### Grinding and polishing:

To reveal the structure with the correct and representative porosity.

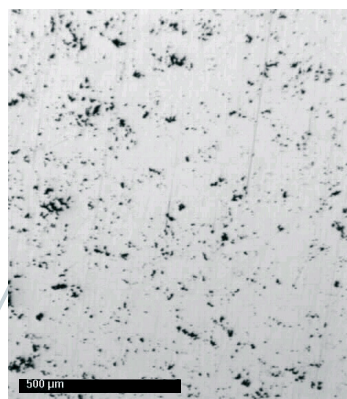


Fig. 1: Porosity of a P/M steel sample after a 4 min. polish with 3 µm diamond

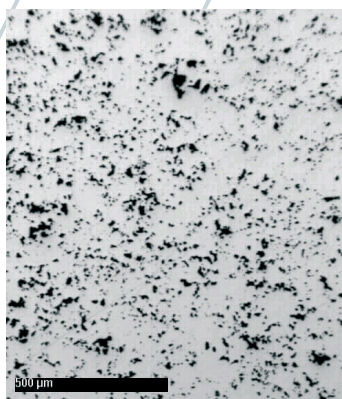


Fig. 2: Same sample as Fig. 1 after 8 min. polish with 3 µm diamond

#### Solution:

Sufficiently long polishing




Different powder metallurgy steel components

# Production of powder metallurgy parts



Synchronising hub

## Manufacturing process of iron and steel powders

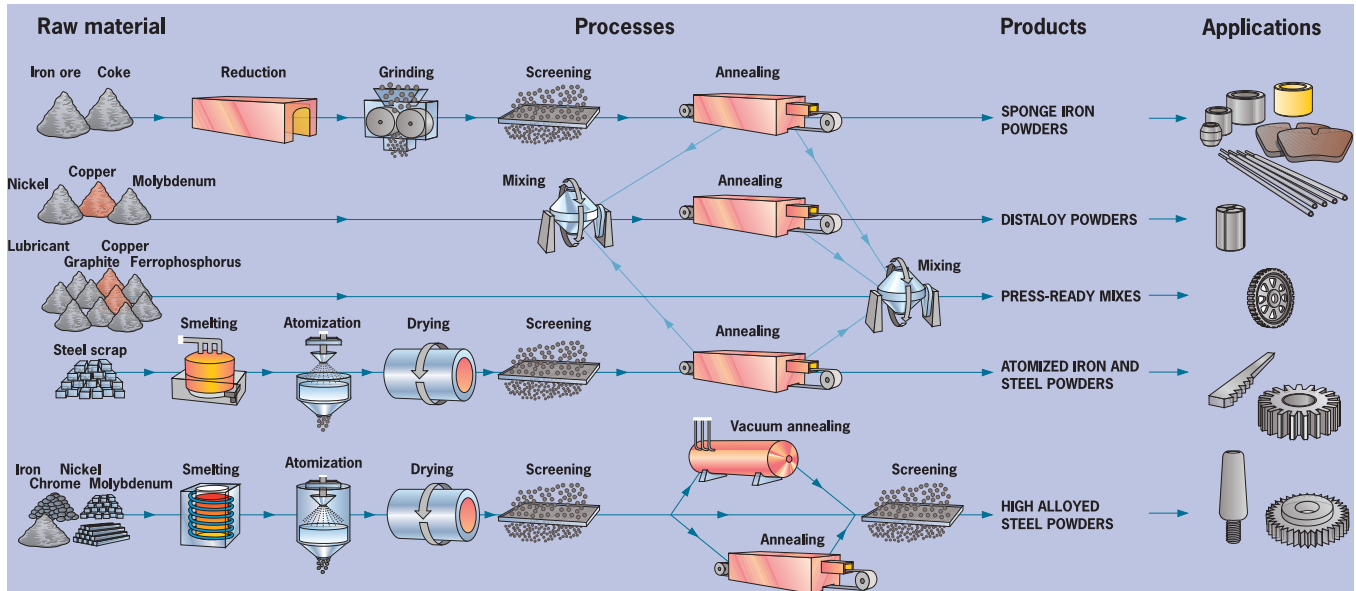

### Production

To achieve the desired structure and near net shape of a powder metal part, stringent process control of the following production steps is required:

- Making the powder
- Mixing the powder with additions such as lubricant, carbon and/or alloying elements
- Compacting the powder in carbide dies
- Sintering at high temperature (1100-1200°C) in a protective atmosphere

Chemical and atomisation methods are the two most common methods for powder production. The chemical method converts metal from ore oxides directly to metal powder at a temperature below the melting point. For example, iron powder is made through direct reduction from iron ore into sponge iron. The sponge iron is then mechanically crushed to powder, which is further refined through annealing under reducing atmosphere to

produce pure iron powder. It is used for alloying and low density applications, such as bearings.

In the atomizing process molten metal of the desired alloy flows through a nozzle and is struck by high pressure water or gas jet. Small droplets are formed which solidify into particles. Atomized powders result in higher densities than mechanically crushed powders, therefore all steel powders are produced by atomizing.

Copper powder is made by atomizing or electrolytic methods. Tungsten carbide powder is produced by adding controlled amounts of carbon to the tungsten powder and carburising it at 1400 - 2650°C


Powder production and mixing is a highly specialized and complex process which produces custom made powder mixes designed to satisfy the needs of a specific application.

A good powder mix not only has the ability to produce the required properties of a specific alloy, but also needs to facilitate handling, compacting and sintering.


For instance, the easy flow of powder and its capability to mix evenly with other powders is important for an even powder distribution before pressing, and ensures uniform properties of the finished part.

For the production of components the mixed powders are first compacted under high pressure in a carbide die. At this stage the part has the geometrical features of the finished component, but not its strength and is called the "green" part. In order to develop the mechanical and physical properties of the material, metallurgical bonding has to take place through sintering at high temperature in a sintering furnace. The bonding occurs through diffusion between adjacent particles. To avoid oxidation, which would impair the inter-particle bonding, the sintering process is conducted in a protective atmosphere.

The bonding increases the density, and pressed and sintered powder metal parts generally contain between 5 and 25% residual porosity. Depending on the application some parts may need an additional hot isostatic pressing for better dimensional accuracy, or surface finish or impregnation with oil. Final treatments, such as surface hardening, plating or coating can be applied.

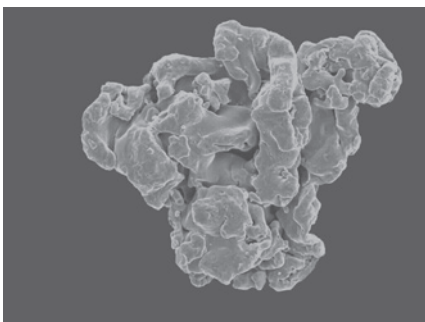

Sponge iron powder, SEM

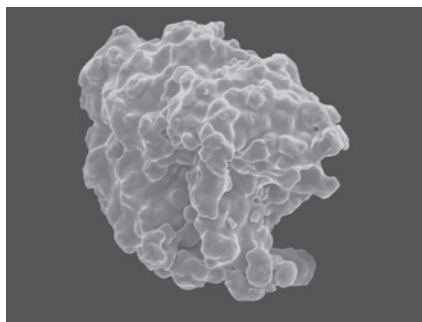

Atomized iron powder, SEM

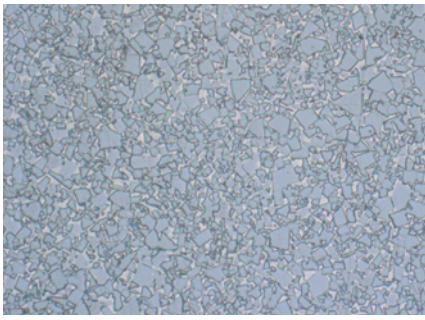


Fig. 3:  
Sintered tungsten carbide (WC/Co), etched with Murakami's reagent.  
1500 x

The sintering of cemented carbides is carried out in a vacuum sintering process. The carbide powder is mixed with 3-25 w% cobalt and small amounts of titanium and tantalum carbides are added to inhibit grain growth. This mixture is pressed and sintered. At 1280 – 1350°C the liquefaction of cobalt takes place and results in the formation of a eutectic-like phase of WC/Co. Densification begins at lower temperatures and reaches a theoretical 100% shortly after the liquefaction has occurred. During liquefaction the part shrinks up to 40% in volume.

### Applications

Components made by powder metallurgy are mainly used for the following applications:

- Mechanical and structural parts, mainly iron based, but also from copper, brass, bronze and aluminium. The largest user of P/M parts is the automotive industry. Component suppliers make connecting rods, synchronizing hubs, chain sprockets, cams and gears.
- Refractory metals which, due to their high melting points, are difficult to produce by melting and casting.
- Porous material in which controlled porosity serves a specific purpose, for instance self lubricating bearings.
- Composite materials that do not form alloys, for instance copper/tungsten for electrical contacts, cemented carbide cutting tools (Fig. 3), materials for brake linings and clutch facings, diamond cutting tools, or metal matrix composites.
- Special high-duty alloys, such as nickel and cobalt based super alloys for jet engine parts, and high speed tool steels, which have an even distribution of carbides and have isotropic qualities (Figs. 4 and 5).

In addition, different powders and powder mixtures for thermal spray coatings are produced and are also subject to metallographic quality control.

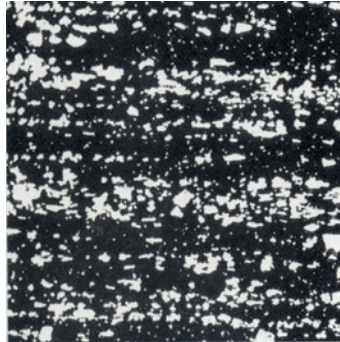


Fig. 4: Carbide distribution in conventionally produced steel

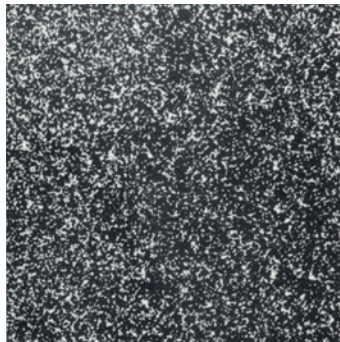


Fig. 5: Carbide distribution in powder metallurgically produced steel



## Difficulties in the metallographic preparation of powder metallurgy parts

The main challenge during the preparation is to show the true porosity after grinding and polishing. Depending on the hardness of the material this can be more or less successful. During the grinding of soft metals, abraded metal is pushed into the pores, which then has to be removed by polishing. Samples from parts in which hard and soft materials are mixed are prone to show pronounced relief. The preparation of green parts needs particular care and patience as they are very fragile.

## Recommendations for the preparation of powder metallurgy parts

### Cutting

For sectioning a powder metallurgy component of a specific metal or alloy one can select an appropriate cut-off wheel using the recommended charts and guidelines. For mixed materials it is recommended to use a cut-off wheel suitable for cutting the material which constitutes the major part. For sintered carbides a resin bonded diamond cut-off wheel is recommended (e.g. BODXX). Green parts need to be mounted in cold mounting resin (see section "Mounting") before cutting, so that they are not crushed by clamping.

### Mounting

In order to assure good adhesion of the mounting resin to the sample material it is essential to degrease the sample thoroughly with acetone, toluene or Isopar C\* before mounting (Use proper safety precautions when handling solvents!).

Sintered parts can be mounted with a resin suitable to the mechanical properties of the sample material: For hot compression mounted either phenolic resin (MultiFast) or resins containing harder filler material (DuroFast, LevoFast). For a cold mounting acrylic resins with filler (DuroCit-3 or LevoCit) can be used. For more information see [www.struers.com](http://www.struers.com) Green parts need to be re-impregnated after sectioning under vacuum with a cold mounting epoxy resin (CaldoFix-2, EpoFix, Specifix-40).

Powders can be mounted by mixing a small amount (about 1/2 teaspoon) of powder with a slow curing epoxy resin and pouring it into a mounting cup. During the approx. 12 hours curing process the particles will form a layer by settling at the bottom of the cup.

\*Available from Exxon

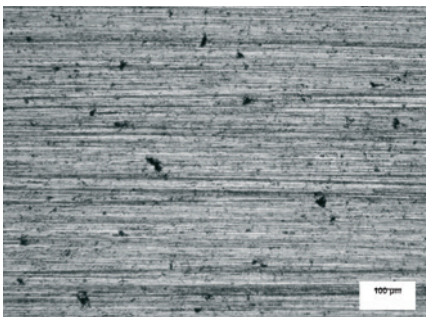


Fig. 6:  
Surface of P/M steel  
after fine grinding  
on MD-Allegro

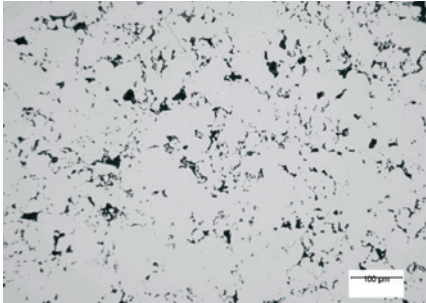


Fig. 7: Same sample as in Fig. 6 showing  
insufficient polish

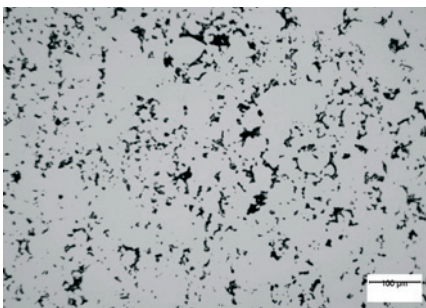


Fig. 8: Same sample as in Fig 7 after longer polish  
showing correct porosity

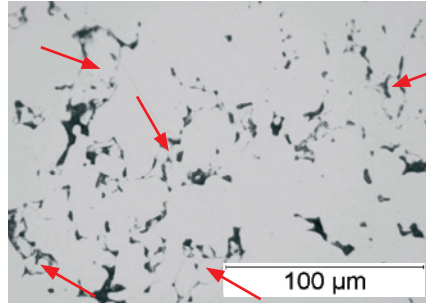


Fig. 9: Larger magnification of surface from Fig. 7,  
showing metal "lids" covering pores

of diamond polishing to remove the metal pushed into the pores during grinding and reveal the correct porosity. For soft metals the final polishing time should not be prolonged unnecessarily, as this leads to rounding of the pore edges. Starting with 500# or 800#, green parts are ground manually on silicon carbide Foil/Paper to 4000#. The surface is re-impregnated as needed. Polishing can be done on a semi-automatic polishing machine for individual samples.

In order to establish the exact polishing time for specific alloys and parts, it is recommended to check the structure during polishing every two minutes with the microscope, and only proceed to the next polishing step when all residual metal has been removed from the pores.

Hard metal powders can be hot compression mounted by mixing with one measuring spoon of the fine grained mounting resin DuroFast.

The mixture is then poured into the cylinder of the mounting press and is topped with phenolic resin.

### Grinding and polishing

The grinding and polishing procedure for powder metals follows the same rules which are applied to prepare ingot based samples of the same material.

Plane grinding high volumes of samples of materials >150 HV can be carried out on an aluminium oxide grinding stone or a diamond grinding disc (MD-Piano). Materials <150 HV can be plane ground on silicon carbide Foil/Paper.

For fine grinding with diamond the fine grinding disc MD-Allegro is suitable for materials >150 HV and MD-Largo for materials <150 HV. This is followed by a thorough diamond polish with 3 µm and a brief final polish with 1 µm or with oxide polishing suspensions. As one of the major goals of preparing a powder metallurgical sample is to show the true porosity, it is important that the diamond polishing step is carried out long enough to achieve this goal (see Figs. 6-9). For large or soft samples it can take up to 10-15 minutes

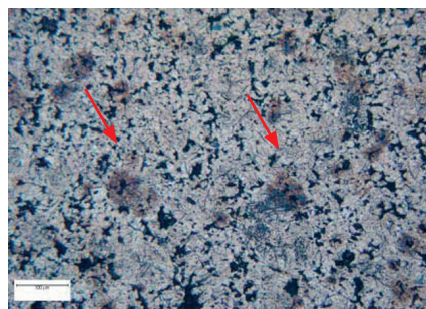


Fig.10: Water stains from cleaning can lead to  
misinterpretation of structure

### Cleaning and drying

After polishing it is essential to clean the sample with a water/detergent mixture to remove remnants of the polishing suspension and lubricant from the pores. The sample is then rinsed with water, followed by a thorough rinse with isopropanol and dried with a warm stream of air, holding the sample at an angle. Do not blow the stream of air directly

### Grinding

FG	Step	PG	
		Foil/Paper	MD-Largo
	Abrasive	Type	SiC
		Size	#320
	Suspension/ Lubricant	Water	DiaPro Allegro/Largo 9
	rpm	300	150
	Force [N]/ specimen	35	25
	Time (min)	As needed	4

### Polishing

Step	DP	OP	
	Surface	MD-Mol	
	Abrasive	Type	Diamond
		Size	3 µm
	Suspension/ Lubricant	DiaPro Mol R3	OP-S NonDry
	rpm	150	
	Force [N]/ specimen	25	
	Time (min)	4	

Table 1: Preparation method for 6 samples of P/M bronze, mounted, 30 mm dia., using the semi-automatic Tegramin, 300 mm dia.

As an alternative to DiaPro polycrystalline diamond suspension P, 9 µm, 3 µm and 1 µm can be used together with red and green or blue lubricant.

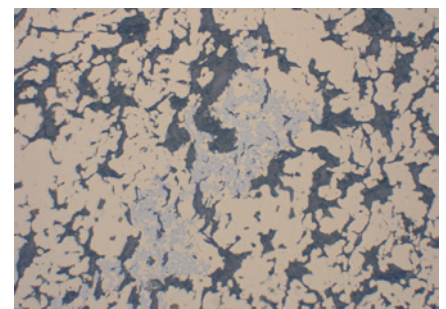


Fig.11: P/M bronze, unetched, containing  
graphite (grey), and α-δ eutectoid (blue)

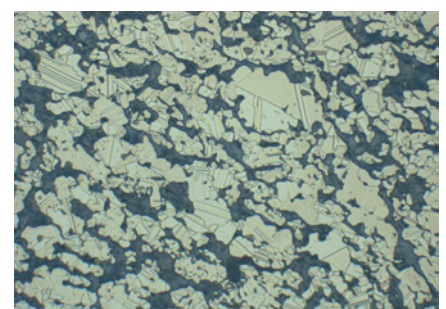










Fig.12: Same as Fig.11, etched with  
iron-III-chloride, showing grain structure  
of bronze

## Grinding

Step		PG 	FG 
	Surface	Foil/Paper	MD-Allegro
	Abrasive	Type	SiC
		Size	#220
	Suspension/ Lubricant	Water	DiaPro Allegro/Largo 9
	rpm	300	150
	Force [N]/ specimen	35	30
	Time (min)	As needed	5

## Polishing









Step		DP 1 	DP 2 
	Surface	MD-Mol*	MD-Nap
	Abrasive	Type	Diamond
		Size	3 µm
	Suspension/ Lubricant	DiaPro Mol B 3	DiaPro Nap B 1
	rpm	150	150
	Force [N]/ specimen	30	20
	Time (min)	6-8	0.5 - 1

Table 2: Preparation method for 6 P/M steel samples, mounted, 30 mm dia., using the semi-automatic Tegramin, 300 mm dia.

As an alternative to DiaPro polycrystalline diamond suspension P, 9 µm, 3 µm and 1 µm can be used together with green or blue lubricant.

\*Alternatively MD-Dac/DiaPro Dac 3 can also be used.

from the top onto the sample surface as it forces the liquids out of the pores, which will leave stains on the surface (Fig.10). It is important to use a high quality alcohol for rinsing to keep staining to a minimum.

## Etching

It is recommended to examine the unetched sample first to check the density, shape and size of the pores, oxidation and inclusions, sintering necks and free graphite (see Figs.11 and 12). It is important to know the theoretical density in order to compare it to the porosity.

After this initial inspection it is recommended to etch the sample immediately to avoid drying stains, which form when cleaning and drying liquids are gradually released by the pores. For very dense sintered carbides this is not as important as for any powder metal with a certain amount of porosity.

## Preparation method for P/M steel

To reveal the structure, the common chemical etching solutions for metals and their alloys, that are recommended in the literature, can be used. The following procedure is recommended for etching: wet the surface with isopropanol, immerse sample face up into etchant and slightly agitate. When the appropriate etching time has elapsed, take the sample out of the etchant and rinse either with isopropanol or water depending on the etchant (see below) and dry with a stream of warm air. Interpretation and photographic documentation should take place immediately after drying.

Etching time depends on the alloy and it needs some experience for the correct timing. Etching too short will not give sufficient contrast of the different phases. If the sample is over-etched, it is difficult to distinguish the various phases (see Figs.15-17). When working with unfamiliar material it is recommended to etch shorter rather than longer and to check the result with the microscope first. More etching can be carried out if necessary, but if the sample is overetched, it needs to be repolished. Following are some of the common etchants. Follow standard safety precautions when mixing and working with chemical reagents!

## Copper and copper alloys:









- 100 ml water  
20 ml hydrochloric acid  
5 g Iron-III-chloride  
10-20 sec  
Rinse with water followed by isopropanol
- 100 ml water  
10 g ammonium persulfate. Use fresh only!  
Rinse with water followed by isopropanol

## Steel:

- 1-3% Nital for iron-carbon alloys, iron-carbon-copper alloys and pre-alloyed iron-molybdenum:  
100 ml ethanol  
1-3 ml nitric acid  
10-60 sec depending on carbon content.  
Rinse with isopropanol (Fig.14)
- Picral to develop difference between martensite and austenite:  
100 ml ethanol  
4 g picric acid  
10-60 sec. depending on carbon content.  
Rinse with water followed by isopropanol (Fig.13)

## Preparation method for sintered carbides

## Grinding

Step		PG 	FG 
	Surface	MD-Piano	MD-Allegro
	Abrasive	Type	Diamond
		Size	#220
	Suspension/ Lubricant	Water	DiaPro Allegro/Largo 9
	rpm	300	150
	Force [N]/ specimen	30	30
	Time (min)	As needed	4

## Polishing


Step		DP 	OP 
	Surface	MD-Dac	MD-Chem*
	Abrasive	Type	Diamond
		Size	3 µm
	Suspension/ Lubricant	DiaPro Dac 3	OP-U NonDry
	rpm	150	150
	Force [N]/ specimen	30	20
	Time (min)	4	1

Table 3: Preparation method for 6 samples of sintered carbides, mounted, 30 mm dia., using the semi-automatic Tegramin, 300 mm dia.

As an alternative to DiaPro polycrystalline diamond suspension P, 9 µm, 3 µm and 1 µm can be used together with green/blue lubricant.

\*Optional step

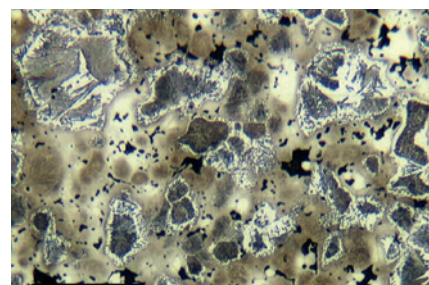


Fig.13: P/M steel with 0.5% C, diffusion alloyed with Ni, Cu and Mo. Etched in picral, showing areas of fine pearlite surrounded by ferrite, martensite, bainite and Ni-rich austenite.

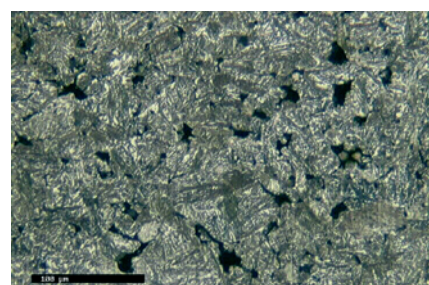


Fig.14: P/M steel with 0.8% C, pre alloyed with 1.5% Mo. Etched with nital, showing dense bainite

**Struers ApS**

 Pederstrupvej 84  
 DK-2750 Ballerup, Denmark  
 Phone +45 44 600 800  
 Fax +45 44 600 801  
 struers@struers.dk  
 www.struers.com

**NETHERLANDS**  
**Struers GmbH Nederland**  
 Zomerdijk 34 A  
 3143 CT Maassluis  
 Telefoon +31 (10) 599 7209  
 Fax +31 (10) 5997201  
 netherlands@struers.de

**NORWAY**  
**Struers ApS, Norge**  
 Sjøskogenveien 44C  
 1407 Vinterbro  
 Telefon +47 970 94 285  
 info@struers.no

**AUSTRIA**  
**Struers GmbH**  
 Zweigniederlassung Österreich  
 Betriebsgebiet Puch Nord 8  
 5412 Puch  
 Telefon +43 6245 70567  
 Fax +43 6245 70567-78  
 austria@struers.de

**POLAND**  
**Struers Sp. z o.o.**  
 Oddział w Polsce  
 ul. Jasnogórska 44  
 31-358 Kraków  
 Phone +48 12 661 20 60  
 Fax +48 12 626 01 46  
 poland@struers.de

**ROMANIA**  
**Struers GmbH, Sucursala Bucuresti**  
 Str. Preciziei nr. 6R  
 062203 sector 6, Bucuresti  
 Phone +40 (31) 101 9548  
 Fax +40 (31) 101 9549  
 romania@struers.de

**SWITZERLAND**  
**Struers GmbH**  
 Zweigniederlassung Schweiz  
 Weissenbrunnstraße 41  
 CH-8903 Birmsdorf  
 Telefon +41 44 777 63 07  
 Fax +41 44 777 63 09  
 switzerland@struers.de

**SINGAPORE**  
**Struers Singapore**  
 627A Aljunied Road,  
 #07-08 BizTech Centre  
 Singapore 389842  
 Phone +65 6299 2268  
 Fax +65 6299 2661  
 struers.sg@struers.dk

**SPAIN**  
**Struers España**  
 Camino Cerro de los Gamos 1  
 Building 1 - Pozuelo de Alarcón  
 CP 28224 Madrid  
 Teléfono +34 917 901 204  
 Fax +34 917 901 112  
 struers.es@struers.de

**FINLAND**  
**Struers ApS, Suomi**  
 Hietalahdenranta 13  
 00180 Helsinki  
 Puhelin +358 (0)207 919 430  
 Faksi +358 (0)207 919 431  
 finland@struers.fi

**SWEDEN**  
**Struers Sverige**  
 Box 20038  
 161 02 Bromma  
 Telefon +46 (0)8 447 53 90  
 Telefaks +46 (0)8 447 53 99  
 info@struers.se

**UNITED KINGDOM**  
**Struers Ltd.**  
 Unit 11 Evolution @ AMP  
 Whittle Way, Catcliffe  
 Rotherham S60 5BL  
 Tel. +44 0845 604 6664  
 Fax +44 0845 604 6651  
 info@struers.co.uk

**USA**  
**Struers Inc.**  
 24766 Detroit Road  
 Westlake, OH 44145-1598  
 Phone +1 440 871 0071  
 Fax +1 440 871 8188  
 info@struers.com

**AUSTRALIA & NEW ZEALAND**  
**Struers Australia**  
 27 Mayneview Street  
 Milton QLD 4064  
 Australia  
 Phone +61 7 3512 9600  
 Fax +61 7 3369 8200  
 info.au@struers.dk

**BELGIUM (Wallonie)**  
**Struers S.A.S.**  
 370, rue du Marché Rollay  
 F- 94507 Champigny  
 sur Marne Cedex  
 Téléphone +33 1 5509 1430  
 Télécopie +33 1 5509 1449  
 struers@struers.fr

**BELGIUM (Flanders)**  
**Struers GmbH Nederland**  
 Zomerdijk 34 A  
 3143 CT Maassluis  
 Telefoon +31 (10) 599 7209  
 Fax +31 (10) 5997201  
 netherlands@struers.de

**CANADA**  
**Struers Ltd.**  
 7275 West Credit Avenue  
 Mississauga, Ontario L5N 5M9  
 Phone +1 905-814-8855  
 Fax +1 905-814-1440  
 info@struers.com

**CHINA**  
**Struers Ltd.**  
 No. 1696 Zhang Heng Road  
 Zhang Jiang Hi-Tech Park  
 Shanghai 201203, P.R. China  
 Phone +86 (21) 6035 3900  
 Fax +86 (21) 6035 3999  
 struers@struers.cn

**CZECH REPUBLIC & SLOVAKIA**  
**Struers GmbH Organizační složka vědeckotechnický park**  
 Přílepská 1920,  
 CZ-252 63 Roztoky u Prahy  
 Phone +420 233 312 625  
 Fax +420 233 312 640  
 czechrepublic@struers.de  
 slovakia@struers.de

**GERMANY**  
**Struers GmbH**  
 Carl-Friedrich-Benz-Straße 5  
 D- 47877 Willich  
 Telefon +49 (0) 2154 486-0  
 Fax +49 (0) 2154 486-222  
 verkauf@struers.de

**FRANCE**  
**Struers S.A.S.**  
 370, rue du Marché Rollay  
 F-94507 Champigny  
 sur Marne Cedex  
 Téléphone +33 1 5509 1430  
 Télécopie +33 1 5509 1449  
 struers@struers.fr

**HUNGARY**  
**Struers GmbH**  
 Magyarországi Fióktelepe  
 2040 Budaörs  
 Szabadság utca 117  
 Phone +36 2380 6090  
 Fax +36 2380 6091  
 Email: hungary@struers.de

**IRELAND**  
**Struers Ltd.**  
 Unit 11 Evolution @ AMP  
 Whittle Way, Catcliffe  
 Rotherham S60 5BL  
 Tel. +44 0845 604 6664  
 Fax +44 0845 604 6651  
 info@struers.co.uk

**ITALY**  
**Struers Italia**  
 Via Monte Grappa 80/4  
 20020 Arese (MI)  
 Tel. +39-02/38236281  
 Fax +39-02/38236274  
 struers.it@struers.it

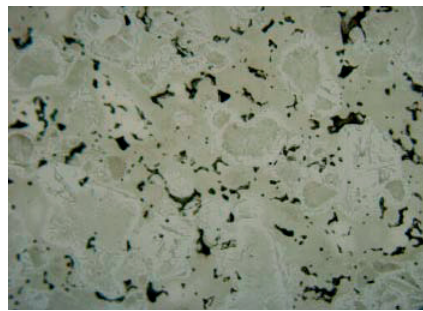
**JAPAN**  
**Marumoto Struers K.K.**  
 Takawawa Muse Bldg. 1F  
 3-14-13 Higashi-Gotanda,  
 Shinagawa  
 Tokyo  
 141-0022 Japan  
 Phone +81 3 5488 6207  
 Fax +81 3 5488 6237  
 struers@struers.co.jp


Fig.15: Etched too short

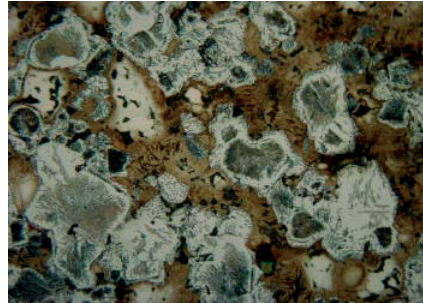


Fig.16: Etched too long, over etched

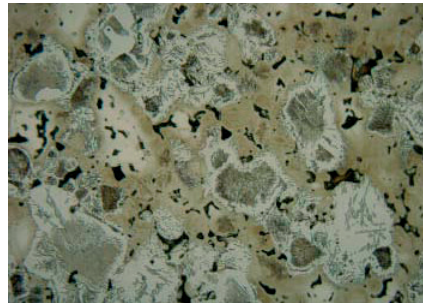


Fig.17: Correct etching

- For contrast between pearlite, bainite and martensite:  
 200 ml ethanol  
 4 g picric acid  
 1-2 ml nitric acid  
 20-100 sec, depending on carbon content and alloying elements.  
 Rinse with water followed by isopropanol

**Stainless steels:**

- Vilella's reagent:  
 45 ml glycerol  
 15 ml nitric acid  
 30 ml hydrochloric acid  
 30 sec to 5 min  
 Rinse thoroughly with water followed by isopropanol

**Sintered tungsten carbides:**

- Murakami's reagent  
 100 ml water  
 10 g potassium or sodium hydroxide  
 10 g potassium ferricyanide  
 Etching by immersion or swab etch  
 Rinse thoroughly with water followed by isopropanol

**Summary**

Powder metallurgy is a method of producing parts from metals which normally would not alloy easily, or from mixtures of metals and non-metals, taking advantage of their combined properties. Although the powder metallurgical process is costly, it has the advantage of economically producing large amounts of small and intricately shaped parts with homogenous structures. The density of compacted and sintered parts affects their strength, ductility and hardness. Therefore metallographic control of porosity is an integral part of quality control. During metallographic grinding, metal is pushed into the pores and if the subsequent polishing steps are not carried out properly, residual metal "lids" covering the pores will obstruct the evaluation of the correct porosity. Thorough grinding and polishing with diamond, and microscopic checks between polishing steps, assure a true representation of the structure. The given procedures for automatic preparation and the chemical etchants have been used successfully in day to day practical laboratory application and give reproducible results.

**Author**


Elisabeth Weidmann, Birgitte Nielsen,  
 Struers A/S, Copenhagen, Denmark  
 Judy Arner, Struers, Inc., Westlake, OH, USA

**Acknowledgements**

We wish to thank Höganäs AB, SE 26383 Höganäs, Sweden, for the co-operation and supplying information and material for reproduction: all photos of P/M components, diagram and SEM photos on page 2, Figs. 1, 2, 6, 7, 8, 9, 10, and 13-17. Special thanks go to M. Carlsson who generously shared her professional experience with us.  
 Figs. 4 and 5 courtesy Böhler Edelstahl GmbH, Kapfenberg, Austria

**Bibliography**

- Höganäs Handbook for sintered components, 6. Metallography, Höganäs AB, 1999  
 Höganäs iron and steel powders for sintered components, Höganäs AB, 1998  
 Structure 24, Metallographic examination of P/M stainless steel matrix composites, N. Frees, Institute for Product Development, Technical University, Lyngby, Denmark, 1991  
 Metallographic etching, G. Petzow, ASM, Metals Park, Ohio, 1978

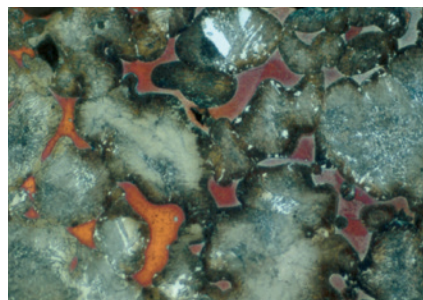

P/M steel with copper infiltration